



# Agilent 1200 Infinity Series Variable Wavelength Detectors

User Manual



**Agilent Technologies**

# Notices

© Agilent Technologies, Inc. 2008, 2010-2012, 2013

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

## Manual Part Number

G1314-90034 Rev B

## Edition

11/2013

Printed in Germany

Agilent Technologies  
Hewlett-Packard-Strasse 8  
76337 Waldbronn

**This product may be used as a component of an in vitro diagnostic system if the system is registered with the appropriate authorities and complies with the relevant regulations. Otherwise, it is intended only for general laboratory use.**

## Warranty

**The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.**

## Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

## Restricted Rights Legend

If software is for use in the performance of a U.S. Government prime contract or subcontract, Software is delivered and licensed as “Commercial computer software” as defined in DFAR 252.227-7014 (June 1995), or as a “commercial item” as defined in FAR 2.101(a) or as “Restricted computer software” as defined in FAR 52.227-19 (June 1987) or any equivalent agency regulation or contract clause. Use, duplication or disclosure of Software is subject to Agilent Technologies’ standard commercial license terms, and non-DOD Departments and Agencies of the U.S. Government will

receive no greater than Restricted Rights as defined in FAR 52.227-19(c)(1-2) (June 1987). U.S. Government users will receive no greater than Limited Rights as defined in FAR 52.227-14 (June 1987) or DFAR 252.227-7015 (b)(2) (November 1995), as applicable in any technical data.

## Safety Notices

### CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

### WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

# In This Guide

This manual covers

- the Agilent 1290 Infinity Variable Wavelength Detector (G1314E),
- the Agilent 1260 Infinity Variable Wavelength Detector (G1314F) and
- the Agilent 1200 series Variable Wavelength Detector (G1314D) (obsolete).

Find information on other Agilent Variable Wavelength Detectors in separate manuals.

## **1 Introduction to the Variable Wavelength Detector**

This chapter gives an introduction to the detector, instrument overview and internal connectors.

## **2 Site Requirements and Specifications**

This chapter gives information on environmental requirements, physical and performance specifications.

## **3 Installing the Detector**

This chapter describes the installation of the detector.

## **4 Using the Detector**

This chapter provides information on how to set up the detector for an analysis and explains the basic settings.

## **5 How to optimize the detector**

This chapter gives hints on how to select the detector parameters and the flow cell.

## **6 Troubleshooting and Diagnostics**

Overview about the troubleshooting and diagnostic features.

### **7 Error Information**

This chapter describes the meaning of detector error messages, and provides information on probable causes and suggested actions how to recover from error conditions.

### **8 Test Functions**

This chapter describes the detector's built in test functions.

### **9 Maintenance and Repair**

This chapter provides general information on maintenance and repair of the detector.

### **10 Parts and Materials for Maintenance**

This chapter provides information on parts for maintenance.

### **11 Identifying Cables**

This chapter provides information on cables used with the Agilent 1200 Infinity Series modules.

### **12 Hardware Information**

This chapter describes the detector in more detail on hardware and electronics.

### **13 LAN Configuration**

This chapter provides information on connecting the detector to the Agilent ChemStation PC.

### **14 Appendix**

This chapter provides addition information on safety, legal and web.

# Contents


<b>1</b>	<b>Introduction to the Variable Wavelength Detector</b>	<b>9</b>
	Introduction to the Detector	10
	Optical System Overview	11
	System Overview	16
<b>2</b>	<b>Site Requirements and Specifications</b>	<b>19</b>
	Site Requirements	20
	Physical Specifications	23
	Performance Specifications G1314D	24
	Performance Specifications G1314E	27
	Performance Specifications G1314F	30
<b>3</b>	<b>Installing the Detector</b>	<b>33</b>
	Unpacking the Detector	34
	Optimizing the Stack Configuration	36
	Installation Information on Leak and Waste Handling	45
	Installing the Detector	49
	Flow Connections to the Detector	52
<b>4</b>	<b>Using the Detector</b>	<b>55</b>
	Leak and Waste Handling	56
	Setting up an Analysis	57
	Special Settings of the Detector	71
<b>5</b>	<b>How to optimize the detector</b>	<b>83</b>
	Optimizing the Detector Performance	84
	Match the Flow Cell to the Column	85
	Set the Detector Parameters	88

<b>6</b>	<b>Troubleshooting and Diagnostics</b>	<b>89</b>
	Overview of the Detector's Indicators and Test Functions	90
	Status Indicators	91
	Available Tests versus Interfaces	93
	Agilent Lab Advisor Software	94
<b>7</b>	<b>Error Information</b>	<b>95</b>
	What Are Error Messages	97
	General Error Messages	98
	Detector Error Messages	104
<b>8</b>	<b>Test Functions</b>	<b>115</b>
	Intensity Test	116
	Cell Test	118
	Wavelength Verification-Calibration	120
	ASTM Drift and Noise Test	122
	Quick Noise Test	123
	Dark Current Test	124
	Holmium Oxide Test	126
<b>9</b>	<b>Maintenance and Repair</b>	<b>129</b>
	Introduction to Maintenance	130
	Warnings and Cautions	131
	Overview of Maintenance	133
	Cleaning the Module	134
	Exchanging a Lamp	135
	Exchanging a Flow Cell	138
	Repairing the Flow Cells	140
	Using the Cuvette Holder	143
	Correcting Leaks	145
	Replacing Leak Handling System Parts	146
	Replacing the Module's Firmware	147

<b>10</b>	<b>Parts and Materials for Maintenance</b>	<b>149</b>
	Overview of Maintenance Parts	150
	Standard Flow Cell 10 mm / 14 $\mu$ L	152
	Micro Flow Cell 3 mm / 2 $\mu$ L	154
	Semi-micro Flow Cell 6 mm / 5 $\mu$ L	156
	High Pressure Flow Cell 10 mm / 14 $\mu$ L	158
	Cuvette Holder	160
	Kits	151
	Leak Parts	161
<b>11</b>	<b>Identifying Cables</b>	<b>163</b>
	Cable Overview	164
	Analog Cables	166
	Remote Cables	168
	BCD Cables	171
	CAN/LAN Cables	173
	RS-232 Cables	174
<b>12</b>	<b>Hardware Information</b>	<b>175</b>
	Firmware Description	176
	Electrical Connections	179
	Interfaces	182
	Setting the 8-bit Configuration Switch	189
	Instrument Layout	193
	Early Maintenance Feedback (EMF)	194
<b>13</b>	<b>LAN Configuration</b>	<b>197</b>
	What you have to do first	198
	TCP-IP parameter configuration	199
	Configuration Switches	200
	Initialization mode selection	201
	Dynamic Host Configuration Protocol (DHCP)	205
	Link configuration selection	208
	Automatic configuration with Bootp	209
	Manual Configuration	219
	PC and User Interface Software Setup	225

**14 Appendix 227**

General Safety Information	228
Radio Interference	231
Sound Emission	232
UV Radiation	233
Solvent Information	234
Declaration of Conformity for HOX2 Filter	236
Agilent Technologies on Internet	237

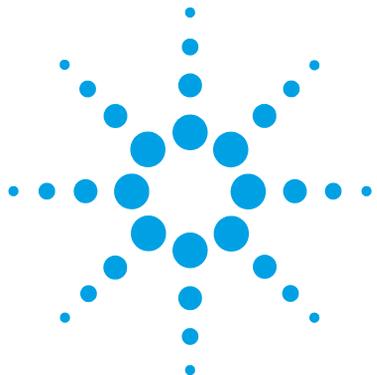

# 1 Introduction to the Variable Wavelength Detector

Introduction to the Detector	10
Optical System Overview	11
Flow Cell	12
Lamp	13
Source Lens Assembly	13
Entrance Slit Assembly	13
Filter Assembly	14
Mirror Assemblies M1 and M2	15
Grating Assembly	15
Beam Splitter Assembly	15
Photo Diodes Assemblies	15
Photo Diode ADC (analog-to-digital converter)	15
System Overview	16
Leak and Waste Handling	16

This chapter gives an introduction to the detector, instrument overview and internal connectors.



## Introduction to the Detector


The Agilent variable wavelength detectors described in this manual are designed for highest optical performance, GLP compliance and easy maintenance with:

- data rate up to 20 Hz for standard-HPLC (G1314D), see [Table 10](#) on page 78
- higher data rate up to 80 Hz for fast-HPLC (G1314F), see [Table 11](#) on page 78
- higher data rate up to 160 Hz for ultra-fast-HPLC (G1314E), see [Table 12](#) on page 79
- data recovery card DRC (G1314E) gives unique "data-never-lost" insurance "[Run Recovery Settings](#)" on page 80,
- deuterium lamp for highest intensity and lowest detection limit over a wavelength range of 190 to 600 nm,
- optional flow-cell cartridges (standard 10 mm, 14  $\mu$ L; high pressure 10 mm, 14  $\mu$ L; micro 3 mm, 2  $\mu$ L; semi-micro 6 mm, 5  $\mu$ L) are available and can be used depending on the application needs (other types may be introduced later),
- easy front access to lamp and flow cell for fast replacement,
- electronic identification of flow cell and lamp with RFID (Radio Frequency Identification) tag for unambiguous identification,
  - lamp information: part number, serial number, production date, ignitions, burn time
  - cell information: part number, serial number, production date, nominal path length, volume, maximum pressure
- built-in electronic temperature control (ETC) for improved baseline stability, and
- built-in holmium oxide filter for fast wavelength accuracy verification.

For specifications refer to [Table 3](#) on page 24.

## Optical System Overview

The optical system of the detector is shown in the figure below. Its radiation source is a deuterium-arc discharge lamp for the ultraviolet (UV) wavelength range from 190 to 600 nm. The light beam from the deuterium lamp passes through a lens, a filter assembly, an entrance slit, a spherical mirror (M1), a grating, a second spherical mirror (M2), a beam splitter, and finally through a flow cell to the sample diode. The beam through the flow cell is absorbed depending on the solutions in the cell, in which UV absorption takes place, and the intensity is converted to an electrical signal by means of the sample photodiode. Part of the light is directed to the reference photodiode by the beam splitter to obtain a reference signal for compensation of intensity fluctuation of the light source. A slit in front of the reference photodiode cuts out light of the sample bandwidth. Wavelength selection is made by rotating the grating, which is driven directly by a stepper motor. This configuration allows fast change of the wavelength. The cutoff filter is moved into the lightpath above 370 nm to reduce higher order light.

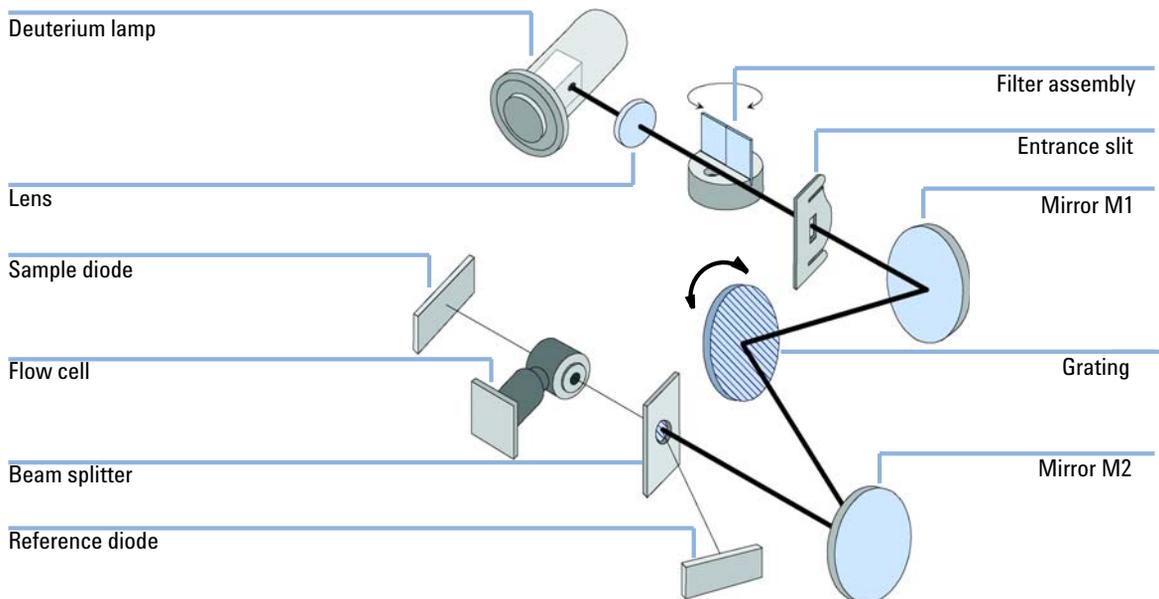


**Figure 1** Optical Path of the Variable Wavelength Detector

## Flow Cell

A variety of flow-cell cartridges can be inserted using the same quick and simple mounting system.

The flow cells have an integrated RFID tag that contains the flow cell specific information (e.g. part number, cell volume, path length, ...). A RFID tag reader reads out this information and transfers it to the user interface.

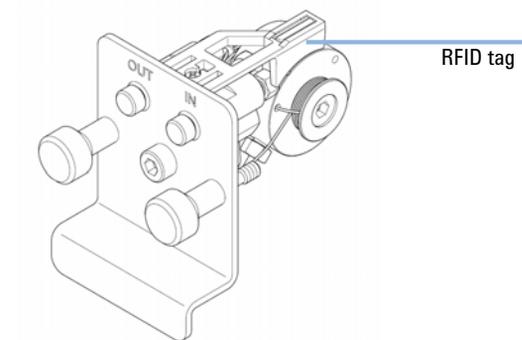

**Figure 2** Flow Cell with RFID tag

**Table 1** Flow Cell Data

	<b>STD</b>	<b>Semi-micro</b>	<b>Micro</b>	<b>High Pressure</b>	
Maximum pressure	40 (4)	40 (4)	120 (12)	400 (40)	bar
Path length	10 (conical)	6 (conical)	3 (conical)	10 (conical)	mm
Volume	14	5	2	14	μL
Inlet i.d.	0.25	0.17	0.12	0.25	mm
Inlet length	750	250	310	750	mm
Outlet i.d.	0.30	0.17	0.17	0.17	mm
Outlet length	120	120	120	120	mm
Total volume	60.77	14.49	14.00	60.77	μL
Materials in contact with solvent	SST, quartz, PTFE, PEEK	SST, quartz, PTFE	SST, quartz, PTFE	SST, quartz, Kapton	

## Lamp

The light source for the UV wavelength range is a deuterium lamp. As a result of plasma discharge in a low pressure deuterium gas, the lamp emits light over the 190 – 600 nm wavelength range.

The lamp has an integrated RFID tag that contains the lamp specific information (e.g. part number, burn time, ...). A RFID tag reader reads out this information and transfers it to the user interface.

## Source Lens Assembly


The source lens receives the light from the deuterium lamp and focuses it onto the entrance slit.

## Entrance Slit Assembly

The entrance slit assembly has an exchangeable slit. The standard one has a 1-mm slit. For replacement and calibration purposes to optimize the alignment, a slit with a hole is needed.

## Filter Assembly

The filter assembly is electromechanically actuated. During wavelength calibrations it moves into the light path.

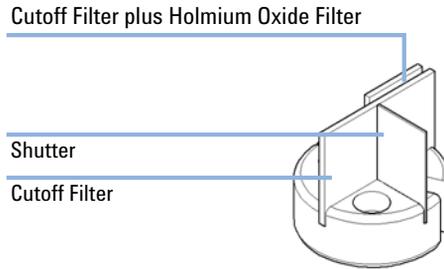

**Figure 3** Filter Assembly

The filter assembly has two filters installed and is processor-controlled.

<b>OPEN</b>	nothing in light path at $\lambda < 370$ nm
<b>CUTOFF</b>	cut off filter in light path at $\lambda > 370$ nm
<b>HOLMIUM</b>	holmium oxide filter for wavelength check
<b>M</b>	
<b>SHUTTER</b>	for measurement of dark current of photo diodes

A photo sensor determines the correct position.

## Mirror Assemblies M1 and M2

The instrument contains two spherical mirrors (M1 and M2). The beam adjustable is vertically and horizontally. Both mirrors are identical.

## Grating Assembly

The grating separates the light beam into all its component wavelengths and reflects the light onto mirror #2.

The stepper motor reference position is determined by a plate fitted onto the motor shaft, interrupting the beam of a photo sensor. The wavelength calibration of the grating is done at the zero order light position and at 656 nm, which is the emission line of the deuterium lamp.

## Beam Splitter Assembly

The beam splitter splits the light beam. One part goes directly to the sample diode. The other part of the light beam goes to the reference diode.

## Photo Diodes Assemblies

Two photo diode assemblies are installed in the optical unit. The sample diode assembly is located on the left side of the optical unit. The reference diode assembly is located in the front of the optical unit.


## Photo Diode ADC (analog-to-digital converter)

The photo diode current is directly converted to digital data direct photo current digitalization. The data is transferred to the detector main board . The photo diode ADC boards are located close to the photo diodes.

## **System Overview**

### **Leak and Waste Handling**

The 1200 Infinity Series has been designed for safe leak and waste handling. It is important that all security concepts are understood and instructions are carefully followed.

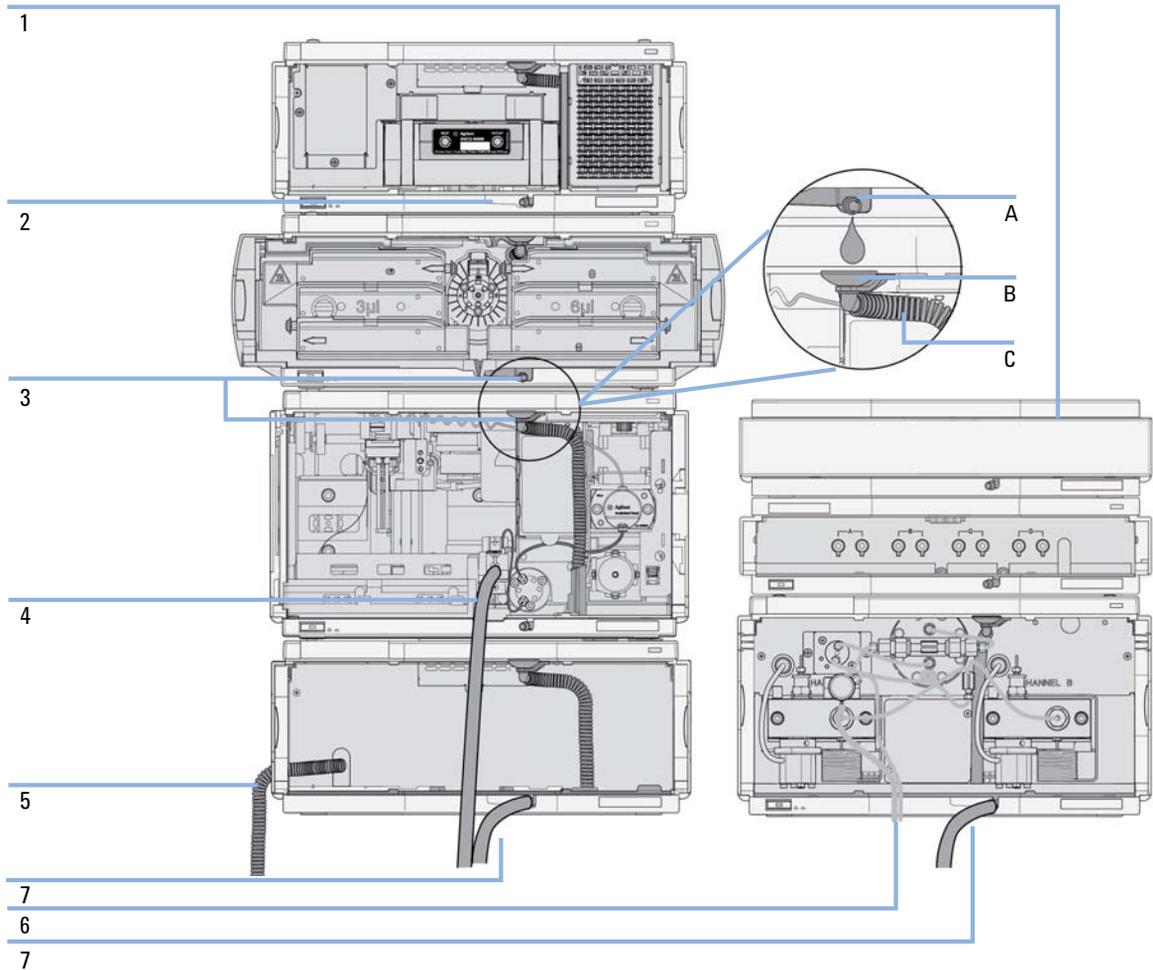

**Figure 4** Leak and waste handling concept (overview - typical stack configuration as an example)

The solvent cabinet (1) is designed to store a maximum volume of 6 L solvent. The maximum volume for an individual bottle stored in the solvent cabinet should not exceed 2.5 L. For details, see the usage guideline for the Agilent 1200 Infinity Series Solvent Cabinets (a printed copy of the guideline has been shipped with the solvent cabinet, electronic copies are available on the Internet).

## 1 Introduction to the Variable Wavelength Detector

### System Overview

The leak pan (2) (individually designed in each module) guides solvents to the front of the module. The concept covers also leakages on internal parts (e.g. the detector's flow cell). The leak sensor in the leak pan stops the running system as soon as the leak detection level is reached.


The leak pan's outlet port (3, A) guides excessive overflow from one module to the next, as the solvent flows into the next module's leak funnel (3, B) and the connected corrugated waste tube (3, C). The corrugated waste tube guides the solvent to the next lower positioned module's leak tray and sensor.

The waste tube of the sampler's needle wash port (4) guides solvents to waste.

The condense drain outlet of the autosampler cooler (5) guides condensate to waste.

The waste tube of the purge valve (6) guides solvents to waste.

The waste tube connected to the leak pan outlet on each of the bottom instruments (7) guides the solvent to a suitable waste container.

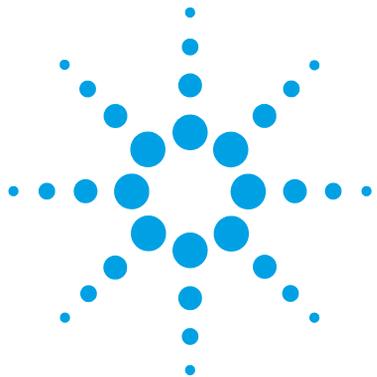

## 2 Site Requirements and Specifications

Site Requirements	20
Physical Specifications	23
Performance Specifications G1314D	24
Performance Specifications G1314E	27
Performance Specifications G1314F	30

This chapter gives information on environmental requirements, physical and performance specifications.



## Site Requirements

A suitable environment is important to ensure optimal performance of the instrument.

### Power Consideration

The detector power supply has wide ranging capabilities, see “[Physical Specifications](#)” on page 23. It accepts any line voltage in the above mentioned range. Consequently, there is no voltage selector in the rear of the detector. There are also no externally accessible fuses, because automatic electronic fuses are implemented in the power supply.

#### WARNING

##### **Instrument is partially energized when switched off**

**The power supply still uses some power, even when the power switch on the front panel is turned OFF. Repair work at the detector can lead to personal injuries, e. g. shock hazard, when the detector cover is opened and the instrument is connected to power.**

→ To disconnect the detector from the power line, unplug the power cord.

---

#### WARNING

##### **Hazard of electrical shock or damage of your instrumentation**

**can result, if the devices are connected to a line voltage higher than specified.**

→ Connect your instrument to the specified line voltage only.

---

#### CAUTION

Inaccessible power plug.

In case of emergency it must be possible to disconnect the instrument from the power line at any time.

→ Make sure the power connector of the instrument can be easily reached and unplugged.

→ Provide sufficient space behind the power socket of the instrument to unplug the cable.

---

## Power Cords

Different power cords are offered as options with the module. The female end of all power cords is identical. It plugs into the power-input socket at the rear. The male end of each power cord is different and designed to match the wall socket of a particular country or region.

### **WARNING**

#### **Absence of ground connection or use of unspecified power cord**

**The absence of ground connection or the use of unspecified power cord can lead to electric shock or short circuit.**

- Never operate your instrumentation from a power outlet that has no ground connection.
  - Never use a power cord other than the Agilent Technologies power cord designed for your region.
- 

### **WARNING**

#### **Use of unsupplied cables**

**Using cables not supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.**

- Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.
- 

### **WARNING**

#### **Unintended use of supplied power cords**

**Using power cords for unintended purposes can lead to personal injury or damage of electronic equipment.**

- Never use the power cords that Agilent Technologies supplies with this instrument for any other equipment.
-

## Bench Space

The detector dimensions and weight (see “Physical Specifications” on page 23) allows you to place the detector on almost any desk or laboratory bench. It needs an additional 2.5 cm (1.0 inch) of space on either side and approximately 8 cm (3.1 inch) in the rear for air circulation and electric connections.

If the bench should carry an Agilent 1200 Infinity Series system, make sure that the bench is designed to bear the weight of all modules.

The detector should be operated in a horizontal position.

## Environment

Your detector will work within specifications at ambient temperatures and relative humidity as described in “Physical Specifications” on page 23.

Better drift performance depends on better control of the temperature fluctuations. To realize the highest performance, minimize the frequency and the amplitude of the temperature changes to below 1 °C/hour (1.8 °F/hour). Turbulences around one minute or less can be ignored.

### CAUTION

Condensation within the module

Condensation will damage the system electronics.

- Do not store, ship or use your module under conditions where temperature fluctuations could cause condensation within the module.
  - If your module was shipped in cold weather, leave it in its box and allow it to warm slowly to room temperature to avoid condensation.
-

## Physical Specifications

**Table 2** Physical Specifications

Type	Specification	Comments
Weight	11 kg (25 lbs)	
Dimensions (height × width × depth)	140 x 345 x 435 mm (5.5 x 13.5 x 17 inches)	
Line voltage	100 – 240 VAC, ± 10 %	Wide-ranging capability
Line frequency	50 or 60 Hz, ± 5 %	
Power consumption	220 VA, 85 W / 290 BTU	Maximum
Ambient operating temperature	0–55 °C (32–131 °F)	
Ambient non-operating temperature	-40 – 70 °C (-40 – 158 °F)	
Humidity	< 95 % r.h. at 40 °C (104 °F)	Non-condensing
Operating altitude	Up to 2000 m (6562 ft)	
Non-operating altitude	Up to 4600 m (15091 ft)	For storing the module
Safety standards: IEC, CSA, UL	Installation category II, Pollution degree 2	For indoor use only.

## Performance Specifications G1314D

### Performance Specifications G1314D

**Table 3** Performance Specifications

Type	Specification	Comments
Detection type	Double-beam photometer	
Light source	Deuterium lamp	
Wavelength range	190 – 600 nm	The UV-lamp is equipped with RFID tag that holds lamp typical information.
Short term noise	$\pm 0.15 \cdot 10^{-5}$ AU at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314D</a> ” on page 26 below the table.
Drift	$< 1 \cdot 10^{-4}$ AU/h at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314D</a> ” on page 26 below the table.
Linearity	$> 2.5$ AU (5 %) at 265 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314D</a> ” on page 26 below the table.
Wavelength accuracy	$\pm 1$ nm	Self-calibration with deuterium lines, verification with holmium oxide filter
Maximum sampling rate	20 Hz	
Band width	6.5 nm typical	
Flow cells	Standard: 14 $\mu$ L volume, 10 mm cell path length and 40 bar (588 psi) pressure maximum High pressure: 14 $\mu$ L volume, cell path length and 400 bar (5880 psi) pressure maximum Micro: 2 $\mu$ L volume, 3 mm cell path length and 120 bar (1760 psi) pressure maximum Semi-micro: 5 $\mu$ L volume, 6 mm cell path length and 40 bar (588 psi) pressure maximum	All flow cells have RFID tags for unambiguous identification. Can be repaired on component level

**Table 3** Performance Specifications

Type	Specification	Comments
Electronic Temperature Control (ETC)	For improved baseline stability in instable environment.	
Control and data evaluation	Agilent ChemStation B.03.02 SR1 or above Instant Pilot (G4208A) with firmware B.02.07 or above	Control and data evaluation Control only
Time programmable	Wavelength, Reference and Sample scan, balance, steps, lamp on/off	
Spectral tools	Stop-flow wavelength scan	
Analog outputs	Recorder/integrator: 100 mV or 1 V, output range 0.001 – 2 AU, one output	
Communications	LAN card integrated on main board, Controller-area network (CAN), RS-232C, APG Remote: ready, start, stop and shut-down signals	
Safety and maintenance	Extensive diagnostics, error detection and display (through Instant Pilot and Data System), leak detection, safe leak handling, leak output signal for shutdown of pumping system. Low voltages in major maintenance areas.	
GLP features	Early maintenance feedback (EMF) for continuous tracking of instrument usage in terms of lamp burn time with user-settable limits and feedback messages. Electronic records of maintenance and errors. Verification of wavelength accuracy with built-in holmium oxide filter. RFID for electronics records of flow cell and UV lamp conditions (path length, volume, product number, serial number, test passed, usage)	
Housing	All materials recyclable.	

## Specification Conditions G1314D

ASTM: “Standard Practice for Variable Wavelength Photometric Detectors Used in Liquid Chromatography”.

Reference conditions: Standard flow cell, path length 10 mm, flow 1 mL/min LC-grade methanol.

**Noise:**

$\pm 0.15 \cdot 10^{-5}$  AU at 230 nm, TC 2 s

RT = 2.2 \* TC

**Linearity:**

Linearity is measured with caffeine at 265 nm.

**NOTE**

The specification are based on the the standard RFID tag lamp (G1314-60101) and may be not achieved when other lamp types or aged lamps are used.

ASTM drift tests require a temperature change below 2°C/hour (3.6°F/hour) over one hour period. Our published drift specification is based on these conditions. Larger ambient temperature changes will result in larger drift.

Better drift performance depends on better control of the temperature fluctuations. To realize the highest performance, minimize the frequency and the amplitude of the temperature changes to below 1°C/hour (1.8°F/hour). Turbulences around one minute or less can be ignored.

Performance tests should be done with a completely warmed up optical unit (> one hour). ASTM measurements require that the detector should be turned on at least 24 hours before start of testing.

## Performance Specifications G1314E

**Table 4** Performance Specifications G1314E

Type	Specification	Comments
Detection type	Double-beam photometer	
Light source	Deuterium lamp	
Wavelength range	190 – 600 nm	The UV-lamp is equipped with RFID tag that holds lamp typical information.
Short term noise	$\pm 0.15 \cdot 10^{-5}$ AU at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314E</a> ” on page 29 below the table.
Drift	$< 1 \cdot 10^{-4}$ AU/h at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314E</a> ” on page 29 below the table.
Linearity	$> 2.5$ AU (5 %) at 265 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314E</a> ” on page 29 below the table.
Wavelength accuracy	$\pm 1$ nm	Self-calibration with deuterium lines, verification with holmium oxide filter
Maximum data rate	160 Hz	
Band width	6.5 nm typical	
Flow cells	Standard: 14 $\mu$ L volume, 10 mm cell path length and 40 bar (588 psi) pressure maximum High pressure: 14 $\mu$ L volume, cell path length and 400 bar (5880 psi) pressure maximum Micro: 2 $\mu$ L volume, 3 mm cell path length and 120 bar (1760 psi) pressure maximum Semi-micro: 5 $\mu$ L volume, 6 mm cell path length and 40 bar (588 psi) pressure maximum	All flow cells have RFID tags for unambiguous identification. Can be repaired on component level
Electronic Temperature Control (ETC)	For improved baseline stability in instable environment.	

## 2 Site Requirements and Specifications

### Performance Specifications G1314E

**Table 4** Performance Specifications G1314E

Type	Specification	Comments
Control and data evaluation	Agilent ChemStation B.03.02 SR1 or above Instant Pilot (G4208A) with firmware B.02.07 or above	Control and data evaluation Control only
Time programmable	Wavelength, Reference and Sample scan, balance, steps, lamp on/off	
Spectral tools	Stop-flow wavelength scan	
Analog outputs	Recorder/integrator: 100 mV or 1 V, output range 0.001 – 2 AU, one output	
Communications	LAN card integrated on main board, Controller-area network (CAN), RS-232C, APG Remote: ready, start, stop and shut-down signals	
Safety and maintenance	Extensive diagnostics, error detection and display (through Instant Pilot and Data System), leak detection, safe leak handling, leak output signal for shutdown of pumping system. Low voltages in major maintenance areas.	
GLP features	Early maintenance feedback (EMF) for continuous tracking of instrument usage in terms of lamp burn time with user-settable limits and feedback messages. Electronic records of maintenance and errors. Verification of wavelength accuracy with built-in holmium oxide filter. RFID for electronics records of flow cell and UV lamp conditions (path length, volume, product number, serial number, test passed, usage)	
Housing	All materials recyclable.	

## Specification Conditions G1314E

ASTM: “Standard Practice for Variable Wavelength Photometric Detectors Used in Liquid Chromatography”.

Reference conditions: Standard flow cell, path length 10 mm, flow 1 mL/min LC-grade methanol.

**Noise:**

$\pm 0.15 \cdot 10^{-5}$  AU at 230 nm, TC 2 s

RT = 2.2 \* TC

**Linearity:**

Linearity is measured with caffeine at 265 nm.

**NOTE**

The specification are based on the the standard RFID tag lamp (G1314-60101) and may be not achieved when other lamp types or aged lamps are used.

ASTM drift tests require a temperature change below 2°C/hour (3.6°F/hour) over one hour period. Our published drift specification is based on these conditions. Larger ambient temperature changes will result in larger drift.

Better drift performance depends on better control of the temperature fluctuations. To realize the highest performance, minimize the frequency and the amplitude of the temperature changes to below 1°C/hour (1.8°F/hour). Turbulences around one minute or less can be ignored.

Performance tests should be done with a completely warmed up optical unit (> one hour). ASTM measurements require that the detector should be turned on at least 24 hours before start of testing.

## Performance Specifications G1314F

**Table 5** Performance Specifications G1314F

Type	Specification	Comments
Detection type	Double-beam photometer	
Light source	Deuterium lamp	
Wavelength range	190 – 600 nm	The UV-lamp is equipped with RFID tag that holds lamp typical information.
Short term noise	$\pm 0.25 \cdot 10^{-5}$ AU at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314F</a> ” on page 32 below the table.
Drift	$< 1 \cdot 10^{-4}$ AU/h at 230 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314F</a> ” on page 32 below the table.
Linearity	$> 2.5$ AU (5 %) at 265 nm	Under specified conditions. See “ <a href="#">Specification Conditions G1314F</a> ” on page 32 below the table.
Wavelength accuracy	$\pm 1$ nm	Self-calibration with deuterium lines, verification with holmium oxide filter
Maximum data rate	80 Hz	
Band width	6.5 nm typical	
Flow cells	Standard: 14 $\mu$ L volume, 10 mm cell path length and 40 bar (588 psi) pressure maximum High pressure: 14 $\mu$ L volume, cell path length and 400 bar (5880 psi) pressure maximum Micro: 2 $\mu$ L volume, 3 mm cell path length and 120 bar (1760 psi) pressure maximum Semi-micro: 5 $\mu$ L volume, 6 mm cell path length and 40 bar (588 psi) pressure maximum	All flow cells have RFID tags for unambiguous identification. Can be repaired on component level
Electronic Temperature Control (ETC)	For improved baseline stability in instable environment.	

**Table 5** Performance Specifications G1314F

Type	Specification	Comments
Control and data evaluation	Agilent ChemStation B.04.02 SP2 or above Instant Pilot (G4208A) with firmware B.02.11 or above	Control and data evaluation Control only
Time programmable	Wavelength, Reference and Sample scan, balance, steps, lamp on/off	
Spectral tools	Stop-flow wavelength scan	
Analog outputs	Recorder/integrator: 100 mV or 1 V, output range 0.001 – 2 AU, one output	
Communications	LAN card integrated on main board, Controller-area network (CAN), RS-232C, APG Remote: ready, start, stop and shut-down signals	
Safety and maintenance	Extensive support for troubleshooting and maintenance is provided by the Instant Pilot, Agilent Lab Advisor, and the Chromatography Data System. Safety-related features are leak detection, safe leak handling, leak output signal for shutdown of pumping system, and low voltages in major maintenance areas.	
GLP features	Early maintenance feedback (EMF) for continuous tracking of instrument usage in terms of lamp burn time with user-settable limits and feedback messages. Electronic records of maintenance and errors. Verification of wavelength accuracy with built-in holmium oxide filter. RFID for electronics records of flow cell and UV lamp conditions (path length, volume, product number, serial number, test passed, usage)	
Housing	All materials recyclable.	

## Specification Conditions G1314F

ASTM: “Standard Practice for Variable Wavelength Photometric Detectors Used in Liquid Chromatography”.

Reference conditions: Standard flow cell, path length 10 mm, flow 1 mL/min LC-grade methanol.

**Noise:**

$\pm 0.25 \cdot 10^{-5}$  AU at 230 nm, TC 2 s

RT = 2.2 \* TC

**Linearity:**

Linearity is measured with caffeine at 265 nm.


**NOTE**

The specification are based on the the standard RFID tag lamp (G1314-60101) and may be not achieved when other lamp types or aged lamps are used.

ASTM drift tests require a temperature change below 2°C/hour (3.6°F/hour) over one hour period. Our published drift specification is based on these conditions. Larger ambient temperature changes will result in larger drift.

Better drift performance depends on better control of the temperature fluctuations. To realize the highest performance, minimize the frequency and the amplitude of the temperature changes to below 1°C/hour (1.8°F/hour). Turbulences around one minute or less can be ignored.

Performance tests should be done with a completely warmed up optical unit (> one hour). ASTM measurements require that the detector should be turned on at least 24 hours before start of testing.

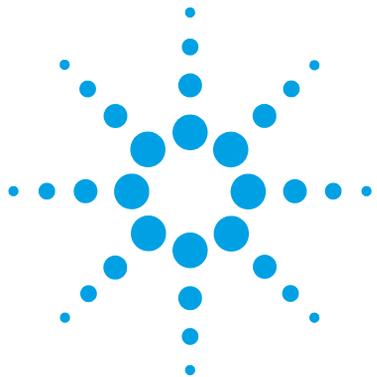

## 3 Installing the Detector

Unpacking the Detector	34
Damaged Packaging	34
Delivery Checklist	35
Detector Accessory Kit Contents	35
Optimizing the Stack Configuration	36
One Stack Configuration	37
Two Stack Configuration	41
Installation Information on Leak and Waste Handling	45
Installing the Detector	49
Flow Connections to the Detector	52

This chapter describes the installation of the detector.



## Unpacking the Detector

### CAUTION

Condensation within the detector

Condensation will damage the system electronics.

- Do not store, ship or use your detector under conditions where temperature fluctuations could cause condensation within the detector.
  - If your detector was shipped in cold weather, leave it in its box and allow it to warm up slowly to room temperature to avoid condensation.
- 

## Damaged Packaging

If the delivery packaging shows signs of external damage, please call your Agilent Technologies sales and service office immediately. Inform your service representative that the instrument may have been damaged during shipment.

### CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
  - An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.
-

## Delivery Checklist

Ensure all parts and materials have been delivered with the detector. The delivery checklist is shown below. Please report missing or damaged parts to your local Agilent Technologies sales and service office.

**Table 6** Variable Wavelength Detector Checklist

Description	Quantity
Variable wavelength detector	1
Power cable	1
Flow cell	As ordered
<i>User Manual</i> on Documentation CD (part of the shipment - not module specific)	1 per order
Accessory kit	1
CompactFlash Card (G1314E)	1

## Detector Accessory Kit Contents

The module is shipped with Accessory kit (G1314-68755) (see [“Accessory Kit”](#) on page 151).


## **Optimizing the Stack Configuration**

If your detector is part of a complete Agilent 1200 Infinity Series system, you can ensure optimum performance by installing the following configuration. This configuration optimizes the system flow path, ensuring minimum delay volume.

## One Stack Configuration

### One Stack Configuration for Agilent 1260 Infinity LC


Ensure optimum performance by installing the modules of the Agilent 1260 Infinity LC System in the following configuration (See [Figure 5](#) on page 37 and [Figure 6](#) on page 38). This configuration optimizes the flow path for minimum delay volume and minimizes the bench space required.

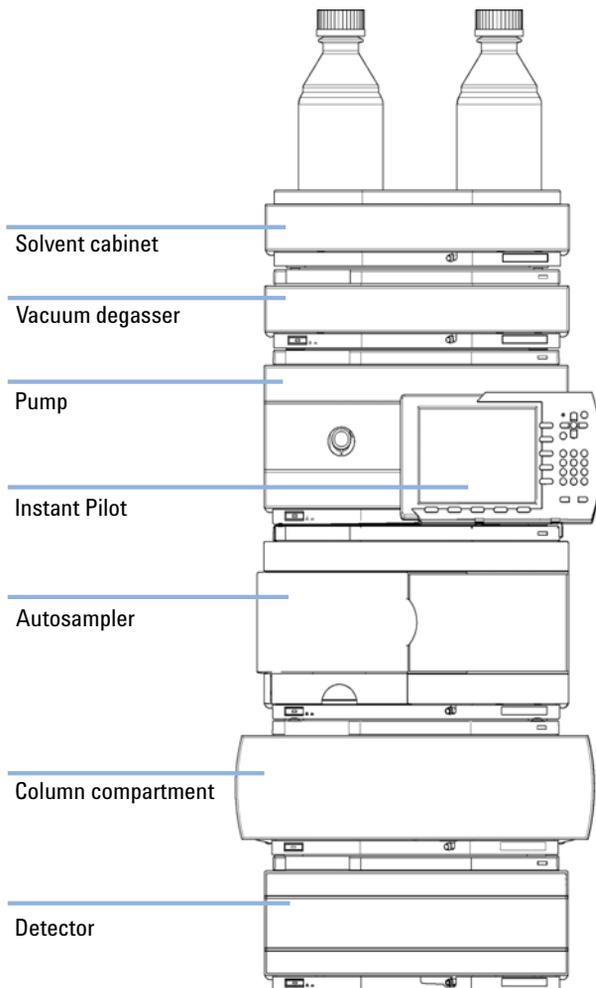

**Figure 5** Recommended Stack Configuration for 1260 Infinity (Front View)

### 3 Installing the Detector

#### Optimizing the Stack Configuration

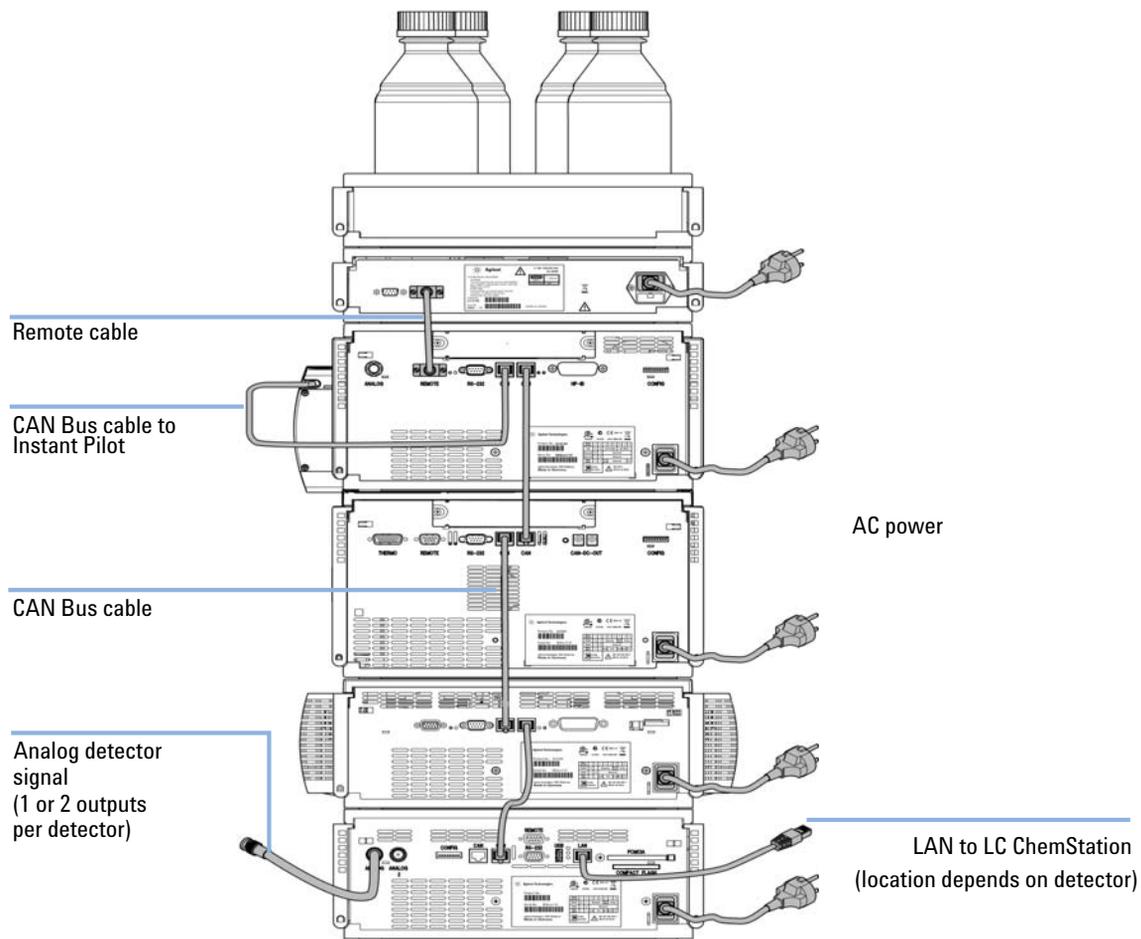


**Figure 6** Recommended Stack Configuration for 1260 Infinity (Rear View)

### One Stack Configuration for Agilent 1290 Infinity LC

Ensure optimum performance by installing the modules of the Agilent 1290 Infinity Binary LC System in the following configuration (See [Figure 7](#) on page 39 and [Figure 8](#) on page 40). This configuration optimizes the flow path for minimum delay volume and minimizes the bench space required.


The Agilent 1290 Infinity Binary Pump should always be installed at the bottom of the stack.

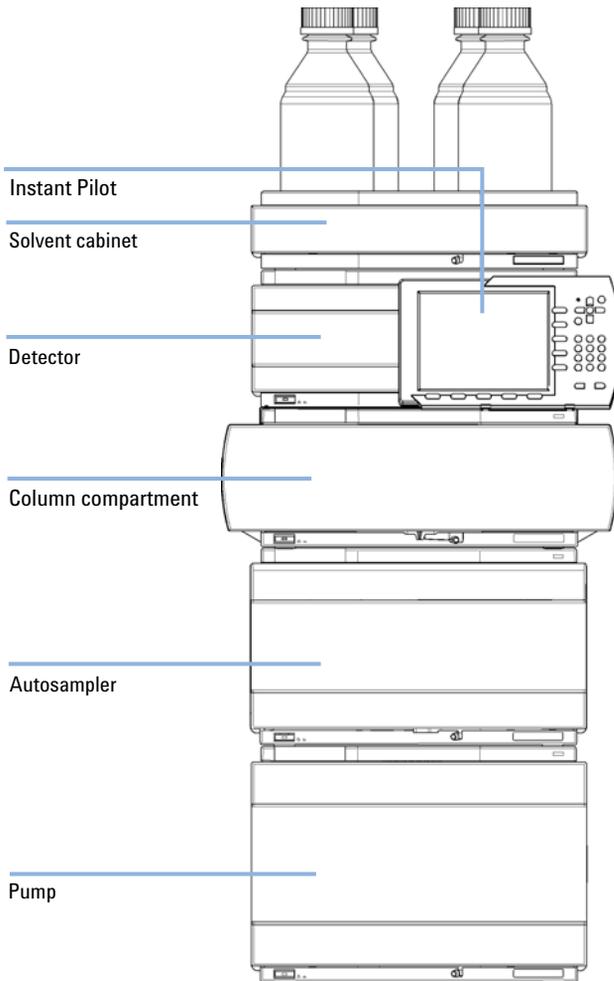

**Figure 7** Recommended stack configuration for 1290 Infinity with binary pump (front view)

### 3 Installing the Detector

#### Optimizing the Stack Configuration

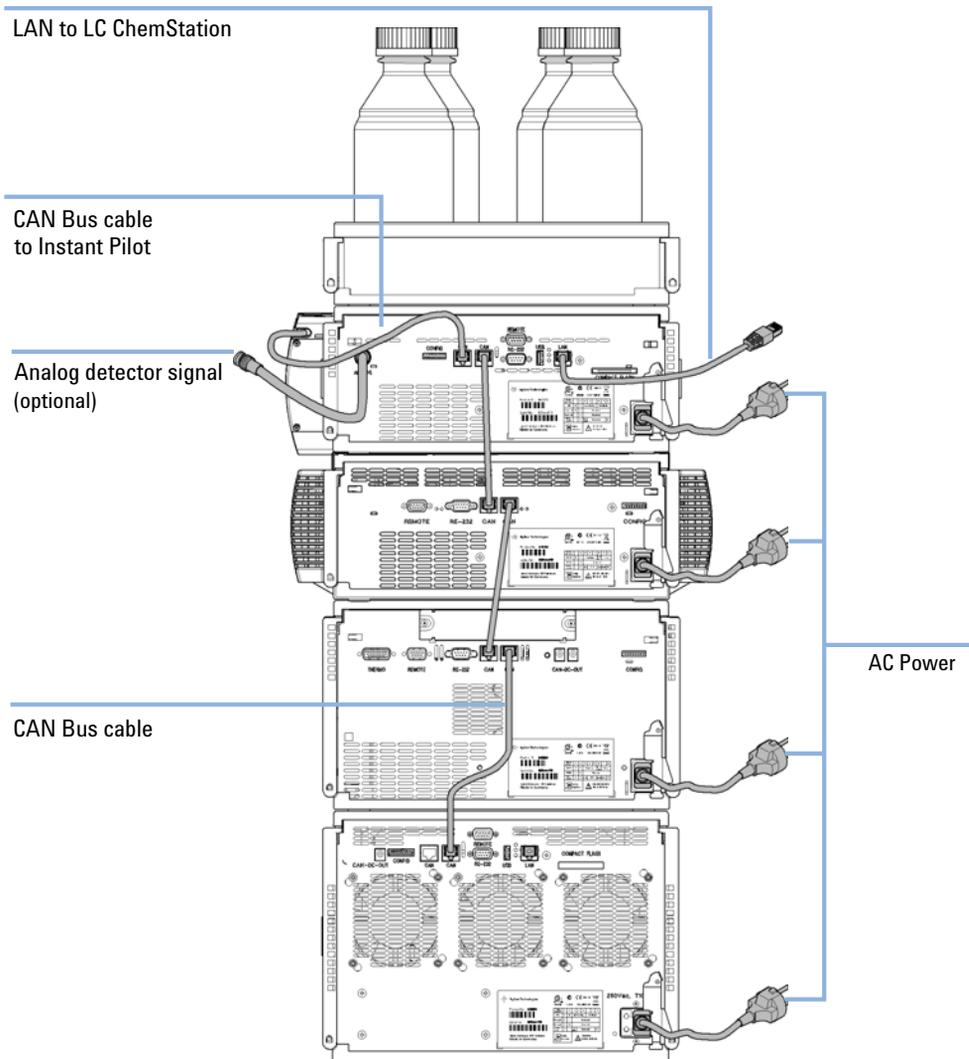


**Figure 8** Recommended stack configuration 1290 Infinity with binary pump (rear view)

## Two Stack Configuration

### Two Stack Configuration for Agilent 1260 Infinity LC


To avoid excessive height of the stack when the autosampler thermostat is added to the system it is recommended to form two stacks. Some users prefer the lower height of this arrangement even without the autosampler thermostat. A slightly longer capillary is required between the pump and autosampler. (See [Figure 9](#) on page 41 and [Figure 10](#) on page 42).

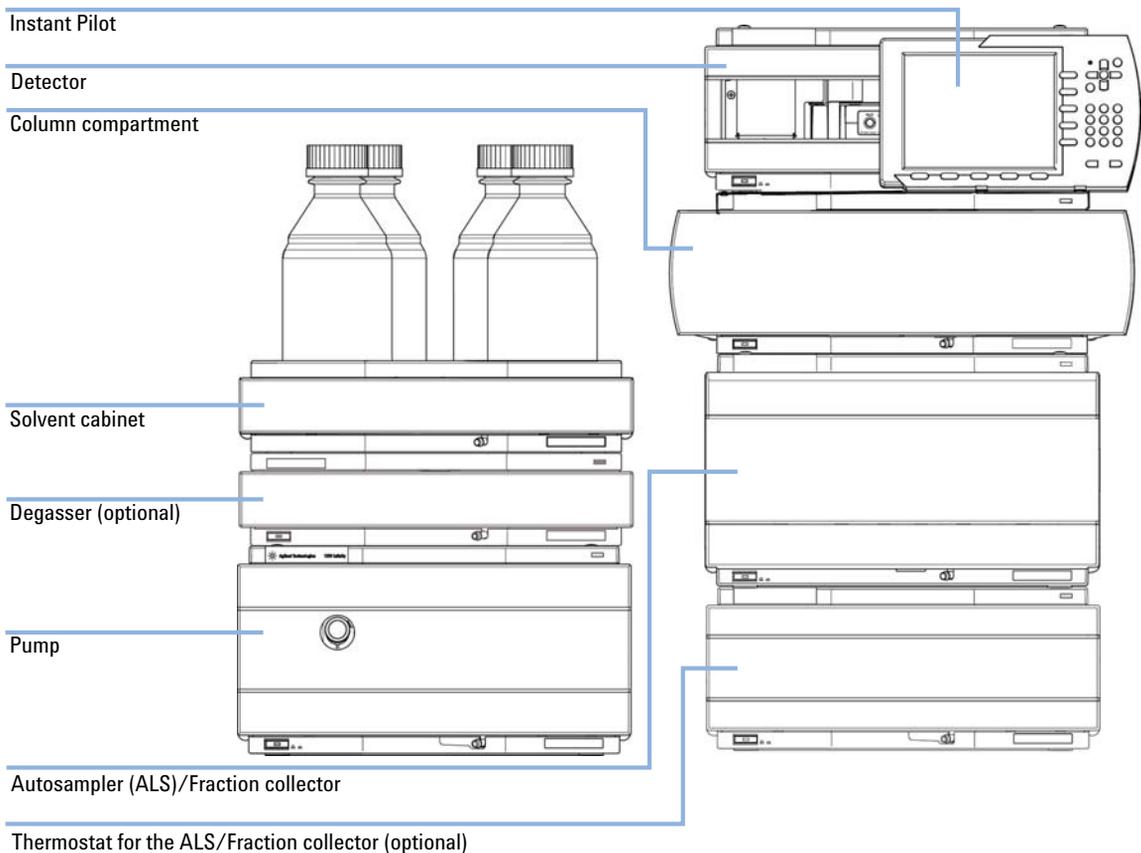

**Figure 9** Recommended Two Stack Configuration for 1260 Infinity (Front View)

### 3 Installing the Detector


#### Optimizing the Stack Configuration

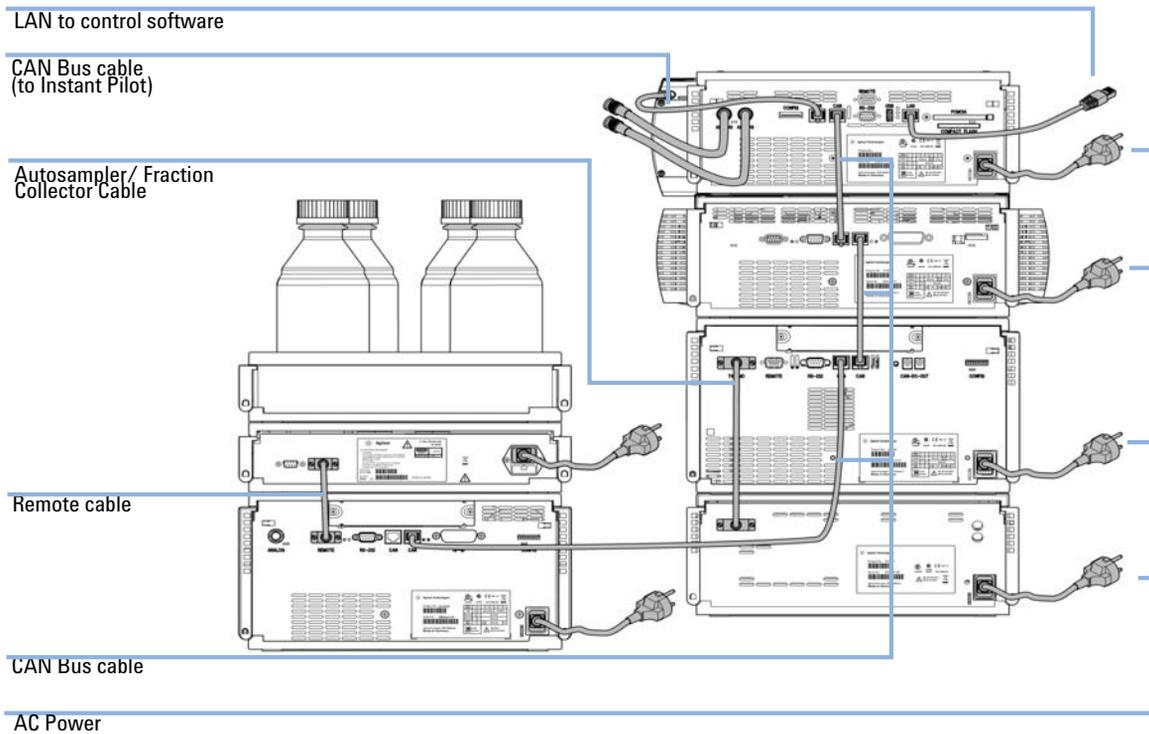

**Figure 10** Recommended Two Stack Configuration for 1260 Infinity (Rear View)

### Two Stack Configuration for Agilent 1290 Infinity LC

In case the autosampler thermostat is added to the system, a two-stack configuration is recommended, which places both heavy modules (1290 Infinity pump and thermostat) at the bottom of each stack and avoids high stacks. Some users prefer the lower height of this arrangement even without the autosampler thermostat. A slightly longer capillary is required between the pump and autosampler. (See [Figure 11](#) on page 43 and [Figure 12](#) on page 44).

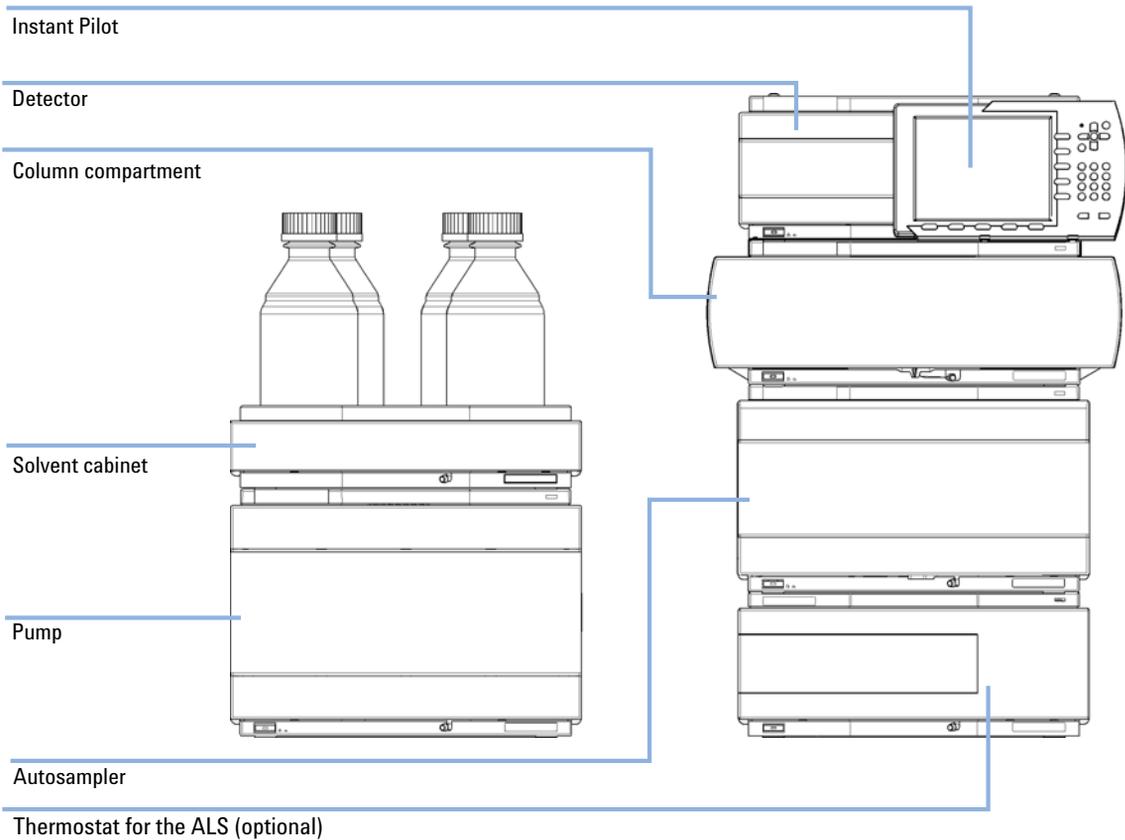

**Figure 11** Recommended two stack configuration for 1290 Infinity with binary pump (front view)



## Installation Information on Leak and Waste Handling

The Agilent 1200 Infinity Series has been designed for safe leak and waste handling. It is important that all security concepts are understood and instructions are carefully followed.

### WARNING

**Toxic, flammable and hazardous solvents, samples and reagents**

**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
- The volume of substances should be reduced to the minimum required for the analysis.
- Never exceed the maximal permissible volume of solvents (6 L) in the solvent cabinet.
- Do not use bottles that exceed the maximum permissible volume as specified in the usage guideline for the Agilent 1200 Infinity Series Solvent Cabinets.
- Arrange the bottles as specified in the usage guideline for the solvent cabinet.
- A printed copy of the guideline has been shipped with the solvent cabinet, electronic copies are available on the Internet.


### NOTE

**Recommendations for Solvent Cabinet**

For details, see the usage guideline for the Agilent 1200 Infinity Series Solvent Cabinets.

### 3 Installing the Detector

#### Installation Information on Leak and Waste Handling

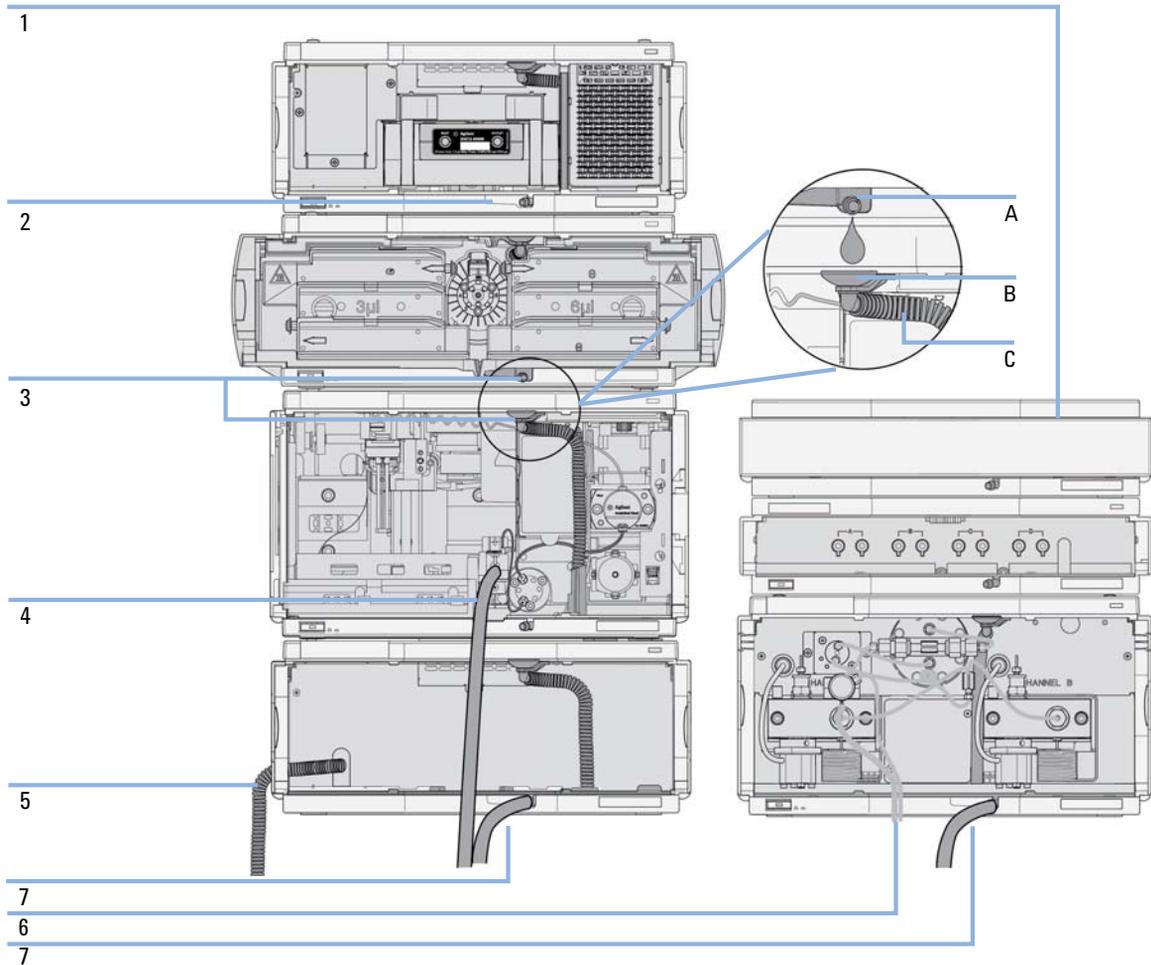

**Figure 13** Leak and waste handling (overview - typical stack configuration as an example)

1	Solvent cabinet
2	Leak pan
3	Leak pan's outlet port (A), leak funnel (B) and corrugated waste tube (C)
4	Waste tube of the sampler's needle wash
5	Condense drain outlet of the autosampler cooler
6	Waste tube of the purge valve
7	Waste tube

- 1 Stack the modules according to the adequate stack configuration.  
The leak pan outlet of the upper module must be vertically positioned above the leak tray of the lower module, see [Figure 13](#) on page 46.
- 2 Connect data and power cables to the modules, see section *Installing the Module* below.
- 3 Connect capillaries and tubes to the modules, see section *Flow Connections to the module* below or the relevant system manual.

**WARNING****Toxic, flammable and hazardous solvents, samples and reagents**

- Keep solvent path free from blockages.
- Keep the flow path closed (in case the pump in the system is equipped with a passive inlet valve, solvent may leak out due to hydrostatic pressure, even if your instrument is off).
- Avoid loops.
- Tubes must not sag.
- Do not bend tubes.
- Do not immerse tube end in waste liquid.
- Do not intubate tubes in other tubes.
- For correct tubing follow instructions on label attached to the module.

### 3 Installing the Detector

#### Installation Information on Leak and Waste Handling



**Figure 14** Warning label (illustration for correct waste tubing)

## Installing the Detector

Parts required	#	p/n	Description
	1		Detector
	1		Power cord
	1		LAN cable (cross-over or twisted pair network cable)
	1		Agilent ChemStation or other control software
	1	G4208A	Instant Pilot
	1		CompactFlash Card (G1314E only)

For other cables see below and section “[Cable Overview](#)” on page 164.

Instant Pilot (G4208A) is optional.

Preparations	
	Other LC modules must have appropriate firmware installed to work with the detector.
	Locate bench space
	Provide power connections
	Unpack the detector

### NOTE

Before adding the detector into an existing system assure that the existing modules have been updated to firmware revision that is supported by the control software.


### NOTE

If you want to use the "data-never-lost" feature, then assure that the CompactFlash Card is installed in the rear of the G1314E VWD.

### 3 Installing the Detector

#### Installing the Detector

- 1 Note the MAC address of the LAN interface (rear of the module, under the configuration switch, see figure below). It's required for LAN Configuration, see chapter *LAN Configuration*.

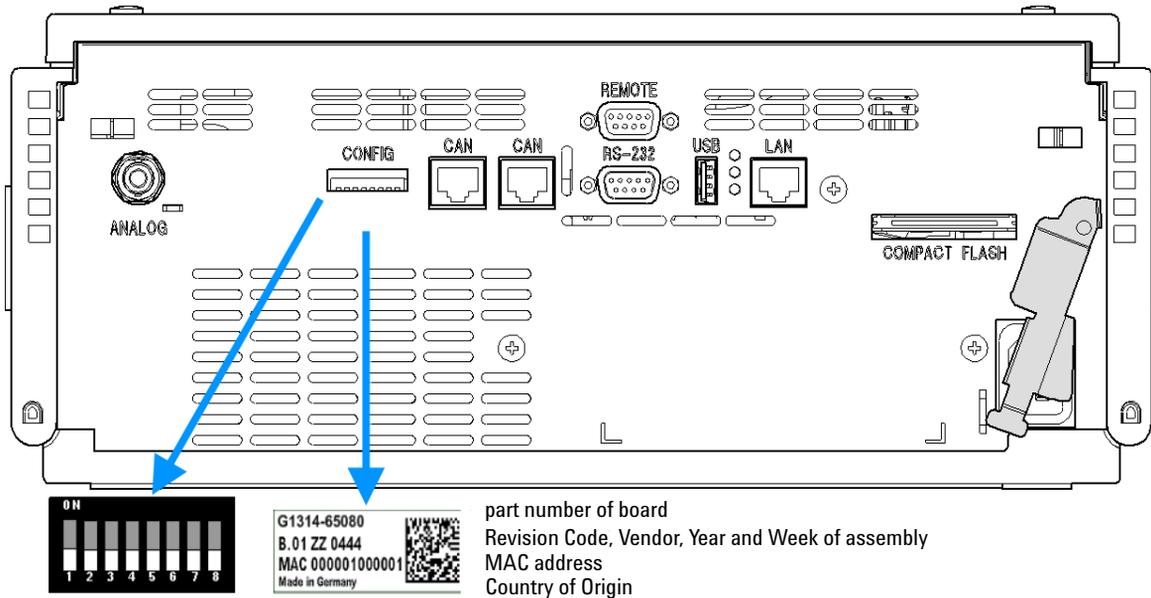

**Figure 15** Rear View of Detector

- 2 Check the setting of the DIP switch at the rear of the detector. All switches should be in down position (BOOTP mode). If another boot mode is required refer to “[Link configuration selection](#)” on page 208.

#### NOTE

The detector was shipped with default configuration settings (all switches down).

- 3 Place the detector in the stack or on the bench in a horizontal position.
- 4 Ensure the line power switch at the front of the detector is OFF.
- 5 Connect the power cable to the power connector at the rear of the detector.
- 6 Connect the CAN cable to other modules.


- 7 Connect the LAN cable (e.g. from a Agilent ChemStation as controller) to the detector's LAN connector.

**NOTE**

In multi-detector configurations the LAN of the Agilent detector with the highest data rate must be used due to its higher data load.

- 8 Connect the analog cable(s) (optional).
- 9 Connect the APG remote cable (optional) for non-Agilent 1200 Infinity Series instruments.
- 10 Turn on power by pushing the button at the lower left hand side of the detector. The status LED should be green.

Status indicator  
green/yellow/red

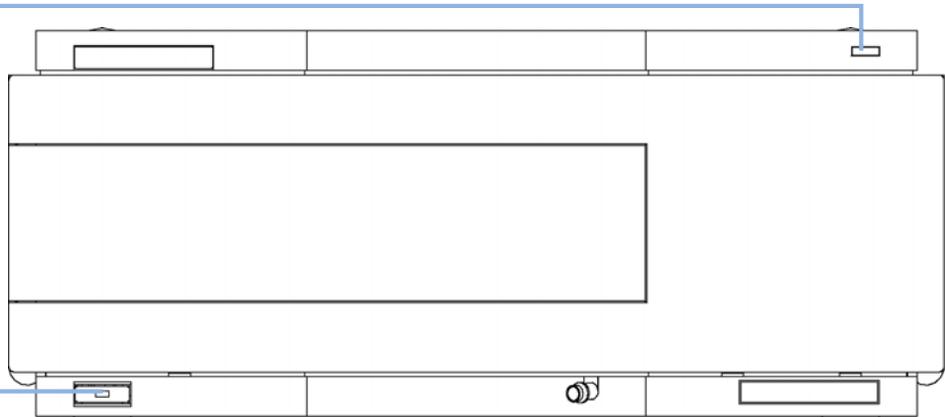

Line power switch  
with green light

**NOTE**

The detector is turned on when the line power switch is pressed and the green indicator lamp is illuminated. The detector is turned off when the line power switch is protruding and the green light is off.

**NOTE**

To disconnect the detector from line, unplug the power cord. The power supply still uses some power, even if the power switch at the front panel is turned off.

## Flow Connections to the Detector

<b>Tools required</b>	<b>Description</b>
	Wrench, 1/4 – 5/16 inch (for capillary connections)
<b>Parts required</b>	<b>#</b> <b>p/n</b> <b>Description</b>
	1    G1314-68755    Accessory kit
<b>Hardware required</b>	Other modules depend on system setup
<b>Preparations</b>	Detector is installed in the LC system.

### WARNING

#### **Toxic, flammable and hazardous solvents, samples and reagents**


**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
- The volume of substances should be reduced to the minimum required for the analysis.
- Do not operate the instrument in an explosive atmosphere.

### NOTE


The flow cell is shipped with a filling of isopropanol (also recommended when the instrument and/or flow cell is shipped to another location). This is to avoid breakage due to subambient conditions.

- 1** Press the release buttons and remove the front cover to have access to the front area.

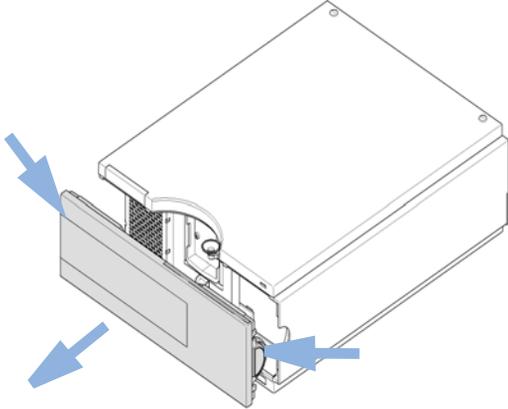

- 2** Loosen the screws of the flow cell dummy plate by turning each screw one turn. Then unscrew the screws completely. This is required to avoid any problem with helicoil insert in the casting.

- 3** Press the flow cell completely into the slot and tighten the cell screws (both parallel) until the mechanical stop.

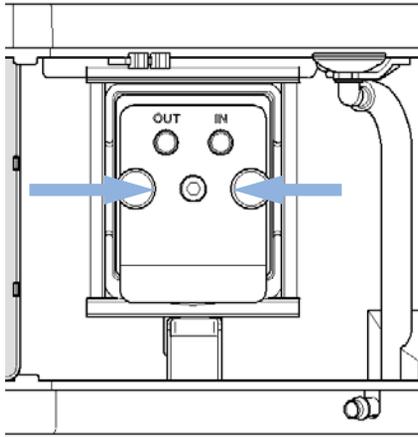


- 4** Assemble the column-detector capillary. Depending on the flow cell type it may be a PEEK or SST capillary.


### 3 Installing the Detector

#### Flow Connections to the Detector


- 5 Connect the newly assembled fitting of the capillary to the inlet connector and connect the other end of the capillary to the column.

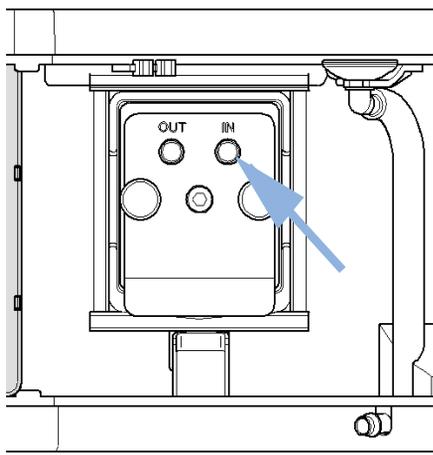

- 6 Connect the PEEK waste capillary to the outlet connector.

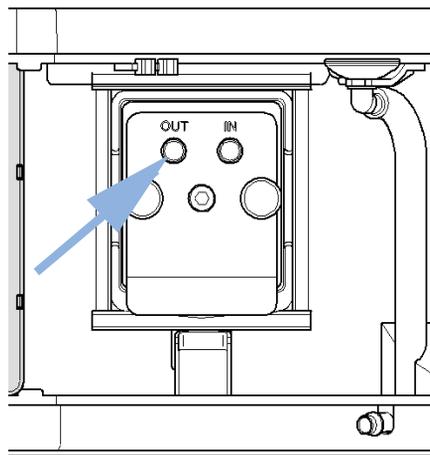

- 7 Establish a flow and observe for leakage.

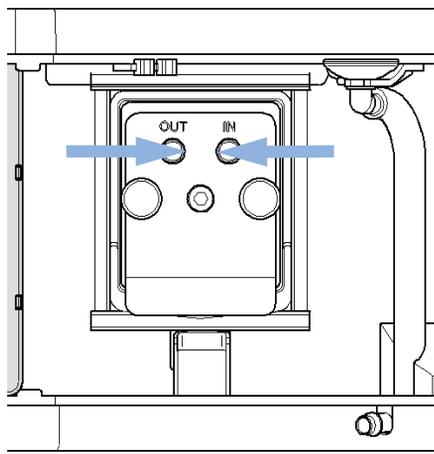


- 8 Replace the front cover.

The installation of the detector is now complete.

#### NOTE

The detector should be operated with the front cover in place to protect the flow cell area against strong drafts from the outside.

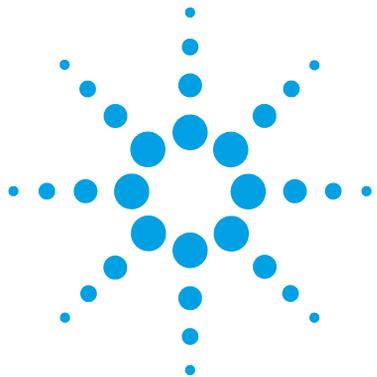

## 4 Using the Detector

Leak and Waste Handling	56
Setting up an Analysis	57
Before Using the System	57
Requirements and Conditions	59
Optimization of the System	61
Preparing the HPLC System	61
Running the Sample and Verifying the Results	70
Special Settings of the Detector	71
Control Settings	71
Configuration Settings	72
Online Spectra	73
Scanning with the VWD	74
Analog Output Settings	75
Special Setpoints	76
Run Recovery (G1314E)	80

This chapter provides information on how to set up the detector for an analysis and explains the basic settings.



## Leak and Waste Handling

### WARNING

**Toxic, flammable and hazardous solvents, samples and reagents**

**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
- The volume of substances should be reduced to the minimum required for the analysis.
- Do not operate the instrument in an explosive atmosphere.
- Never exceed the maximal permissible volume of solvents (6 L) in the solvent cabinet.
- Do not use bottles that exceed the maximum permissible volume as specified in the usage guideline for the Agilent 1200 Infinity Series Solvent Cabinets.
- Arrange the bottles as specified in the usage guideline for the solvent cabinet.
- A printed copy of the guideline has been shipped with the solvent cabinet, electronic copies are available on the Internet.
- The residual free volume in the appropriate waste container must be large enough to collect the waste liquid.
- Check the filling level of the waste container regularly.
- To achieve maximal safety, check the correct installation regularly.

### NOTE

**Recommendations for Solvent Cabinet**

For details, see the usage guideline for the Agilent 1200 Infinity Series Solvent Cabinets.

For details on correct installation, see [“Installation Information on Leak and Waste Handling”](#) on page 45.

## Setting up an Analysis

This chapter can be used for

- preparing the system,
- to learn the set up of an HPLC analysis and
- to use it as an instrument check to demonstrate that all modules of the system are correctly installed and connected. It is not a test of the instrument performance.
- Learn about special settings

## Before Using the System

### Solvent Information

Observe recommendations on the use of solvents in chapter “Solvents” in the pump’s reference manual.

### Priming and Purging the System

When the solvents have been exchanged or the pumping system has been turned off for a certain time (for example, overnight) oxygen will re-diffuse into the solvent channel between the solvent reservoir, vacuum degasser (when available in the system) and the pump. Solvents containing volatile ingredients will slightly lose these. Therefore priming of the pumping system is required before starting an application.

## 4 Using the Detector

### Setting up an Analysis

**Table 7** Choice of Priming Solvents for Different Purposes

Activity	Solvent	Comments
After an installation	Isopropanol	Best solvent to flush air out of the system
When switching between reverse phase and normal phase (both times)	Isopropanol	Best solvent to flush air out of the system
After an installation	Ethanol or Methanol	Alternative to Isopropanol (second choice) if no Isopropanol is available
To clean the system when using buffers	Bidistilled water	Best solvent to re-dissolve buffer crystals
After a solvent change	Bidistilled water	Best solvent to re-dissolve buffer crystals
After the installation of normal phase seals (P/N 0905-1420)	Hexane + 5% Isopropanol	Good wetting properties

#### NOTE

The pump should never be used for priming empty tubings (never let the pump run dry). Use a syringe to draw enough solvent for completely filling the tubings to the pump inlet before continuing to prime with the pump.

- 1 Open the purge valve of your pump (by turning it counterclockwise) and set flow rate to 3 – 5 mL/min.
- 2 Flush all tubes with at least 30 mL of solvent.
- 3 Set flow to required value of your application and close the purge valve.

#### NOTE

Pump for approximately 10 minutes before starting your application.

## Requirements and Conditions

### What You Will Need

The table below lists the items you need to have for the set up of the analysis. Some of these are optional (not required for the basic system).

**Table 8** What you will need

Agilent 1200 Infinity Series system	Pump (plus degassing)
	Autosampler
	Detector, standard flow cell installed
	Degasser (optional)
	Column Compartment (optional)
	Agilent ChemStation or Instant Pilot G4208, optional for basic operation.
	System should be correctly set up for LAN communication with the Agilent ChemStation
Column:	Zorbax Eclipse XDB-C18, 4.6 x 150 mm, 5 µm (993967-902) or an equivalent column
Standard:	Agilent isocratic checkout sample (01080-68704)

### Conditions


A single injection of the isocratic test standard is made under the conditions given in [Table 9](#) on page 59:

**Table 9** Conditions

Flow	1.5 mL/min
Stoptime	8 min
Solvent	100% (30% water/70% Acetonitrile)
Temperature	Ambient
Wavelength	sample 254 nm
Injection Volume	1 µL
Column Temperature (optional):	25 °C or ambient

### Typical Chromatogram

A typical chromatogram for this analysis is shown in [Figure 16](#) on page 60. The exact profile of the chromatogram will depend on the chromatographic conditions. Variations in solvent quality, column packing, standard concentration and column temperature will all have a potential effect on peak retention and response.

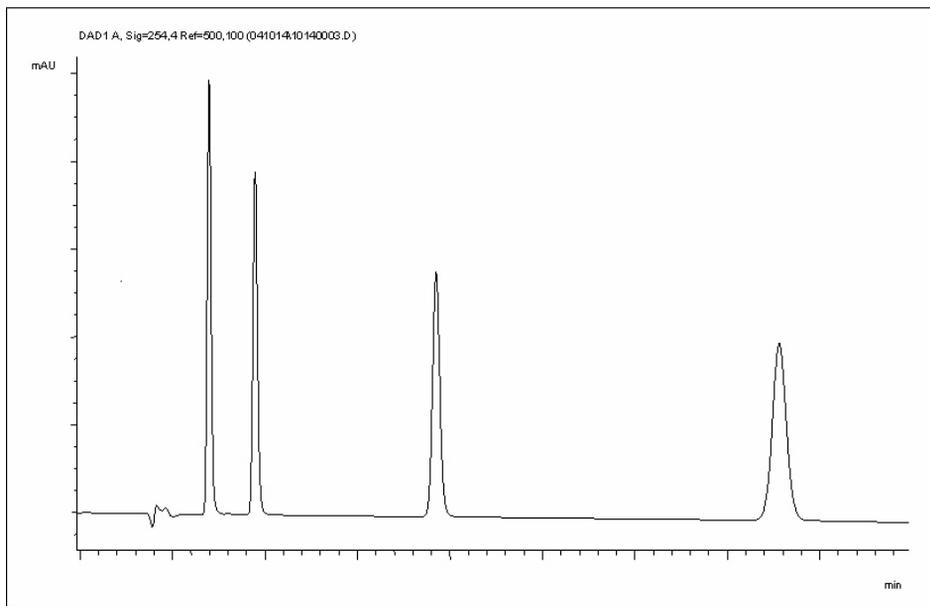


**Figure 16** Typical Chromatogram with UV-detector

## Optimization of the System

The settings used for this analysis are specific for this purpose. For other applications the system can be optimized in various ways. Please refer to the section “[Optimizing the Detector Performance](#)” on page 84.

## Preparing the HPLC System

- 1 Turn on the Agilent ChemStation PC and the monitor.
- 2 Turn on the modules.
- 3 Start the Agilent ChemStation software. If the pump, autosampler, thermostatted column compartment and detector are found, the Agilent ChemStation screen should look like shown in [Figure 17](#) on page 61. The System status is red (**Not Ready**).

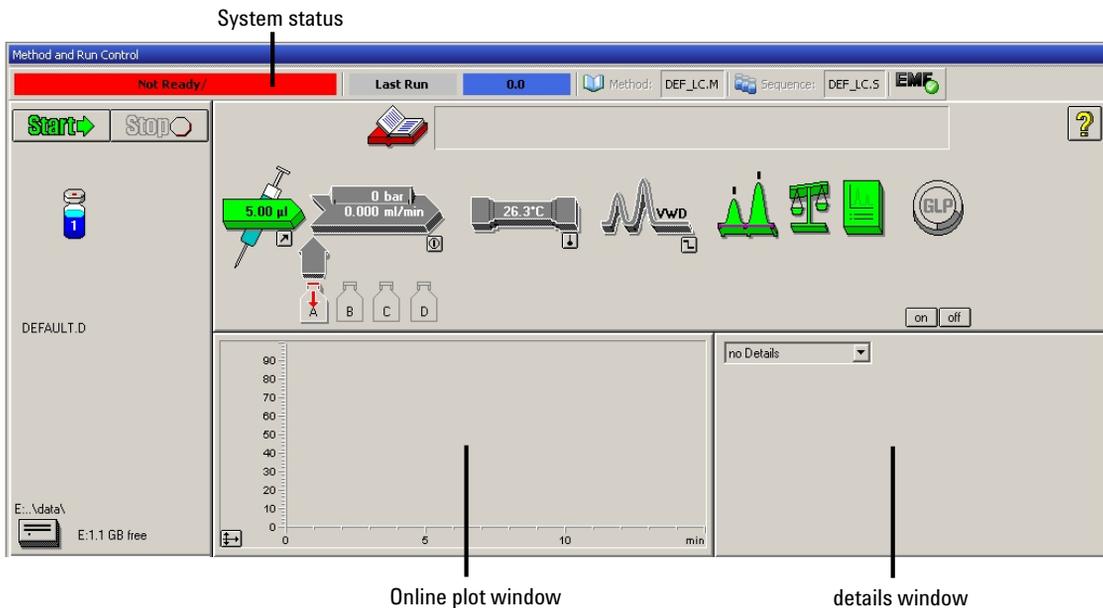

**Figure 17** Initial Agilent ChemStation screen (**Method and Run Control**)

## 4 Using the Detector Setting up an Analysis

- 4 Turn on the detector lamp, pump and autosampler by clicking the **System On** button or the buttons below the module icons on the graphical user interface (GUI).

After some time, the pump, thermostatted column compartment and detector module will turn to green.


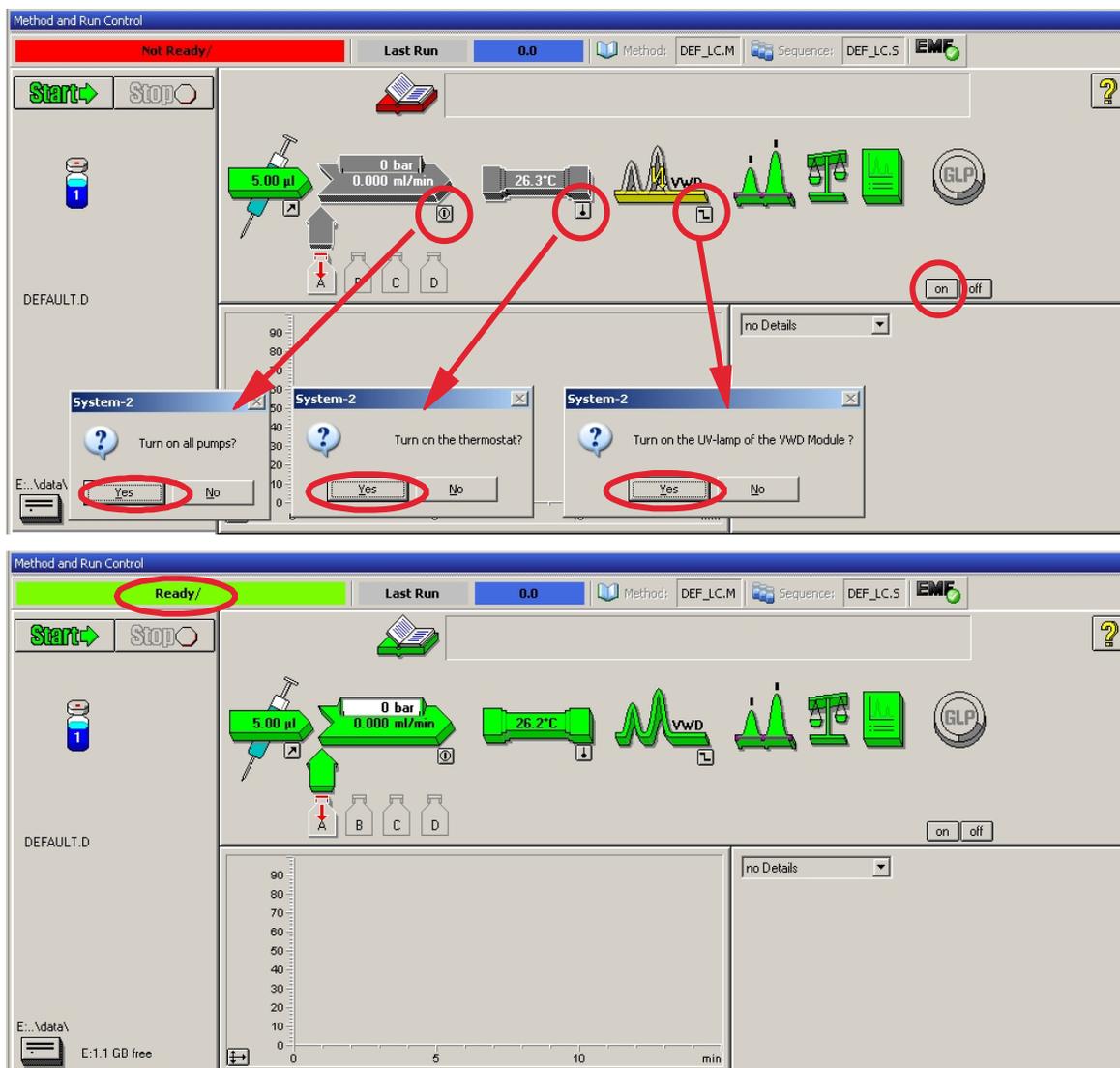


Figure 18 Turning on the HPLC Module

- 5 Purge the pump. For more information “[Priming and Purging the System](#)” on page 57.
- 6 Allow the detector to warm up of at least 60 minutes to provide a stable baseline (example: [Figure 19](#) on page 63).

**NOTE**

For reproducible chromatography, the detector and lamp should be on for at least one hour. Otherwise the detector baseline may still drift (depending on the environment).

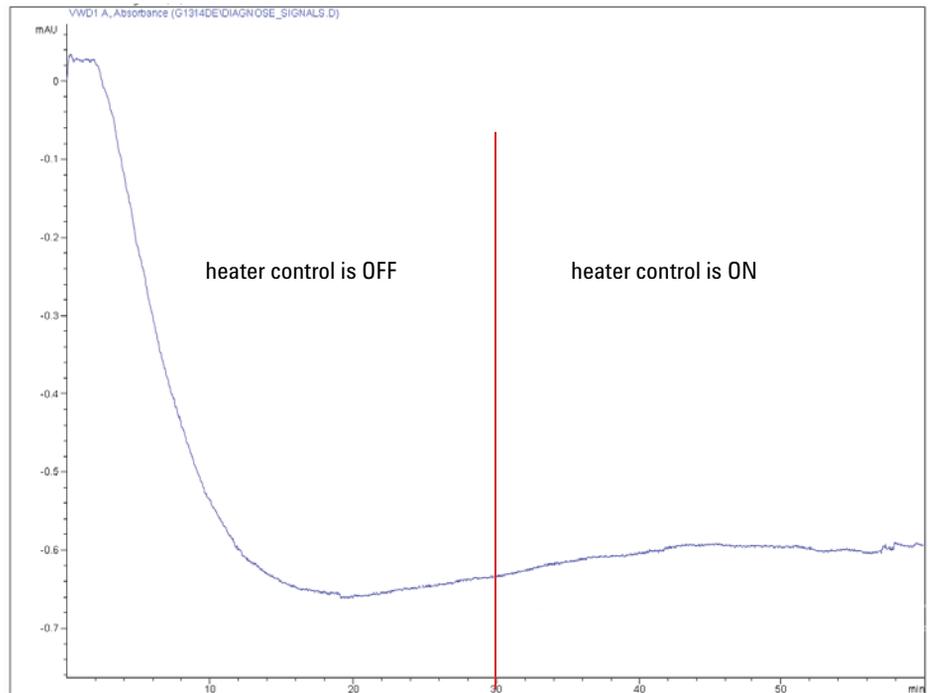


**Figure 19** Stabilization of baseline after turn on of detector

- 7 For the isocratic pump, fill the solvent bottle with the mixture of HPLC-grade bi-distilled water (30 %) and acetonitrile (70 %). For binary- and quaternary pumps you can use separate bottles.

## 4 Using the Detector Setting up an Analysis

- Click on the **Load Method** button, select **DEF\_LC.M** and press **OK**. Alternatively, double-click on the method in the method window. The default LC method parameters are transferred into the modules.

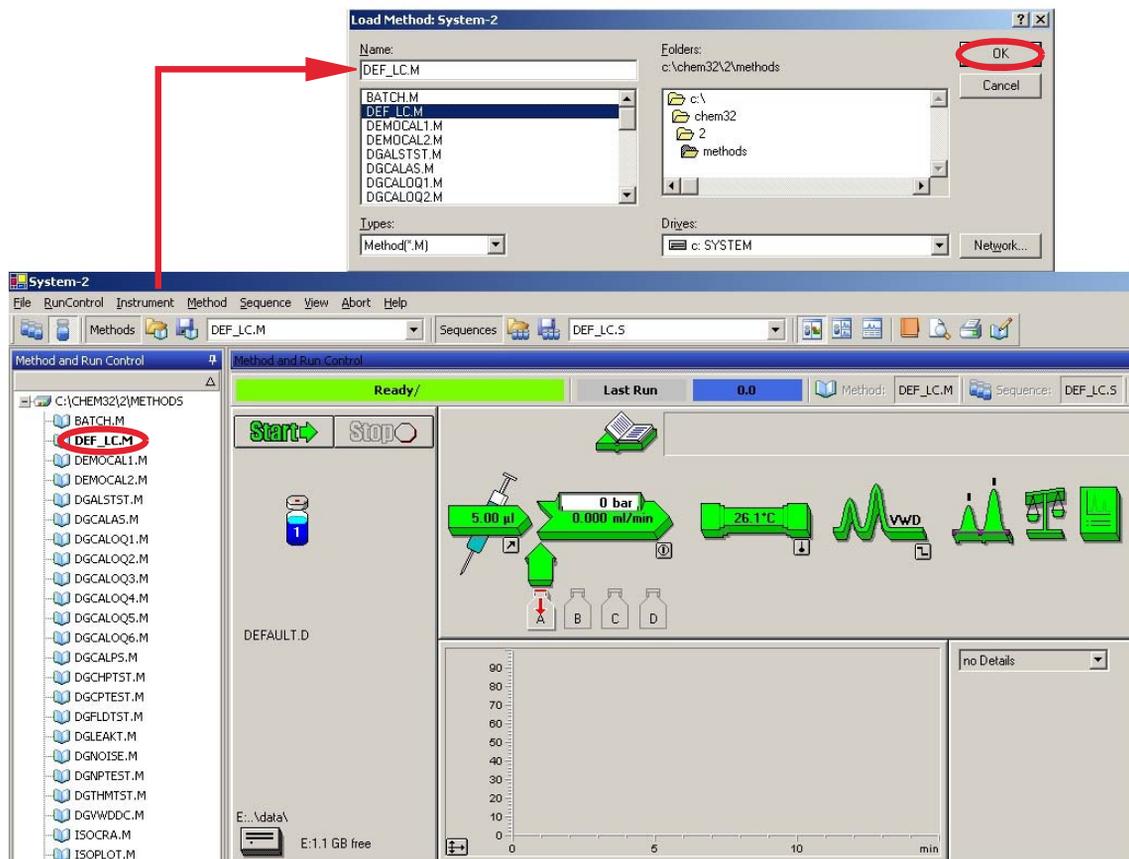

**Figure 20** Loading Default LC Method

- Click on the module icons (Figure 21 on page 65) and open the **Setup** of these modules. Figure 22 on page 66 shows the detector settings (do not change the detector parameters at this time).


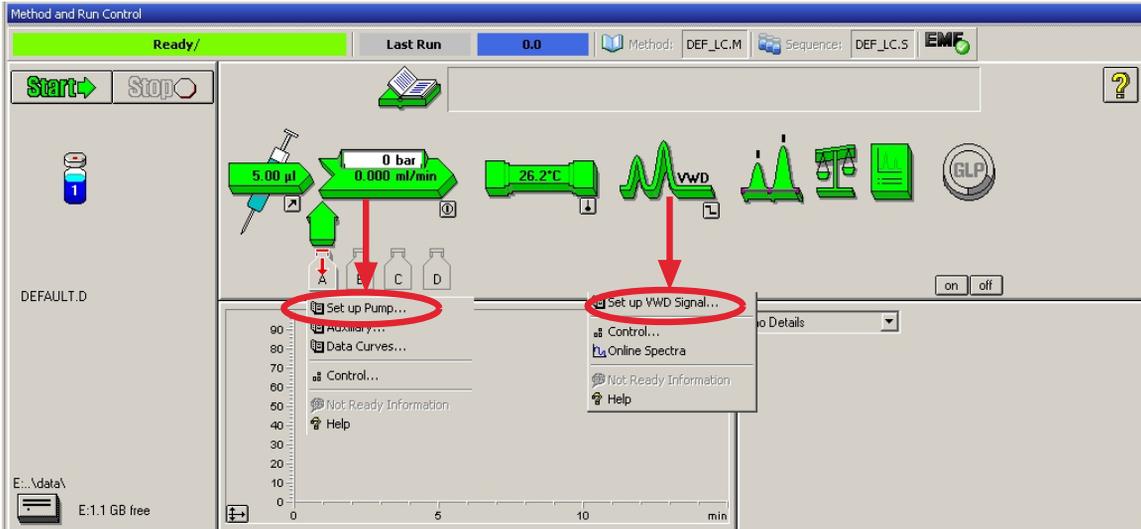


Figure 21 Open the module menu

## 4 Using the Detector

### Setting up an Analysis

10 Enter the pump parameters mentioned under [Table 9](#) on page 59.

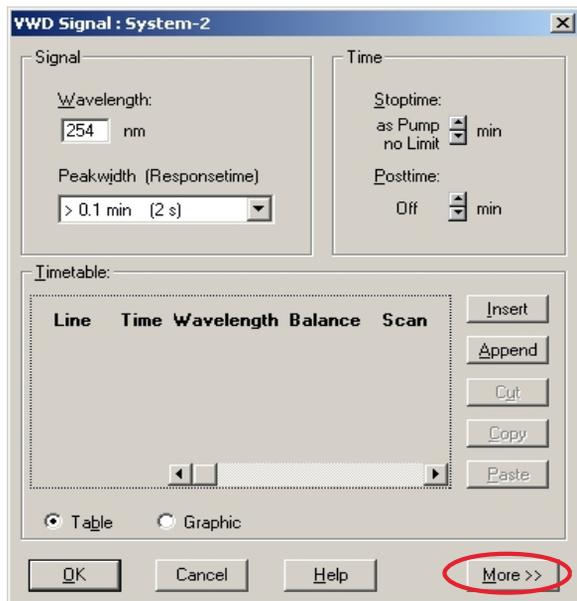

- 1 signal with individual wavelength setting
- stop and post time can be set (if required)
- time table for programmable actions during the run

- Zero Offset Limits: 1 to 99 % in steps of 1 %
- Attenuation Limits: 0.98 to 4000 mAU at discrete values for either 100 mV or 1 V full scale

- additional signals can be stored with the normal signal (for diagnostics)

- autobalance to zero absorbance (on the analog output plus offset) at begin and/or end of run

- see [“Special Setpoints”](#) on page 76.

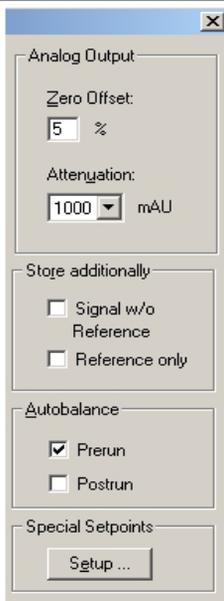


Figure 22 Detector Settings (default)

- 11 Pump the water/acetonitrile (30/70 %) mobile phase through the column for 10 minutes for equilibration.
- 12 Click the button  and select **Change...** to open the Signal Plot information. Select the **Pump: Pressure** and the **VWD A: Signal 254** as signals. Change the Y-range for the VWD to 1 mAU and the offset to 20 % and the pressure offset to 50 %. The X-axis range should be 15 minutes. Press **OK** to exit this screen.


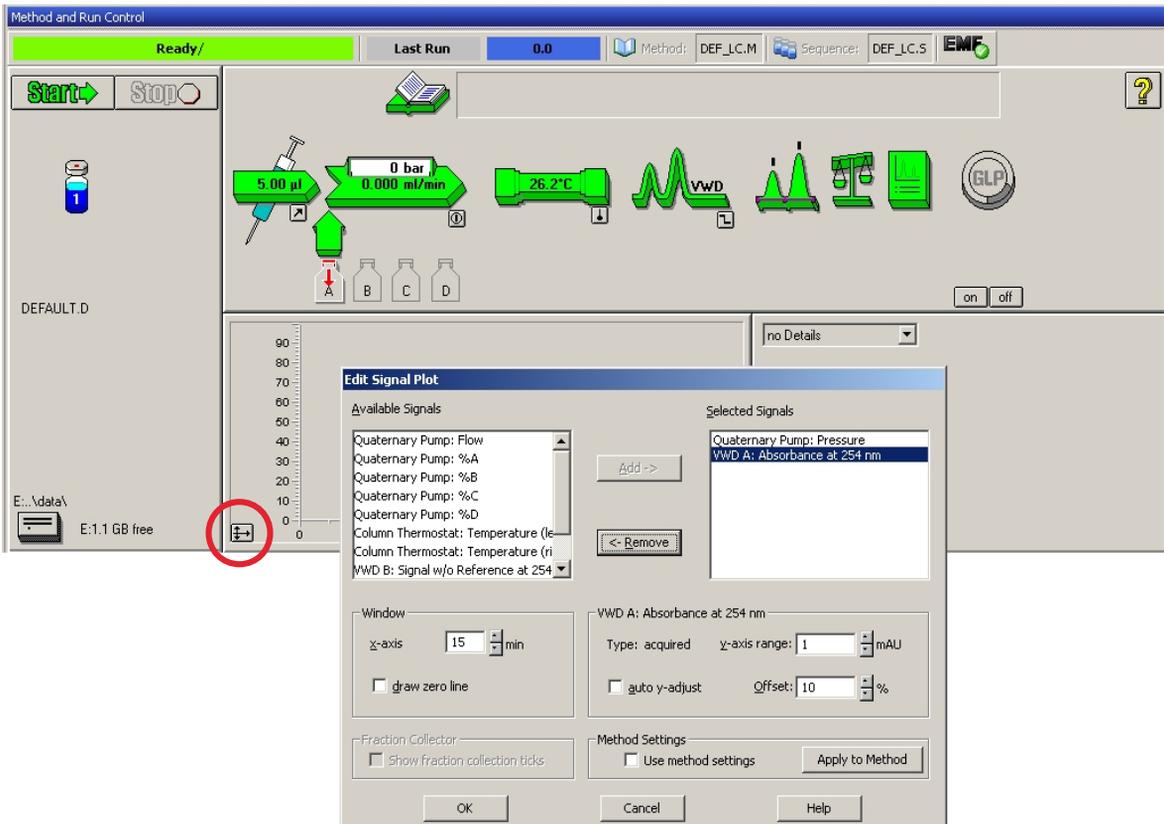


Figure 23 Edit Signal Plot Window

## 4 Using the Detector

### Setting up an Analysis

The Online Plot (Figure 24 on page 68) shows both, the pump pressure and the detector absorbance signals. Pressing **Adjust** the signals can be reset to the offset value and **Balance** would do a balance on the detector.

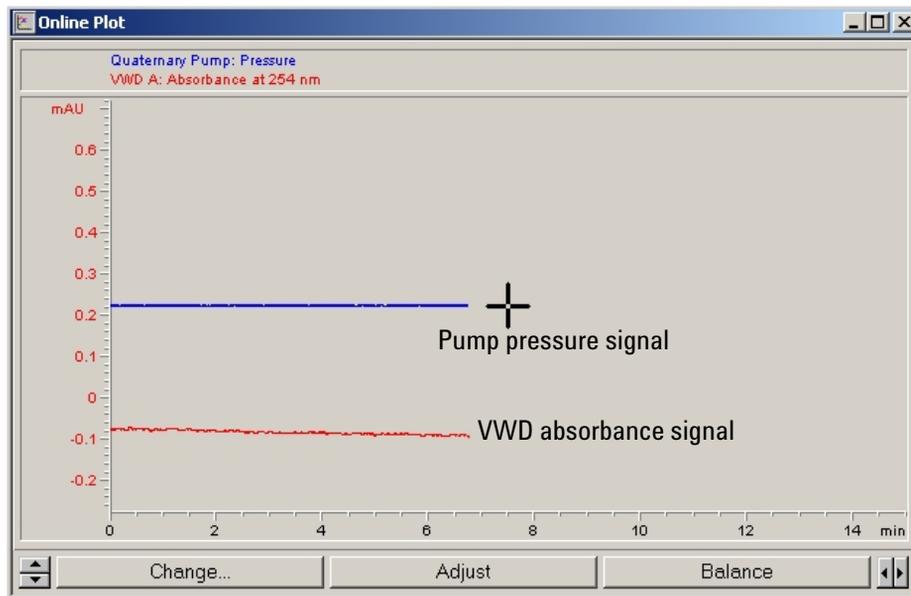

**Figure 24** Online Plot Window

**13** If both baselines are stable, set the Y-range for the detector signal to 100 mAU.

#### NOTE

If you start with a new UV-lamp for the first time, the lamp may show initial drift for some time (burn-in effect).

- 14 Select the menu item **RunControl** > **Sample Info** and enter information about this application (Figure 25 on page 69). Press **OK** to leave this screen.


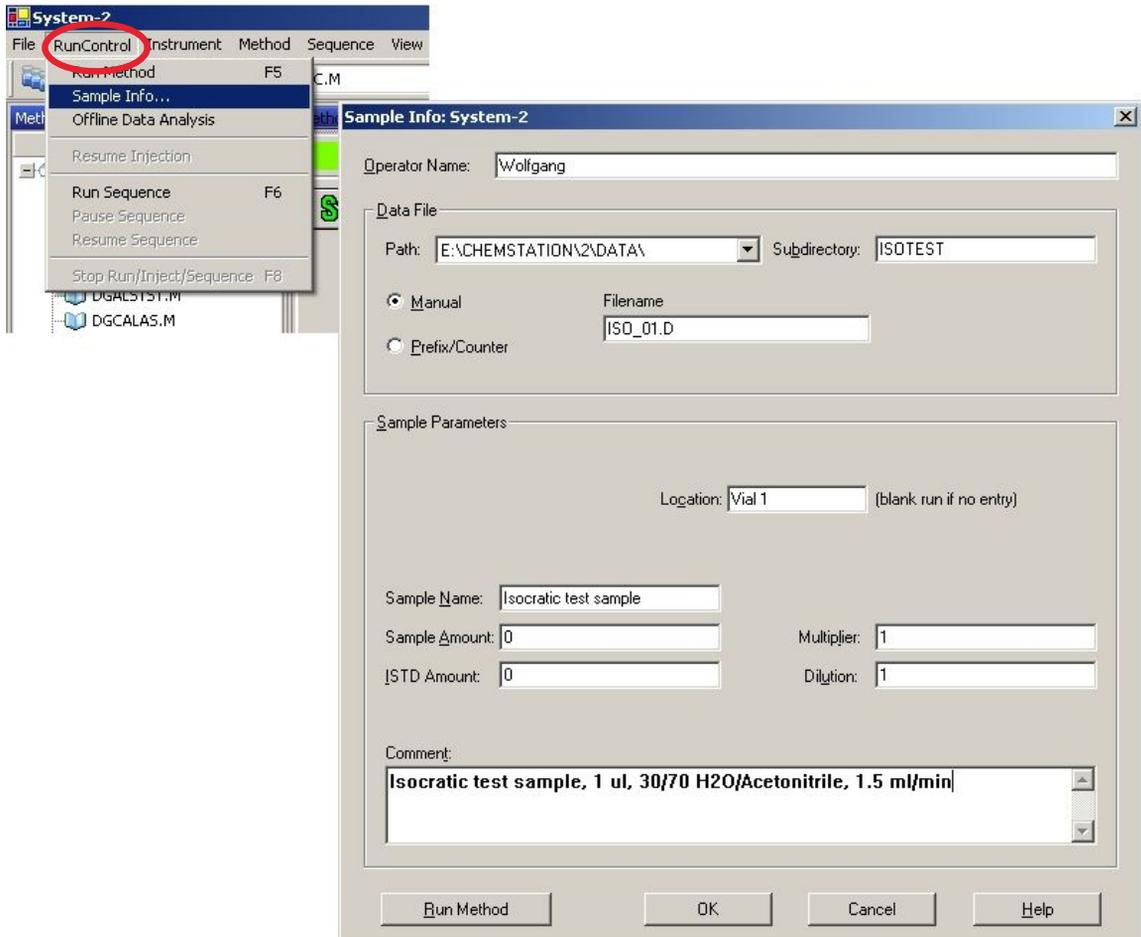


Figure 25 Sample Information

- 15 Fill the content of an isocratic standard sample ampoule into a vial and seal the vial with a cap and place the vial into autosampler tray (position #1).

## Running the Sample and Verifying the Results

- 1 To start a run select the menu item **RunControl > Run Method**.
- 2 This will start the modules and the online plot on the Agilent ChemStation will show the resulting chromatogram.

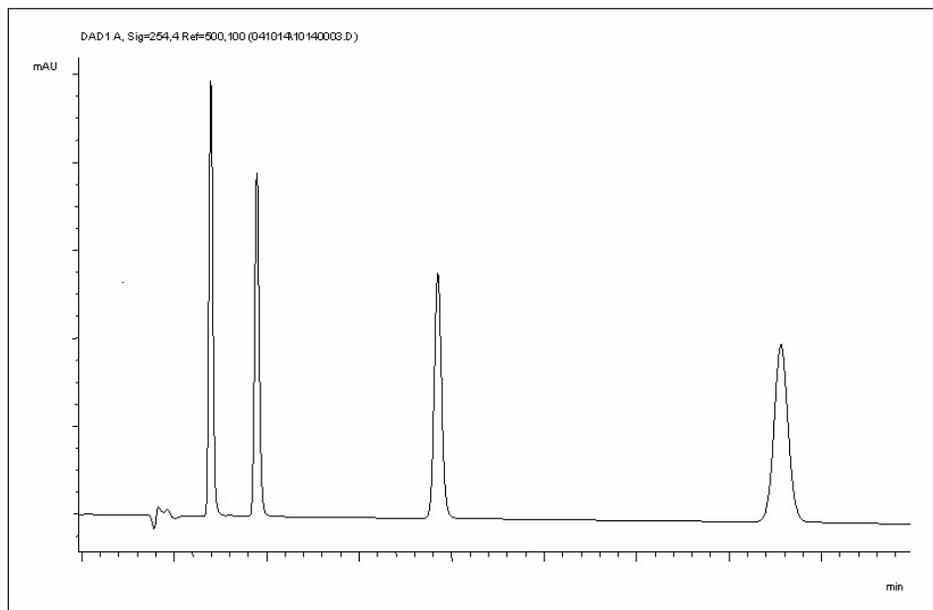

**Figure 26** Chromatogram with Isocratic Test Sample

**NOTE**

Information about using the Data Analysis functions can be obtained from the Using your ChemStation manual supplied with your system.

## Special Settings of the Detector

In this chapter special settings of the detector are described.

### Control Settings

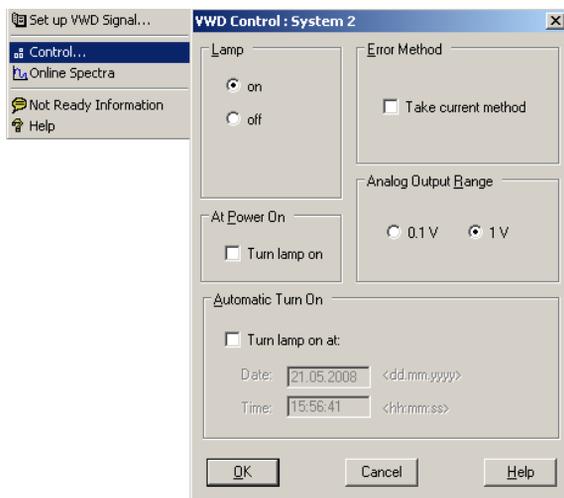




Figure 27 Detector control settings

- **Lamp:** turn on and off of UV-lamp.
- **At Power On:** automatic lamp-on at power on.
- **Error Method:** take error method or current method (in case of an error).
- **Analog Output Range:** can be set to either 100 mV or 1 V full scale, see “Analog Output Settings” on page 75.
- **Automatic Turn On:** lamps can be programmed (detector must be on for this).
- **Help:** online help.

## 4 Using the Detector

### Special Settings of the Detector

# Configuration Settings

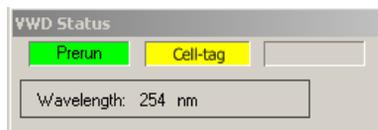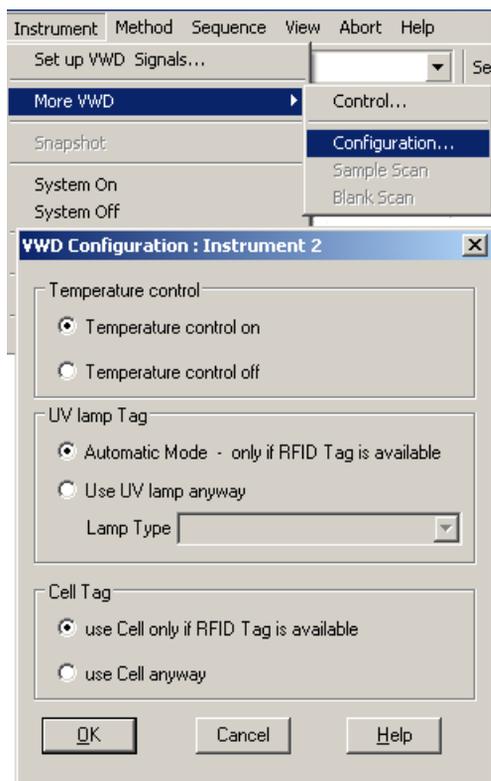

**Figure 28** Detector Configuration Settings

- **Temperature Control:** the optical unit is kept on constant temperature (some degrees above ambient) and improves the baseline stability in unstable environments. See also note below.
- **UV lamp tag:** Automatic mode for Agilent lamps with RFID tags. If no RFID tag lamp is used, detector icon will become gray (lamp tag not ready) and analysis is disabled. The specifications are based on the lamp with RFID tag. Use UV lamp anyway - here you can select non-RFID tag lamps, such as of type VWD lor DAD (have different heating mechanismn). The correct selection is important for optimal performance and lifetime.
- **Cell tag:** for Agilent flow cells with RFID tags. If no RFID tag cell is used, detector icon will become gray (cell tag not ready) and analysis is disabled.
- **Help:** online help.

The detector status shows “Cell-tag” in yellow in case the flow cell with the RFID tag is not inserted. The detector icon is gray and the system is not ready.

## NOTE


If the flow cell temperature is critical for your chromatography or your environment is stable, you may set the Temperature Control to off. This will lower the optical unit and flow cell temperature by some degree.

## Online Spectra

- 1 To view the online spectra select **Online Spectra**.

### NOTE

This online spectrum is taken during a stop-flow condition only while the peak is kept in the flow cell, see “Scanning with the VWD” on page 74.

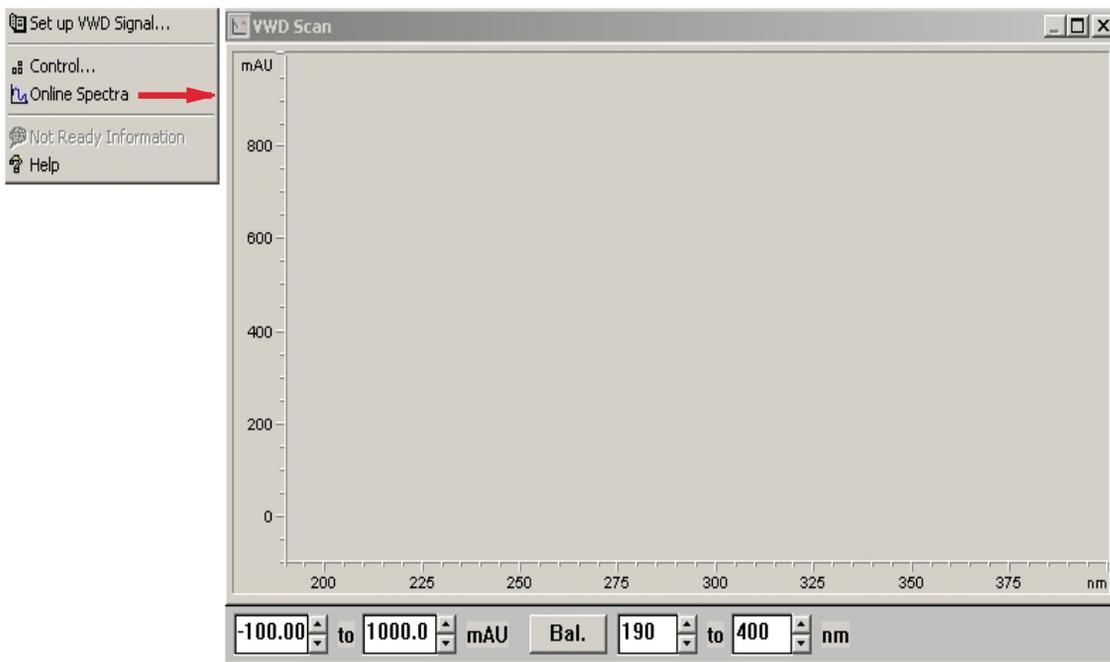

**Figure 29** Online Spectra Window

- 2 Change the absorbance and wavelength range according your needs.


## Scanning with the VWD

### NOTE

Access to the scan feature is only possible during run.

- 1 Set up a run.
- 2 Start a run.
- 3 While running on the baseline, select from the menu **Instrument > More VWD > Blank Scan**.

A background scan is stored in the memory.

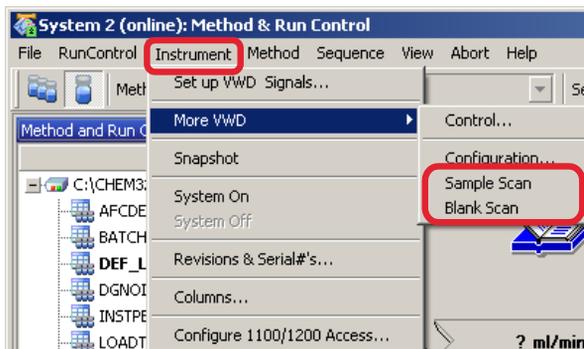

- *Step 1: Blank Scan:* scan of the background (solvent) is stored in the memory.
- *Step 2: Sample Scan:* scan of the peak of interest is taken while the peak stays in the flow cell (stop-flow condition).
- **Online Spectrum: Sample Scan minus Blank Scan.**

- 4 When the peak of interest enters the flow cell, stop the flow (set flow rate to zero or open the purge valve) and wait a few moments to stabilize the concentration.

### NOTE


Turning off the pump would stop the run and no access to the sample scan is possible.

- 5 Select from the menu **Instrument > More VWD > Sample Scan**.

A sample scan is taken in the range defined under “[Special Setpoints](#)” on page 76 and the Online Spectra window (see “[Online Spectra](#)” on page 73) displays the result (Sample Scan minus Blank Scan).

## Analog Output Settings

- 1 To change the Output Range of the analog outputs select **VWD Control**.
- 2 To change the offset and the attenuation select **VWD Signal > More**.

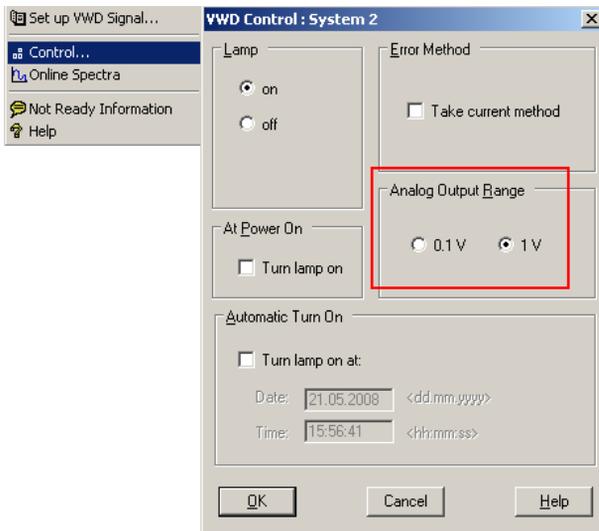

- **Analog Output Range:** can be set to either 100 mV or 1 V full scale.
- **Zero Offset:** can be set to either 100 mV or 1 V full scale.
- **Attenuation Limits:** 0.98 to 4000 mAU at discrete values for either 100 mV or 1 V full scale.

Figure 30 Analog Output Settings

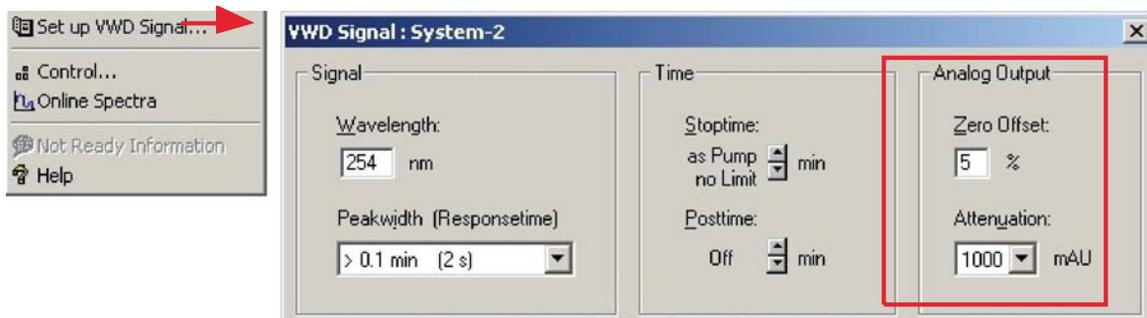


Figure 31 Analog Output Settings

- 3 Change the values if required.

## Special Setpoints

- 1 To change the offset and the attenuation select **VWD Signal > More > Special Setpoints**.

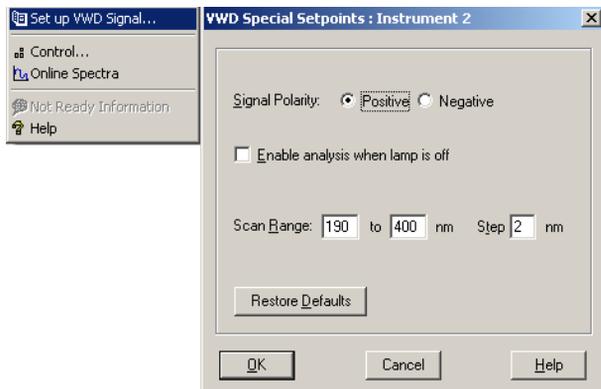


Figure 32 Special Setpoints

- **Signal Polarity:** can be switched to negative (if required).
- **Enable analysis when lamp is off:** if the VWD is not used in a dual detector setup (lamp off), the not-ready condition is not stopping the analysis.
- **Scan Range / Step:** Used for stop-flow scanning, see “Scanning with the VWD” on page 74.

## Peakwidth Settings

### NOTE

Do not use peak width shorter than necessary, see details below.

- 1 To change the Peakwidth settings select **Setup Detector Signals**.
- 2 In the section **Peakwidth (Responsetime)** click on the drop-down list.
- 3 Change the Peakwidth according to your needs.

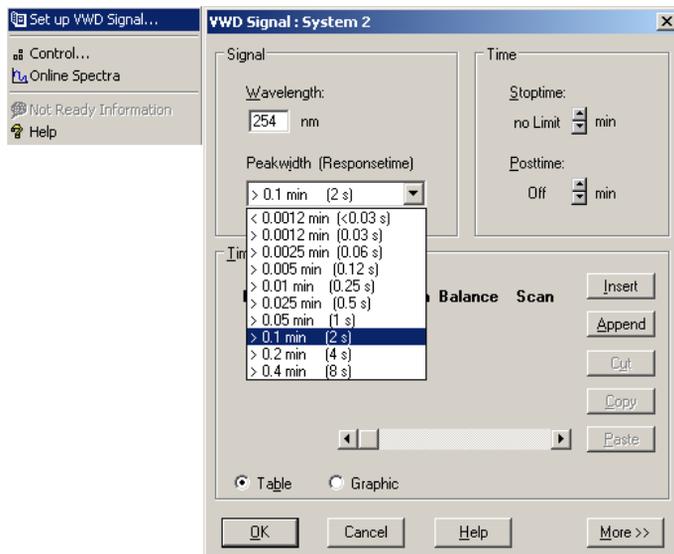


Figure 33 Peakwidth Setting

**Peakwidth** enables you to select the peak width (response time) for your analysis. The peak width is defined as the width of a peak, in minutes, at half the peak height. Set the peak width to the narrowest expected peak in your chromatogram. The peak width sets the optimum response time for your detector. The peak detector ignores any peaks that are considerably narrower, or wider, than the peak width setting. The response time is the time between 10 % and 90 % of the output signal in response to an input step function.

**Limits:** When you set the peak width (in minutes), the corresponding response time is set automatically and the appropriate data rate for signal acquisition is selected as shown in [Table 10](#) on page 78, [Table 11](#) on page 78 and [Table 12](#) on page 79 .

## 4 Using the Detector

### Special Settings of the Detector

**Table 10** Peak Width – Response Time – Data Rate (G1314D)

peakwidth at half heighth [min]	Response Time [s]	Data Rate [Hz]
<0.005	<0.12	20
>0.005	0.12	20
>0.01	0.25	20
>0.025	0.5	20
>0.05	1.0	10
>0.10	2.0	5
>0.20	4.0	2.5
>0.40	8.0	1.25

**Table 11** Peak Width - Response Time - Data Rate (G1314F)

peakwidth at half heighth [min]	Response Time [s]	Data Rate [Hz]
<0.003125	<0.0625	80
>0.003125	0.0625	80
>0.00625	0.125	80
>0.0125	0.25	40
>0.025	0.5	20
>0.05	1	10
>0.1	2	5
>0.2	4	2.5
>0.4	8	1.25

**Table 12** Peak Width – Response Time – Data Rate (G1314E)

<b>peakwidth at half heighth [min]</b>	<b>Response Time [s]</b>	<b>Data Rate [Hz]</b>
<0.0012	<0.03	160
>0.0012	0.03	160
>0.0025	0.06	160
>0.005	0.12	80
>0.01	0.25	40
>0.025	0.5	20
>0.05	1.0	10
>0.1	2.0	5
>0.2	4.0	2.5
>0.4	8.0	1.25

## Run Recovery (G1314E)

### NOTE

This feature is not supported with ChemStation B.03.02 SR1. It will be implemented with ChemStation B.04.01. The figures used in this chapter are from the G1315C DAD VL+. The G1314E VWD figures will look the same.

---

### Run Recovery Settings

### CAUTION

For this recovery mode the CompactFlash Card must be in the detector.

If the LAN communication is interrupted, no data is stored.

→ Have the CompactFlash Card always inserted.

---

The detector supports run buffering, which means that an amount of run data (\*.uv and \*.ch files) is stored in a storage medium (CompactFlash Card) in the detector until either it is overwritten or the detector undergoes a power cycle.

If there is a temporary network failure or the PC is not able to constantly take the data, the stored data is transferred to the ChemStation automatically when the network connection is restored or the PC can take the data, so that no loss of data occurs.

If there is a permanent network failure, the Run Recovery dialog box allows you to restore the stored data to the data directory. From there you can copy the files to the directory where the files are corrupted or not complete.

### NOTE

On very large recovery files it may take a long time to restore it to the Agilent ChemStation. A sequence will be stopped in case of a network problem.

---

### NOTE




When during recovery an error "Method/Sequence stopped" appears, the instrument logbook shows an entry "No Run data available in device".

In this case refer to "[No Run Data Available In Device](#)" on page 112.

---

## Automated Run Recovery in case of temporary communication failures

**Table 13** Automated Run Recovery in case of temporary communication failures

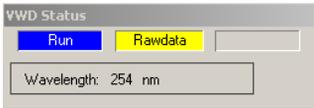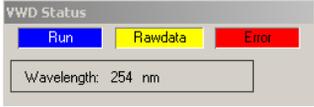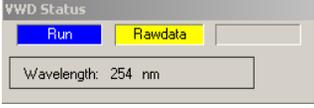Situation	Reaction	On ChemStation
All OK	<ul style="list-style-type: none"> <li>Run in progress - Data Analysis</li> <li>Run / Rawdata</li> <li>Elapsed run time is running</li> <li>Data are stored on PC and on card</li> </ul>	
LAN breaks	<ul style="list-style-type: none"> <li>Run in progress - Data Analysis</li> <li>Run / Rawdata</li> <li>Error Power Fail</li> <li>Elapsed run time stops</li> <li>Data continues to be stored on card</li> </ul>	
LAN recovers	<ul style="list-style-type: none"> <li>Run in progress - Data Analysis</li> <li>Run / Rawdata</li> <li>Error Power Fail cleared</li> <li>Elapsed run time continues at actual time</li> <li>Spectra counter continues</li> <li>Data continues to be stored on PC and on card</li> <li>ChemStation tries already to add missing data (depends on the data load).</li> </ul>	
Stop time elapsed	<ul style="list-style-type: none"> <li>Run in progress - Data Analysis</li> <li>Prerun / Rawdata</li> <li>Elapsed run time stops</li> <li>ChemStation continues to add missing data</li> </ul>	
Run ends	<ul style="list-style-type: none"> <li>Ready</li> <li>Run finished</li> <li>Prerun / Ready</li> </ul>	


**NOTE**

If the detector status window is not opened, you will realize only the Power Fail error and the long Run In Progress information until the data is recovered from disk.


## 4 Using the Detector

### Special Settings of the Detector


### Manual Run Recovery in case of permanent communication failures

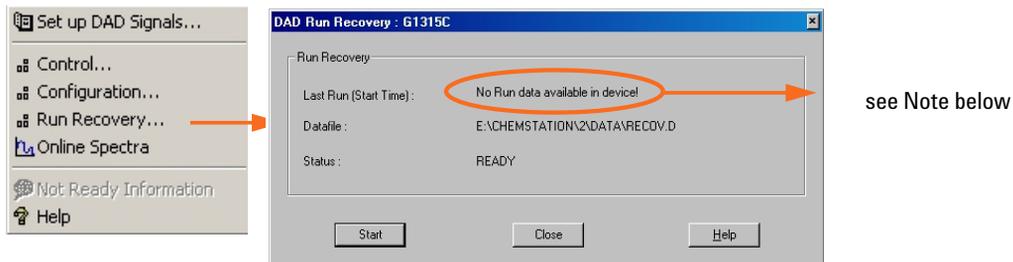

Start a recover

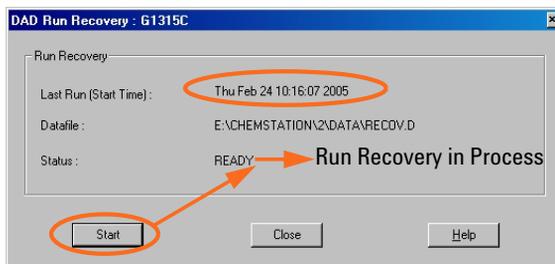


After a recover

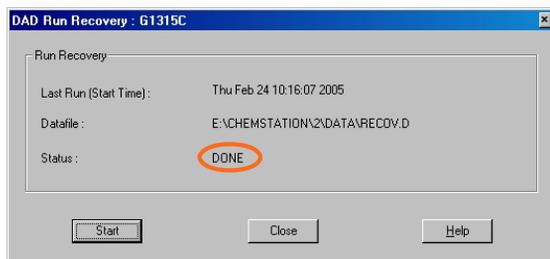

#### NOTE

When during recovery an error “Method/Sequence stopped” appears, the instrument logbook shows an entry “No Run data available in device”.

In this case refer to “No Run Data Available In Device” on page 112.

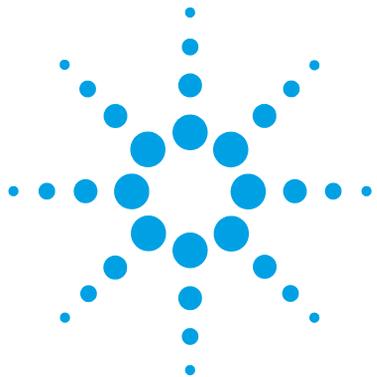

## 5 How to optimize the detector

Optimizing the Detector Performance 84

Match the Flow Cell to the Column 85

Set the Detector Parameters 88

This chapter gives hints on how to select the detector parameters and the flow cell.



## **Optimizing the Detector Performance**

The detector has a variety of parameters that can be used to optimize performance.

The information below will guide you on how to get the best detector performance. Follow these rules as a start for new applications. It gives a rule-of-thumb for optimizing the detector parameters.

## Match the Flow Cell to the Column

The tables below recommend the flow cell that matches the column used. If more than one selection is appropriate, use the larger flow cell to get the best detection limit. Use the smaller flow cell for best peak resolution.

### Standard HPLC Applications

Column length	Typical peak width	Recommended flow cell				
<= 5 cm	0.025 min	Micro flow cell	Semimicro flow cell	Standard flow cell	High Pressure flow cell	High Pressure flow cell
10 cm	0.05 min					
20 cm	0.1 min					
>= 40 cm	0.2 min					
	Typical flow rate	0.05-0.2 ml/min	0.2- 0.4 ml/min	0.4- 0.8 ml/min	1-2 ml/min	0.01- 5 ml/min
	Internal column diameter	1.0 mm	2.1mm	3.0 mm	4.6 mm	

**Figure 34** Choosing a Flow Cell (Standard HPLC Applications)

### Ultra fast separation with RRLC systems

Column ID	2.1 mm	3.0 mm	4.6 mm
Configuration	No damper No mixer ++	Damper Mixer	Damper Mixer
Flow cell	2 µl, 3 mm	5 µl, 6 mm +	14 µl, 10 mm +

**Figure 35** Choosing a Flow Cell for G1314E (for ultra fast separation with RRLC systems)

- (+) For ultra fast analysis with step gradients the micro flow cell (2 µL, 3 mm) gives the best performance
- (++) In high resolution analysis time is not the highest priority. Higher delay volumes are accepted. Therefore we recommend to use the damper plus mixer for a highest signal to noise.
- If longer columns (> 50 mm) for higher resolution are used, then the next larger flow cell is the preferred choice for higher sensitivity.

## Flow Cell Path Length

Lambert-Beer's law shows a linear relationship between the flow cell path length and absorbance.

$$\text{Absorbance} = -\log T = \log \frac{I_0}{I} = \epsilon \times C \times d$$

where

---


T	is the transmission, defined as the quotient of the intensity of the transmitted light I divided by the intensity of the incident light, I <sub>0</sub> ,
e	is the extinction coefficient, which is a characteristic of a given substance under a precisely-defined set of conditions of wavelength, solvent, temperature and other parameters,
C [mol/L]	is the concentration of the absorbing species,
d [m]	is the path length of the cell used for the measurement.

---

Therefore, flow cells with longer path lengths yield higher signals. Although noise usually increases little with increasing path length, there is a gain in the signal-to-noise ratio. For example, in [Figure 36](#) on page 87 the noise increased by less than 10 % but a 70 % increase in signal intensity was achieved by increasing the path length from 6 mm to 10 mm.

When increasing the path length, the cell volume usually increases – in the example from 5 – 14 µL. Typically, this causes more peak dispersion. As demonstrated, this did not affect the resolution in the gradient separation in the example that is shown below.

As a rule-of-thumb, the flow cell volume should be about 1/3 of the peak volume at half height. To determine the volume of your peaks, take the peak width as reported in the integration results multiply it by the flow rate and divide it by 3).

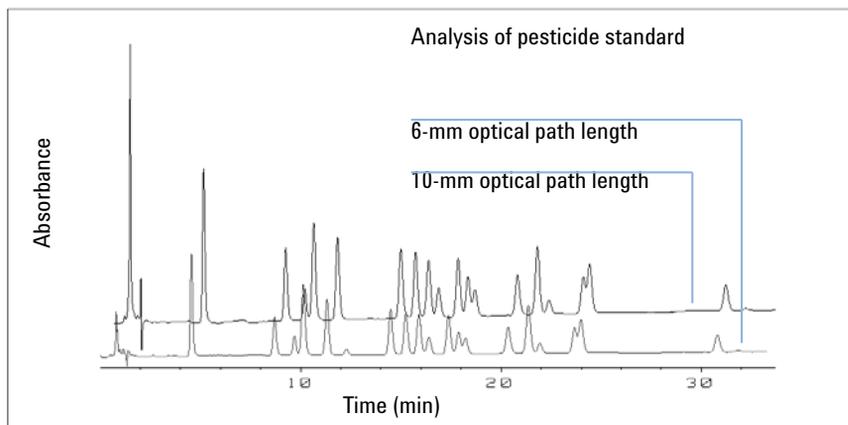

**Figure 36** Influence of Cell Path Length on Signal Height

Traditionally LC analysis with UV detectors is based on comparing measurements with internal or external standards. To check photometric accuracy of the detector, it is necessary to have more precise information on path lengths of the flow cells.

The correct response is:

expected response \* correction factor

Please find below the details of the flow cells:

**Table 14** Correction factors for Agilent VWD flow cells


Part number	Path length (actual)	Correction factor
Standard flow cell 10 mm, 14 µL, 40 bar (G1314-60186)	10.15 ± 0.19 mm	10/10.15
Semi-micro flow cell 6 mm, 5 µL (G1314-60183)	6.10 ± 0.19 mm	6/6.10
Micro flow cell 3 mm, 2 µL, 120 bar (G1314-60187)	2.80 ± 0.19 mm	3/2.8
High pressure flow cell 10 mm, 14 µL, 400 bar (G1314-60182)	10.00 ± 0.19 mm	10/10

**NOTE**

However you have to be aware that there is additional tolerance of gasket thickness and its compression ratio which is supposed to be very small in comparison with the machining tolerance.

## Set the Detector Parameters

- 1** Set peakwidth as close as possible to the width (at half height) of a narrow peak of interest.
  
- 2** Choose the sample wavelength.
  - at a longer wavelength than the cut-off wavelength of the mobile phase,
  - at a wavelength where the analytes have strong absorptivity if you want to get the lowest possible detection limit,
  - at a wavelength with moderate absorptivity if you work with high concentrations, and
  - preferably where the spectrum is flat for better linearity.
  
- 3** Consider to use time-programming to further optimization.

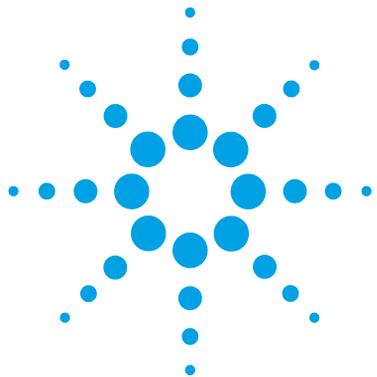

## 6 Troubleshooting and Diagnostics

Overview of the Detector's Indicators and Test Functions	90
Status Indicators	91
Power Supply Indicator	91
Module Status Indicator	92
Available Tests versus Interfaces	93
Agilent Lab Advisor Software	94

Overview about the troubleshooting and diagnostic features.



## Overview of the Detector's Indicators and Test Functions

### Status Indicators

The detector is provided with two status indicators which indicate the operational state (prerun, run, and error states) of the detector. The status indicators provide a quick visual check of the operation of the detector "[Status Indicators](#)" on page 91.

### Error Messages

In the event of an electronic, mechanical or hydraulic failure, the detector generates an error message in the user interface. For each message, a short description of the failure, a list of probable causes of the problem, and a list of suggested actions to fix the problem are provided by the user interface. Refer to the Service Manual for details.

### Test Functions

A series of test functions are available for troubleshooting and operational verification after exchanging internal components. Refer to the User Interface and/or the Service Manual for details.

### Wavelength Verification / Recalibration


Wavelength recalibration is recommended after repair of internal components, and on a regular basis to ensure correct operation of the detector. The detector uses the deuterium alpha and beta emission lines for wavelength calibration, see "[Wavelength Verification-Calibration](#)" on page 120.

### Diagnostic Signals

The detector has several signals (internal temperatures, voltages and currents of lamps) that can be used for diagnosing baseline problems. Refer to the Service Manual for details.

## Status Indicators

Two status indicators are located on the front of the detector. The lower left indicates the power supply status, the upper right indicates the detector status.

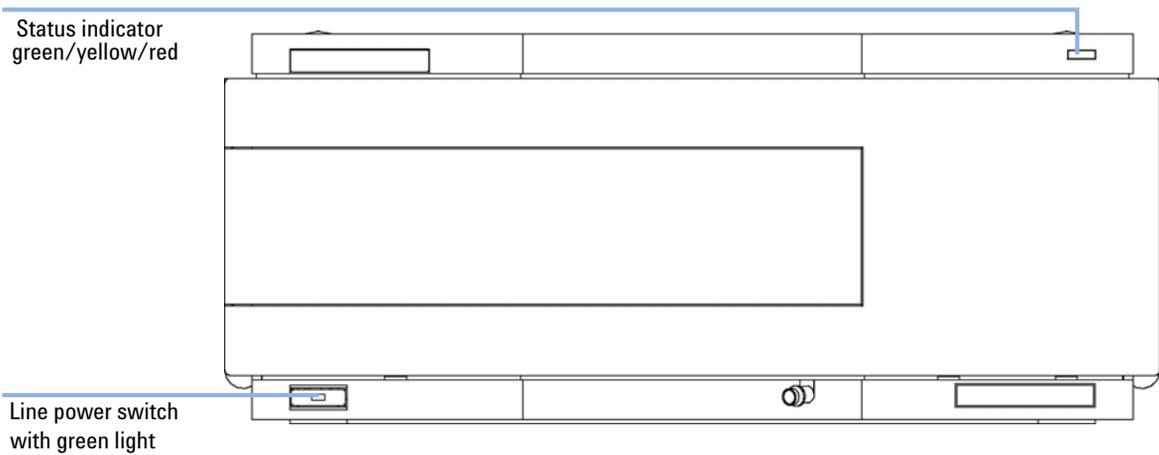

**Figure 37** Location of Status Indicators

## Power Supply Indicator

The power supply indicator is integrated into the main power switch. When the indicator is illuminated (*green*) the power is ON.

## Module Status Indicator

The module status indicator indicates one of six possible module conditions:

- When the status indicator is *OFF* (and power switch light is on), the module is in a *prerun* condition, and is ready to begin an analysis.
- A *green* status indicator, indicates the module is performing an analysis (*run* mode).
- A *yellow* indicator indicates a *not-ready* condition. The module is in a not-ready state when it is waiting for a specific condition to be reached or completed (for example, immediately after changing a set point), or while a self-test procedure is running.
- An *error* condition is indicated when the status indicator is *red*. An error condition indicates the module has detected an internal problem which affects correct operation of the module. Usually, an error condition requires attention (e.g. leak, defective internal components). An error condition always interrupts the analysis.

If the error occurs during analysis, it is propagated within the LC system, i.e. a red LED may indicate a problem of a different module. Use the status display of your user interface for finding the root cause/module of the error.

- A *blinking* indicator indicates that the module is in resident mode (e.g. during update of main firmware).
- A *fast blinking* indicator indicates that the module is in a low-level error mode. In such a case try to re-boot the module or try a cold-start (see “[Special Settings](#)” on page 191). Then try a firmware update (see “[Replacing the Module’s Firmware](#)” on page 147). If this does not help, a main board replacement is required.

## Available Tests versus Interfaces

**NOTE**

Depending on the used interface, the available tests and the screens/reports may vary. Preferred tool should be the Agilent Lab Advisor, see “[Agilent Lab Advisor Software](#)” on page 94.

In future, a user interface may not show the Diagnostics/Tests anymore. Then the Agilent Lab Advisor must be used instead.

The Agilent ChemStation may not include any maintenance/test functions.

**Table 15** Available Tests versus Interfaces

Interface Test	Agilent Lab Advisor	Agilent ChemStation	Instant Pilot G4208A
Wavelength Verification/Re-calibration	yes (*)	Tests (*)	Maintenance (*)
Lamp Intensity	yes (*)	Tests (*)	Diagnosis (*)
Holmium Test	yes (*)	Tests (*)	Diagnosis (*)
Cell Test	yes (*)	Tests (*)	n/a
D/A Converter Test	yes (*)	Tests (*)	n/a
Filter / Grating Motor Test	yes	Tests (*)	command line (***)
Test Chromatogram	yes	command line (**)	command line (***)
Spectrum (Blank, Sample, Holmium)	yes	n/a	Control
Service Dialog	for service only	n/a	for service only

(\*) interface provides passed/fail information or a plot.

(\*\*) requires a comand via comand line

(\*\*\*) requires a comand via comand line in Service Mode


## Agilent Lab Advisor Software

The Agilent Lab Advisor software is a standalone product that can be used with or without data system. Agilent Lab Advisor software helps to manage the lab for high quality chromatographic results and can monitor in real time a single Agilent LC or all the Agilent GCs and LCs configured on the lab intranet.

Agilent Lab Advisor software provides diagnostic capabilities for all Agilent 1200 Infinity Series modules. This includes diagnostic capabilities, calibration procedures and maintenance routines for all the maintenance routines.

The Agilent Lab Advisor software also allows users to monitor the status of their LC instruments. The Early Maintenance Feedback (EMF) feature helps to carry out preventive maintenance. In addition, users can generate a status report for each individual LC instrument. The tests and diagnostic features as provided by the Agilent Lab Advisor software may differ from the descriptions in this manual. For details refer to the Agilent Lab Advisor software help files.

The Instrument Utilities is a basic version of the Lab Advisor with limited functionality required for installation, use and maintenance. No advanced repair, troubleshooting and monitoring functionality is included.

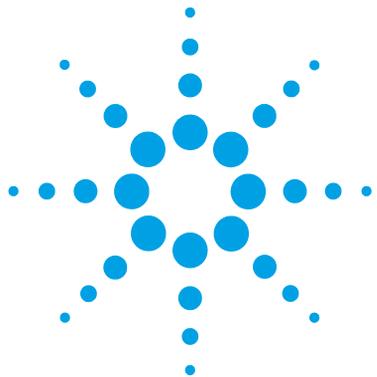

## 7 Error Information

What Are Error Messages	97
General Error Messages	98
Timeout	98
Shutdown	98
Remote Timeout	99
Lost CAN Partner	100
Leak	100
Leak Sensor Open	101
Leak Sensor Short	101
Compensation Sensor Open	102
Compensation Sensor Short	102
Fan Failed	103
Open Cover	103
Detector Error Messages	104
UV lamp: no current	104
UV lamp: no voltage	105
Ignition Failed	105
No heater current	106
Wavelength calibration setting failed	106
Wavelength holmium check failed	107
Grating or Filter Motor Errors	108
Wavelength test failed	109
Cutoff filter doesn't decrease the light intensity at 250 nm	109
ADC Hardware Error	110
Illegal Temperature Value from Sensor on Main Board	110
Illegal Temperature Value from Sensor at Air Inlet	111
Heater at fan assembly failed	111
Heater Power At Limit	112



## **7 Error Information**

Agilent Lab Advisor Software

No Run Data Available In Device [112](#)

Cover Violation [113](#)

This chapter describes the meaning of detector error messages, and provides information on probable causes and suggested actions how to recover from error conditions.

## What Are Error Messages

Error messages are displayed in the user interface when an electronic, mechanical, or hydraulic (flow path) failure occurs which requires attention before the analysis can be continued (for example, repair, or exchange of consumables is necessary). In the event of such a failure, the red status indicator at the front of the module is switched on, and an entry is written into the module logbook.

If an error occurs outside a method run, other modules will not be informed about this error. If it occurs within a method run, all connected modules will get a notification, all LEDs get red and the run will be stopped. Depending on the module type, this stop is implemented differently. For example, for a pump the flow will be stopped for safety reasons. For a detector, the lamp will stay on in order to avoid equilibration time. Depending on the error type, the next run can only be started, if the error has been resolved, for example liquid from a leak has been dried. Errors for presumably single time events can be recovered by switching on the system in the user interface.

Special handling is done in case of a leak. As a leak is a potential safety issue and may have occurred at a different module from where it has been observed, a leak always causes a shutdown of all modules, even outside a method run.

In all cases, error propagation is done via the CAN bus or via an APG remote cable (see documentation for the APG interface).

## General Error Messages

General error messages are generic to all Agilent series HPLC modules and may show up on other modules as well.

### Timeout

**Error ID: 0062**

The timeout threshold was exceeded.

**Probable cause**

- 1** The analysis was completed successfully, and the timeout function switched off the module as requested.
- 2** A not-ready condition was present during a sequence or multiple-injection run for a period longer than the timeout threshold.

**Suggested actions**

- Check the logbook for the occurrence and source of a not-ready condition. Restart the analysis where required.
- Check the logbook for the occurrence and source of a not-ready condition. Restart the analysis where required.

### Shutdown

**Error ID: 0063**

An external instrument has generated a shutdown signal on the remote line.

The module continually monitors the remote input connectors for status signals. A LOW signal input on pin 4 of the remote connector generates the error message.

<b>Probable cause</b>	<b>Suggested actions</b>
<p><b>1</b> Leak detected in another module with a CAN connection to the system.</p>	<p>Fix the leak in the external instrument before restarting the module.</p>
<p><b>2</b> Leak detected in an external instrument with a remote connection to the system.</p>	<p>Fix the leak in the external instrument before restarting the module.</p>
<p><b>3</b> Shut-down in an external instrument with a remote connection to the system.</p>	<p>Check external instruments for a shut-down condition.</p>
<p><b>4</b> The degasser failed to generate sufficient vacuum for solvent degassing.</p>	<p>Check the vacuum degasser for an error condition. Refer to the <i>Service Manual</i> for the degasser or the 1260 pump that has the degasser built-in.</p>

## Remote Timeout

### Error ID: 0070

A not-ready condition is still present on the remote input. When an analysis is started, the system expects all not-ready conditions (for example, a not-ready condition during detector balance) to switch to run conditions within one minute of starting the analysis. If a not-ready condition is still present on the remote line after one minute the error message is generated.

<b>Probable cause</b>	<b>Suggested actions</b>
<p><b>1</b> Not-ready condition in one of the instruments connected to the remote line.</p>	<p>Ensure the instrument showing the not-ready condition is installed correctly, and is set up correctly for analysis.</p>
<p><b>2</b> Defective remote cable.</p>	<p>Exchange the remote cable.</p>
<p><b>3</b> Defective components in the instrument showing the not-ready condition.</p>	<p>Check the instrument for defects (refer to the instrument's documentation).</p>

## Lost CAN Partner

### Error ID: 0071

During an analysis, the internal synchronization or communication between one or more of the modules in the system has failed.

The system processors continually monitor the system configuration. If one or more of the modules is no longer recognized as being connected to the system, the error message is generated.

#### Probable cause

- 1 CAN cable disconnected.
- 2 Defective CAN cable.
- 3 Defective main board in another module.

#### Suggested actions

- Ensure all the CAN cables are connected correctly.
  - Ensure all CAN cables are installed correctly.
- Exchange the CAN cable.
- Switch off the system. Restart the system, and determine which module or modules are not recognized by the system.

## Leak

### Error ID: 0064

A leak was detected in the module.

The signals from the two temperature sensors (leak sensor and board-mounted temperature-compensation sensor) are used by the leak algorithm to determine whether a leak is present. When a leak occurs, the leak sensor is cooled by the solvent. This changes the resistance of the leak sensor which is sensed by the leak-sensor circuit on the main board.

#### Probable cause

- 1 Loose fittings.
- 2 Broken capillary.
- 3 Leaking flow cell.

#### Suggested actions

- Ensure all fittings are tight.
- Exchange defective capillaries.
- Exchange flow cell components.

## Leak Sensor Open

**Error ID: 0083**

The leak sensor in the module has failed (open circuit).

The current through the leak sensor is dependent on temperature. A leak is detected when solvent cools the leak sensor, causing the leak-sensor current to change within defined limits. If the current falls outside the lower limit, the error message is generated.

<b>Probable cause</b>	<b>Suggested actions</b>
<ol style="list-style-type: none"> <li>1 Leak sensor not connected to the main board.</li> </ol>	<p>Please contact your Agilent service representative.</p>
<ol style="list-style-type: none"> <li>2 Defective leak sensor.</li> </ol>	<p>Please contact your Agilent service representative.</p>
<ol style="list-style-type: none"> <li>3 Leak sensor incorrectly routed, being pinched by a metal component.</li> </ol>	<p>Please contact your Agilent service representative.</p>

## Leak Sensor Short

**Error ID: 0082**

The leak sensor in the module has failed (short circuit).

The current through the leak sensor is dependent on temperature. A leak is detected when solvent cools the leak sensor, causing the leak sensor current to change within defined limits. If the current increases above the upper limit, the error message is generated.

<b>Probable cause</b>	<b>Suggested actions</b>
<ol style="list-style-type: none"> <li>1 Defective leak sensor.</li> </ol>	<p>Please contact your Agilent service representative.</p>
<ol style="list-style-type: none"> <li>2 Leak sensor incorrectly routed, being pinched by a metal component.</li> </ol>	<p>Please contact your Agilent service representative.</p>

## Compensation Sensor Open

**Error ID: 0081**

The ambient-compensation sensor (NTC) on the main board in the module has failed (open circuit).

The resistance across the temperature compensation sensor (NTC) on the main board is dependent on ambient temperature. The change in resistance is used by the leak circuit to compensate for ambient temperature changes. If the resistance across the sensor increases above the upper limit, the error message is generated.

**Probable cause**

- 1 Defective main board.

**Suggested actions**

Please contact your Agilent service representative.

## Compensation Sensor Short

**Error ID: 0080**

The ambient-compensation sensor (NTC) on the main board in the module has failed (open circuit).

The resistance across the temperature compensation sensor (NTC) on the main board is dependent on ambient temperature. The change in resistance is used by the leak circuit to compensate for ambient temperature changes. If the resistance across the sensor falls below the lower limit, the error message is generated.

**Probable cause**

- 1 Defective main board.

**Suggested actions**

Please contact your Agilent service representative.

## Fan Failed

**Error ID: 0068**

The cooling fan in the module has failed.

The hall sensor on the fan shaft is used by the main board to monitor the fan speed. If the fan speed falls below a certain limit for a certain length of time, the error message is generated.

This limit is given by 2 revolutions/second for longer than 5 seconds.

Depending on the module, assemblies (e.g. the lamp in the detector) are turned off to assure that the module does not overheat inside.

**Probable cause**

- 1 Fan cable disconnected.
- 2 Defective fan.
- 3 Defective main board.

**Suggested actions**

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## Open Cover

**Error ID: 0205**

The top foam has been removed.

The sensor on the main board detects when the top foam is in place. If the foam is removed, the fan is switched off, and the error message is generated.

**Probable cause**

- 1 The top foam was removed during operation.
- 2 Foam not activating the sensor.
- 3 Defective sensor or main board.

**Suggested actions**

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## Detector Error Messages

These errors are detector specific.

### UV lamp: no current

**Error ID: 7450**

The lamp anode current is missing. The processor continually monitors the anode current drawn by the lamp during operation. If the anode current falls below the lower current limit, the error message is generated.

**Probable cause**

- 1** Lamp disconnected.
- 2** Top foam removed while lamp is on.
- 3** Defective or non-Agilent lamp.
- 4** Defective main board.
- 5** Defective power supply.

**Suggested actions**

- Ensure the lamp connector is seated firmly.
- Please contact your Agilent service representative.
- Exchange the lamp.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## UV lamp: no voltage

**Error ID: 7451**

The lamp anode voltage is missing. The processor continually monitors the anode voltage across the lamp during operation. If the anode voltage falls below the lower limit, the error message is generated.

**Probable cause**

- 1 Defective or non-Agilent lamp.
- 2 Defective power supply.
- 3 Defective main board.

**Suggested actions**

- Exchange the lamp.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## Ignition Failed

**Error ID: 7452**

The lamp failed to ignite. The processor monitors the lamp current during the ignition cycle. If the lamp current does not rise above the lower limit within 2 – 5 s, the error message is generated.

**Probable cause**

- 1 Lamp disconnected.
- 2 Defective or non-Agilent lamp.
- 3 Defective power supply.
- 4 Defective main board.

**Suggested actions**

- Ensure the lamp is connected.
- Exchange the lamp.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## No heater current

**Error ID: 7453**

The lamp heater current in the detector is missing. During lamp ignition, the processor monitors the heater current. If the current does not rise above the lower limit within 1 , the error message is generated.

**Probable cause**

- 1 Lamp disconnected.
- 2 Ignition started without the top foam in place.
- 3 Defective main board.
- 4 Defective or non-Agilent lamp.
- 5 Defective power supply.

**Suggested actions**

- Ensure the lamp is connected.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Exchange the lamp.
- Please contact your Agilent service representative.

## Wavelength calibration setting failed

**Error ID: 7310**

The intensity maximum was not found during wavelength calibration.

**Calibration 0 Failed:**

Zero-order calibration failed.

**Calibration 1 Failed:**

656 nm calibration failed.

**Probable cause**

- 1 Lamp is OFF.
- 2 Incorrect flow cell installation.
- 3 Flow cell contamination or air bubbles.

**Suggested actions**

- Switch on the lamp.
- Ensure the flow cell is installed correctly.
- Clean/replace flow cell windows or remove air bubbles.

<b>Probable cause</b>	<b>Suggested actions</b>
<b>4</b> Intensity too low.	Replace lamp.
<b>5</b> Current step value too far from maximum.	<ul style="list-style-type: none"> <li>• Repeat the calibration.</li> <li>• Please contact your Agilent service representative.</li> </ul>
<b>6</b> Misaligned/defective grating assembly.	Please contact your Agilent service representative.
<b>7</b> Defective main board.	Please contact your Agilent service representative.

## Wavelength holmium check failed

### Error ID: 7318

The holmium oxide test in the detector has failed. During the holmium test, the detector moves the holmium filter into the light path, and compares the measured absorbance maxima of the holmium oxide filter with expected maxima. If the measured maxima are outside the limits, the error message is generated.

<b>Probable cause</b>	<b>Suggested actions</b>
<b>1</b> Misaligned/defective grating assembly.	<ul style="list-style-type: none"> <li>• Ensure the flow cell is inserted correctly, and is free from contamination (cell windows, buffers, and so on).</li> <li>• Run the filter-motor test to determine if the filter motor assembly is defective. If defective, please contact your Agilent service representative.</li> <li>• Run the grating-motor test to determine if the grating assembly is defective. If defective, please contact your Agilent service representative.</li> </ul>

## Grating or Filter Motor Errors

**Error ID:** Grating: 7800, 7801, 7802, 7803, 7804, 7805, 7806, 7808, 7809;  
Filter: 7810, 7811, 7812, 7813, 7814, 7815, 7816

The motor test has failed.

<b>Test 0 Failed:</b>	Filter motor.
<b>Test 1 Failed:</b>	Grating motor.

During the motor tests, the detector moves the motor to the end position while monitoring the end-position sensor. If the end position is not found, the error message is generated.

<b>Probable cause</b>	<b>Suggested actions</b>
<b>1</b> Motor is not connected.	Please contact your Agilent service representative.
<b>2</b> Defective motor.	Please contact your Agilent service representative.
<b>3</b> Defective/missing grating or filter.	Please contact your Agilent service representative.
<b>4</b> Cable/connector defective.	Please contact your Agilent service representative.

## Wavelength test failed

### Error ID: 7890

The automatic wavelength check after lamp ignition has failed. When the lamp is switched on, the detector waits 1 min to warm-up the lamp. Then a check of the deuterium emission line (656 nm) via the reference diode is performed. If the emission line is more than 3 nm away from 656 nm, the error message is generated.

#### Probable cause

- 1 Calibration incorrect.

#### Suggested actions

Recalibrate the detector.

## Cutoff filter doesn't decrease the light intensity at 250 nm

### Error ID: 7813

The automatic filter check after lamp ignition has failed. When the lamp is switched on, the detector moves the cutoff filter into the light path. If the filter is functioning correctly, a decrease in lamp intensity is seen. If the expected intensity decrease is not detected, the error message is generated.

#### Probable cause

- 1 Motor is not connected.
- 2 Defective motor.
- 3 Defective/missing grating or filter.
- 4 Cable/connector defective.

#### Suggested actions

Please contact your Agilent service representative.

## ADC Hardware Error

**Error ID: 7830, 7831**

A/D-Converter hardware is defective.

**Probable cause**

- 1 A/D-Converter hardware is defective.

**Suggested actions**

Please contact your Agilent service representative.

## Illegal Temperature Value from Sensor on Main Board

**Error ID: 1071**

This temperature sensor (located on the detector main board) delivered a value outside the allowed range. The parameter of this event equals the measured temperature in 1/100 centigrade. As a result the temperature control is switched off.

**Probable cause**

- 1 Defective sensor or main board.
- 2 Detector is exposed to illegal ambient conditions.

**Suggested actions**

Please contact your Agilent service representative.

Verify that the ambient conditions are within the allowed range.

## Illegal Temperature Value from Sensor at Air Inlet

**Error ID: 1072**

This temperature sensor delivered a value outside the allowed range. The parameter of this event equals the measured temperature in 1/100 centigrade. As a result the temperature control is switched off.

**Probable cause**

- 1** The temperature sensor is defect.
  
- 2** Detector is exposed to illegal ambient conditions.

**Suggested actions**

- Replace the cable to the main board.
  - Please contact your Agilent service representative.
- Verify that the ambient conditions are within the allowed range.

## Heater at fan assembly failed

**Error ID: 1073**

Every time the deuterium lamp or the tungsten lamp (DAD only) is switched on or off a heater self-test is performed. If the test fails an error event is created. As a result the temperature control is switched off.

**Probable cause**

- 1** Defective connector or cable.
  
- 2** Defective heater.

**Suggested actions**

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## Heater Power At Limit

**Error ID: 1074**

The available power of the heater reached either the upper or lower limit. This event is sent only once per run. The parameter determines which limit has been hit:

0 means upper power limit hit (excessive ambient temperature drop).

1 means lower power limit hit (excessive ambient temperature increase).

**Probable cause**

- 1 Excessive ambient temperature change.

**Suggested actions**

Wait until temperature control equilibrates.

## No Run Data Available In Device

In a very rare case the capacity of the CompactFlash Card is not sufficient. This could happen for example when the interrupt of LAN communication takes longer and the detector uses special settings (e.g full data rate at 80 Hz plus full spectra plus all signals) during data buffering.

**Probable cause**

- 1 CompactFlash Card is full.

**Suggested actions**

- Correct communication problem.
- Reduce data rate.

## Cover Violation

**Error ID: 7461**

The top foam has been removed.

The sensor on the main board detects when the top foam is in place. If the foam is removed while the lamps are on (or if an attempt is made to switch on for example the lamps with the foam removed), the lamps are switched off, and the error message is generated.

**Probable cause**


- 1 The top foam was removed during operation.
- 2 Foam not activating the sensor.

**Suggested actions**

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

## **7 Error Information**

### **Detector Error Messages**

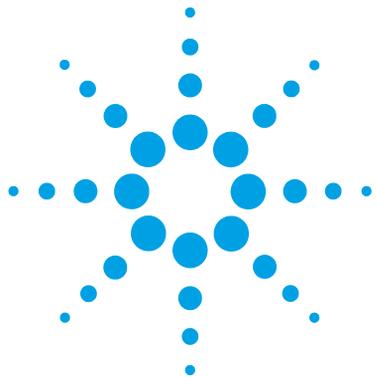

## 8 Test Functions

Intensity Test	116
Intensity Test Failed	117
Cell Test	118
Wavelength Verification-Calibration	120
ASTM Drift and Noise Test	122
Quick Noise Test	123
Dark Current Test	124
Dark Current Test Failed	125
Holmium Oxide Test	126
Holmium Oxide Test Failed	127

This chapter describes the detector's built in test functions.



## Intensity Test

The intensity test measures the intensity of the deuterium lamp over the full VWD wavelength range (190 - 600 nm). The test can be used to determine the performance of the lamp, and to check for dirty or contaminated flow cell windows. When the test is started, the gain is set to zero. To eliminate effects due to absorbing solvents, the test should be done with water in the flow cell. The shape of the intensity spectrum is primarily dependent on the lamp, grating, and diode characteristics. Therefore, intensity spectra will differ slightly between instruments. The figure below shows a typical intensity test spectrum.

The Intensity Test is available in

- Agilent Lab Advisor (preferred tool).
- Agilent Instant Pilot G4208A, via **More-Diagnosis-VWD-Lamp Intensity Test**.

### Intensity Test Evaluation

The Agilent Lab Advisor and the Instant Pilot evaluate three values automatically and display the limits for each value, the average, the minimum and the maximum of all data points and **passed** or **failed** for each value.



## Cell Test

The cell test compares the intensity of the deuterium lamp measured by the sample and reference diodes (unfiltered and not logarithmized) when the grating is in the zero-order position. The resulting intensity ratio (sample:reference) is a measure of the amount of light absorbed by the flow cell.

The test can be used to check for dirty or contaminated flow cell windows. When the test is started, the gain is set to -1. To eliminate effects due to absorbing solvents, the test should be done with water in the flow cell.

Limits: No real limit. The reason is that it depends on the position/alignment of the reference side (beam splitter – reference slit – reference diode). Therefore the reference side value can be higher/smaller than the sample side value.

With a clean cell the counts for sample and reference (photocurrent) are in the same range. If the sample side shows much lower values than the reference side the flow cell might have a problem.

Pre-requisite:

Flush the flow cell with a flow of 1 mL/min for at least 10 minutes.

Probable Cause	Suggested Action
Cell contaminated	Flush flow cell
Cell windows are contaminated	Clean/replace cell windows
Mechanical problem	Check cell position

In the Agilent Instant Pilot G4208A, the photocurrent readings are available via **More > Diagnosis > VWD > LampIntensity Test**, see [Figure 40](#) on page 119.

<b>Test Name</b>	Cell Test	<b>Description</b>	Calculate the ratio of the sample signal and the reference signal, measured in the zero order of the grating.
<b>Module</b>	G1314C:DE60555128		
<b>Status</b>	Passed		
<b>Start Time</b>	7/6/2011 1:24:55 PM		
<b>Stop Time</b>	7/6/2011 1:26:18 PM		

Test Procedure

- ✓ 1. Check Prerequisites...
- ✓ 2. Flush Flow Cell.
- ✓ 3. Measure Sample and Reference Intensity...
- ✓ 4. Evaluate Data...

Result	
Name	Value
Accumulated UV Lamp Burn Time	60.49 h
UV Lamp On-Time	4.36 h
Intensity Sample	241,908 Counts
Intensity Reference	422,625 Counts
Intensity Ratio	0.57

Figure 39 Cell Test with Lab Advisor

### Checking the Photocurrent with the Instant Pilot

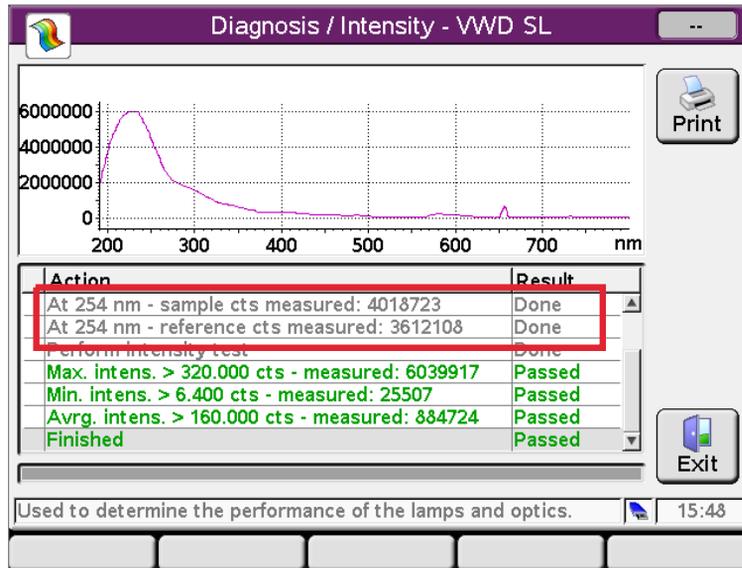


Figure 40 Checking the Photocurrent with the Instant Pilot

## Wavelength Verification-Calibration

Wavelength calibration of the detector is done using the zero-order position and 656 nm emission line position of the deuterium lamp. The calibration procedure involves two steps. First the grating is calibrated on the zero-order position. The stepper-motor step position where the zero-order maximum is detected is stored in the detector. Next, the grating is calibrated against the deuterium emission-line at 656 nm, and the motor position at which the maximum occurs is stored in the detector.

In addition to the zero-order and 656 nm (alpha-emission line) calibration, the beta-emission line at 486 nm and the three holmium lines are used for the complete wavelength calibration process. These holmium lines are at 360.8 nm, 418.5 nm and 536.4 nm.

### NOTE

The wavelength verification/calibration takes about 2.5 min and is disabled within the first 10 min after ignition of the lamp because initial drift may distort the measurement.

When the lamp is turned **ON**, the 656 nm emission line position of the deuterium lamp is checked automatically.

The Wavelength Verification/Calibration is available in

- Agilent Lab Advisor (preferred tool).
- Agilent Instant Pilot G4208A, via **More-Diagnosis-VWD-Calibration**.

### When to Calibrate the Detector

The detector is calibrated at the factory, and under normal operating conditions should not require recalibration. However, it is advisable to recalibrate:

- after maintenance (flow cell or lamp),
- after repair of components in the optical unit,
- after exchange of the optical unit or VWM board,
- at a regular interval, at least once per year (for example, prior to an Operational Qualification/Performance Verification procedure), and
- when chromatographic results indicate the detector may require recalibration.






## Quick Noise Test

The noise test measures the noise of the detector, with HPLC-grade water flowing through the flow cell at 1 mL/min, in one minute intervals over a total of 5 minutes.

The noise of the detector is calculated by using the maximum amplitude for all random variations of the detector signal of frequencies greater than one cycle per hour. The noise is determined for 5 one minute intervals and is based on the accumulated peak-to-peak noise for the intervals. At least seven data points per cycles are used in the calculation.

The cycles in the noise determination are not overlapping.


In order to obtain reliable results, the lamp should be turned on for at least 10 minutes prior to measurement.

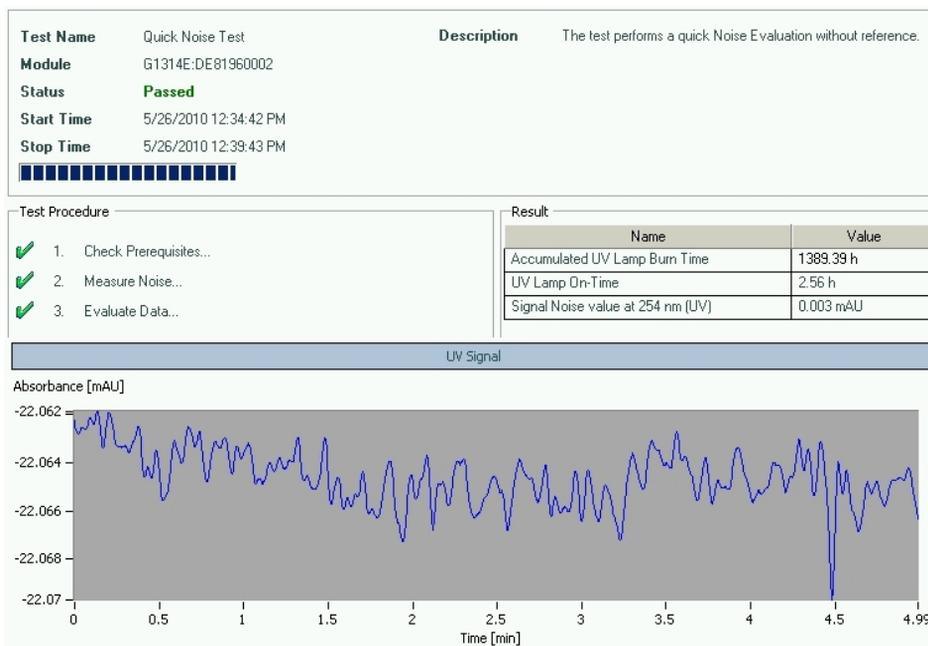

**Figure 43** Quick Noise Test with Agilent Lab Advisor

## Dark Current Test

The dark-current test measures the leakage current from the sample and reference circuits. The test is used to check for defective sample or reference diodes or ADC circuits which may cause non-linearity or excessive baseline noise. During the test, the shutter is moved into the light path. Next, the leakage current from both diodes is measured.

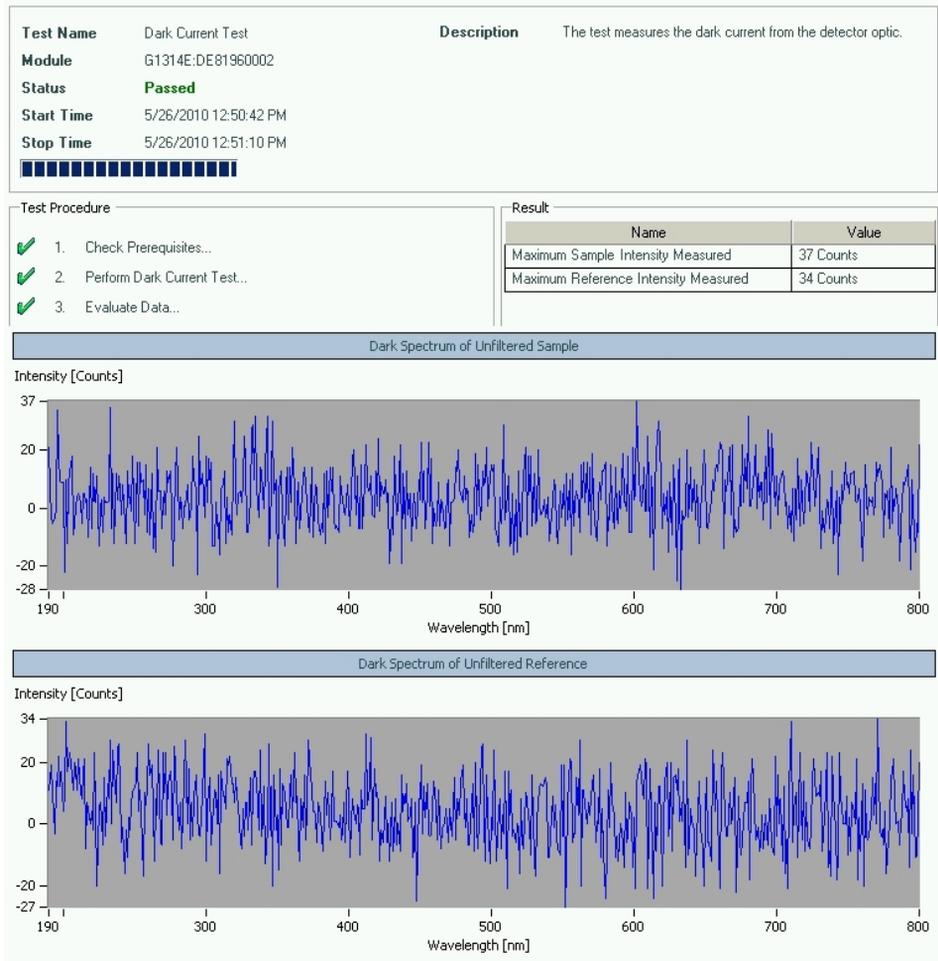

**Figure 44** Dark Current Test with Agilent Lab Advisor

## Dark Current Test Failed

### Probable cause

- 1 Defective sample or reference diode.
- 2 Defective sample or reference ADC board.
- 3 Defective main board.

### Suggested actions

Please contact your Agilent service representative.

Please contact your Agilent service representative.

Please contact your Agilent service representative.

## Holmium Oxide Test

This test verifies the calibration of the detector against the three wavelength maxima of the built-in holmium oxide filter. The test displays the difference between the expected and measured maxima. The figure below shows a holmium test spectrum.

The Holmium Oxide Test is available in

- Agilent Lab Advisor (preferred tool).
- Agilent Instant Pilot G4208A, via **More-Diagnosis-VWD-Holmium Spectrum Test**.

The test uses the following holmium maxima:

- 360.8 nm
- 418.5 nm
- 536.4 nm

### NOTE

See also “[Declaration of Conformity for HOX2 Filter](#)” on page 236.

### When to do the Test

- after recalibration,
- as part of the Operational Qualification/Performance Verification procedure, or
- after flow cell maintenance or repair.


### Interpreting the Results

The test is passed successfully when all three wavelengths are within  $\pm 1$  nm of the expected value. This indicates the detector is calibrated correctly.

## Holmium Oxide Test with Agilent Lab Advisor

<b>Test Name</b>	Holmium Oxide Test	<b>Description</b>	The test measures the Holmium spectrum from the built-in Holmium filter. The spectrum is evaluated for peaks at different wavelengths.
<b>Module</b>	G1314E:DE81960002		
<b>Status</b>	<b>Passed</b>		
<b>Start Time</b>	5/28/2010 10:28:23 AM		
<b>Stop Time</b>	5/28/2010 10:31:45 AM		

<b>Test Procedure</b>		<b>Result</b>	
✓	1. Check Prerequisites...	<b>Name</b>	<b>Value</b>
✓	2. Perform Holmium Oxide Test...	Accumulated UV Lamp Burn Time	1395.27 h
✓	3. Evaluate Data...	UV Lamp On-Time	0.63 h
		Holmium Deviation to 360.8 nm	0.00 nm
		Holmium Deviation to 418.5 nm	-0.50 nm
		Holmium Deviation to 536.4 nm	-0.10 nm

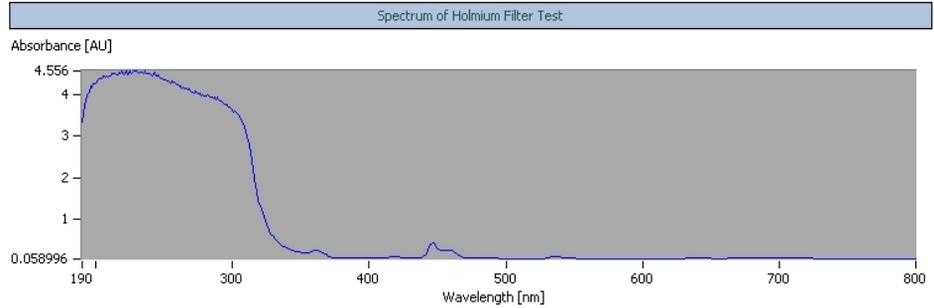

**Figure 45** Holmium Oxide Test with Agilent Lab Advisor

## Holmium Oxide Test Failed

### Probable cause


- 1 Detector not calibrated.
- 2 Dirty or defective flow cell.
- 3 Dirty or defective holmium oxide filter.
- 4 Optical misalignment.

### Suggested actions

- Recalibrate the detector.
- Repeat the test with the flow cell removed. If the test is OK, exchange the flow cell components.
- Run the holmium oxide filter test. If the test fails, contact your Agilent service representative.
- Please contact your Agilent service representative.

## **8 Test Functions**

### **Holmium Oxide Test**

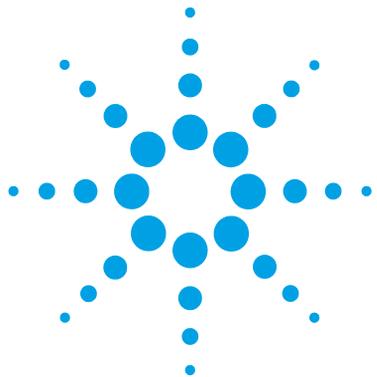

## 9 Maintenance and Repair

Introduction to Maintenance	130
Warnings and Cautions	131
Overview of Maintenance	133
Cleaning the Module	134
Exchanging a Lamp	135
Exchanging a Flow Cell	138
Repairing the Flow Cells	140
Using the Cuvette Holder	143
Correcting Leaks	145
Replacing Leak Handling System Parts	146
Replacing the Module's Firmware	147

This chapter provides general information on maintenance and repair of the detector.



## Introduction to Maintenance

The module is designed for easy maintenance. Maintenance can be done from the front with module in place in the system stack.

**NOTE**

There are no serviceable parts inside.  
Do not open the module.

---

## Warnings and Cautions

### WARNING

#### Toxic, flammable and hazardous solvents, samples and reagents

The handling of solvents, samples and reagents can hold health and safety risks.

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
  - The volume of substances should be reduced to the minimum required for the analysis.
  - Do not operate the instrument in an explosive atmosphere.
- 

### WARNING

#### Eye damage by detector light



Eye damage may result from directly viewing the UV-light produced by the lamp of the optical system used in this product.

- Always turn the lamp of the optical system off before removing it.
- 

### WARNING

#### Electrical shock

Repair work at the module can lead to personal injuries, e.g. shock hazard, when the cover is opened.

- Do not remove the cover of the module.
  - Only certified persons are authorized to carry out repairs inside the module.
-

**WARNING**

**Personal injury or damage to the product**

**Agilent is not responsible for any damages caused, in whole or in part, by improper use of the products, unauthorized alterations, adjustments or modifications to the products, failure to comply with procedures in Agilent product user guides, or use of the products in violation of applicable laws, rules or regulations.**

→ Use your Agilent products only in the manner described in the Agilent product user guides.

---

**CAUTION**

**Safety standards for external equipment**

→ If you connect external equipment to the instrument, make sure that you only use accessory units tested and approved according to the safety standards appropriate for the type of external equipment.

---

## Overview of Maintenance

The following pages describe maintenance (simple repairs) of the detector that can be carried out without opening the main cover.

**Table 16** Simple Repairs

Procedures	Typical Frequency	Notes
Deuterium lamp exchange	If noise and/or drift exceeds your application limits or lamp does not ignite.	A VWD test should be performed after replacement.
Flow cell exchange	If application requires a different flow cell type.	A VWD test should be performed after replacement.
Cleaning flow cell parts cleaning or exchange	If leaking or if intensity drops due to contaminated flow cell windows.	A pressure tightness test should be done after repair.
Leak sensor drying	If leak has occurred.	Check for leaks.
Leak handling system replacement	If broken or corroded.	Check for leaks.

## Cleaning the Module

To keep the module case clean, use a soft cloth slightly dampened with water, or a solution of water and mild detergent.

### **WARNING**

**Liquid dripping into the electronic compartment of your module can cause shock hazard and damage the module**

- Do not use an excessively damp cloth during cleaning.
  - Drain all solvent lines before opening any connections in the flow path.
-

## Exchanging a Lamp

**When** If noise or drift exceeds application limits or lamp does not ignite.

**Tools required** **Description**  
Screwdriver, Pozidriv #1 PT3

Parts required	#	p/n	Description
	1	G1314-60101	Deuterium lamp (with RFID tag)

**Preparations** Turn the lamp OFF.

### NOTE

If you want to use an Agilent DAD lamp instead of the VWD lamp, you have to change the lamp settings in the *VWD Configuration* to the required lamp type. This ensures that the DAD lamp's filament heating is operated like in the DAD.

### NOTE

The specification are based on the the standard RFID tag lamp (G1314-60101) and may be not achieved when other lamp types or aged lamps are used.

### WARNING

#### Injury by touching hot lamp

**If the detector has been in use, the lamp may be hot.**

→ If so, wait for lamp to cool down.

### WARNING

#### Injury by sharp metal edges

→ Be careful when touching the RFI sheet metal at the rear of the fan. There are sharp edges.

### CAUTION


Electronic boards and components are sensitive to electrostatic discharge (ESD).

→ To prevent accidental electrostatic discharge when coming into contact with components inside the instrument, touch one of the metal housing panels at the front of the instrument.


## 9 Maintenance and Repair

### Exchanging a Lamp


- 1** Press the release buttons and remove the front cover to have access to the front area.

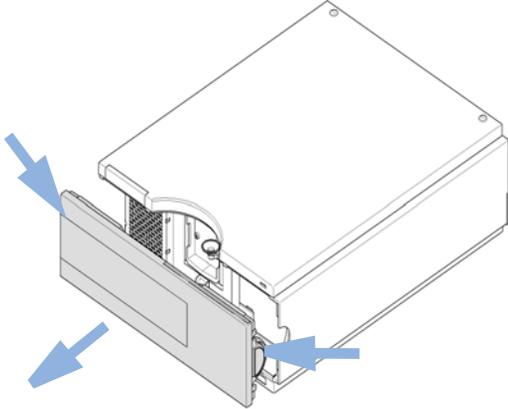


- 2** Unscrew the heater assembly and remove it.

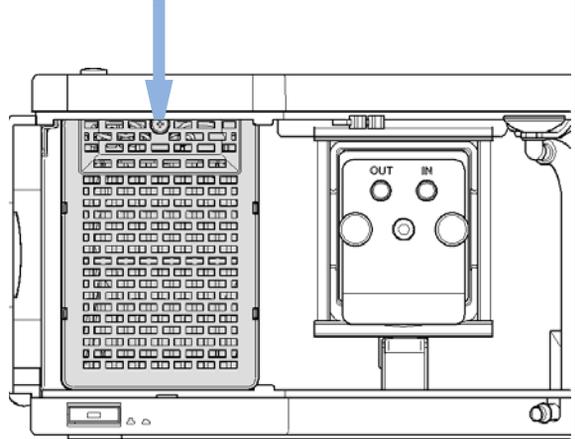

- 3** Unscrew, disconnect and remove the lamp. Insert, fix and reconnect the lamp.

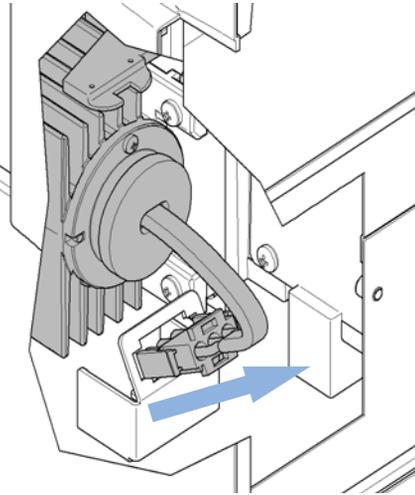

- 4** Replace the heater assembly.

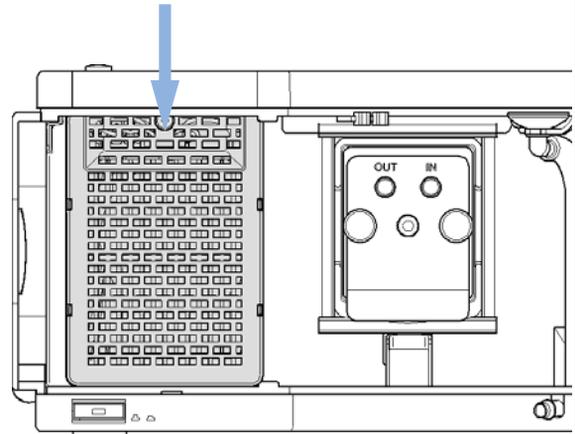

**Next Steps:**

- 5 Replace the front cover.
- 6 Reset the lamp counter as described in the User Interface documentation (required for non-RFID tag lamps only).
- 7 Turn the lamp ON.
- 8 Give the lamp more than 10 minutes to warm-up.
- 9 Perform “[Wavelength Verification-Calibration](#)” on page 120 to check the correct positioning of the lamp.

**NOTE**

If the detector was turned off during the replacement, then the detector requires a warm-up time of 60 minutes. No measurements should be performed during this time.

## Exchanging a Flow Cell

**When** If application needs a different type of flow cell or the flow cell needs repair.

**Tools required** **Description**  
Wrench, 1/4 inch  
for capillary connections

**Parts required**


#	Description
1	Flow cell

For flow cell details see


- “Standard Flow Cell 10 mm / 14  $\mu$ L” on page 152
- “Micro Flow Cell 3 mm / 2  $\mu$ L” on page 154
- “Semi-micro Flow Cell 6 mm / 5  $\mu$ L” on page 156
- “High Pressure Flow Cell 10 mm / 14  $\mu$ L” on page 158

**Preparations** Turn the lamp OFF.


**1** Press the release buttons and remove the front cover to have access to the flow cell area.

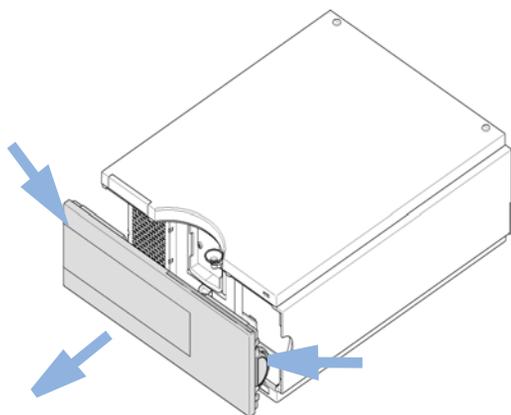

**2** Disconnect the inlet and outlet capillaries.

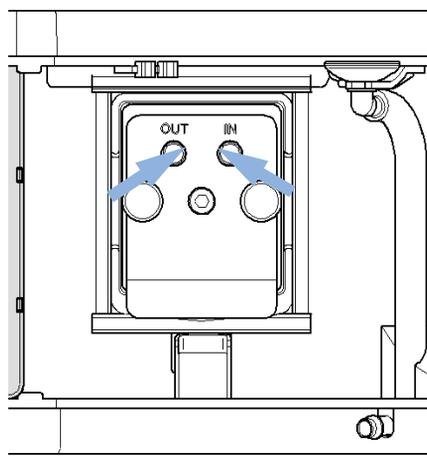


- 3** Unscrew both thumb screws parallel and remove the flow cell.

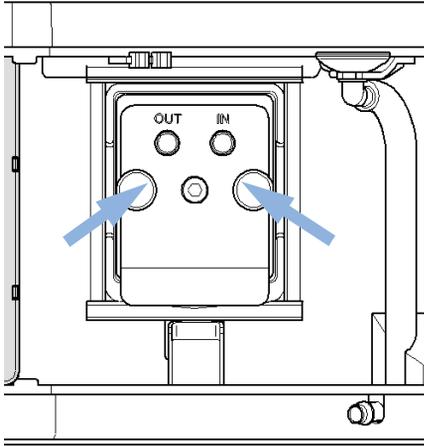


**NOTE**

If you want to maintain flow cell parts, see [“Overview of Maintenance Parts”](#) on page 150 or the information provided with your flow cell.

- 4** Replace the flow cell and fix the thumb screws. Reconnect the inlet and outlet capillaries to the flow cell.

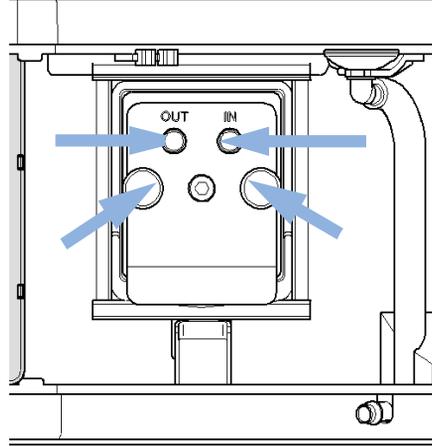

- 5** Replace the front cover.

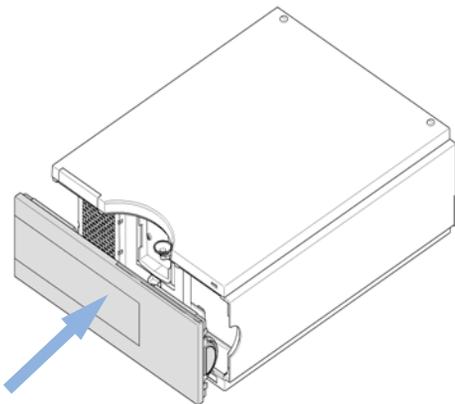

**Next Steps:**

- 6** To check for leaks, establish a flow and observe the flow cell (outside of the cell compartment) and all capillary connections.
- 7** Insert the flow cell.
- 8** Perform [“Wavelength Verification-Calibration”](#) on page 120 to check the correct positioning of the flow cell.
- 9** Replace the front cover.

## Repairing the Flow Cells

Parts required	#	Description
	1	Flow cell


For details on flow cells see

- “Standard Flow Cell 10 mm / 14  $\mu$ L” on page 152
- “Micro Flow Cell 3 mm / 2  $\mu$ L” on page 154
- “Semi-micro Flow Cell 6 mm / 5  $\mu$ L” on page 156
- “High Pressure Flow Cell 10 mm / 14  $\mu$ L” on page 158

### NOTE

The shown cell parts will differ depending upon the flow cell type. For detailed parts schematics, refer to above mentioned pages.

- 1 - Cell screw
- 2 - Conical springs
- 3 - Ring #1 PEEK
- 4 - Gasket #1 (small hole)
- 5 - Window Quartz
- 6 - Gasket #2 (large hole)
- 7 - Ring #2 PEEK
- 8 - RFID tag

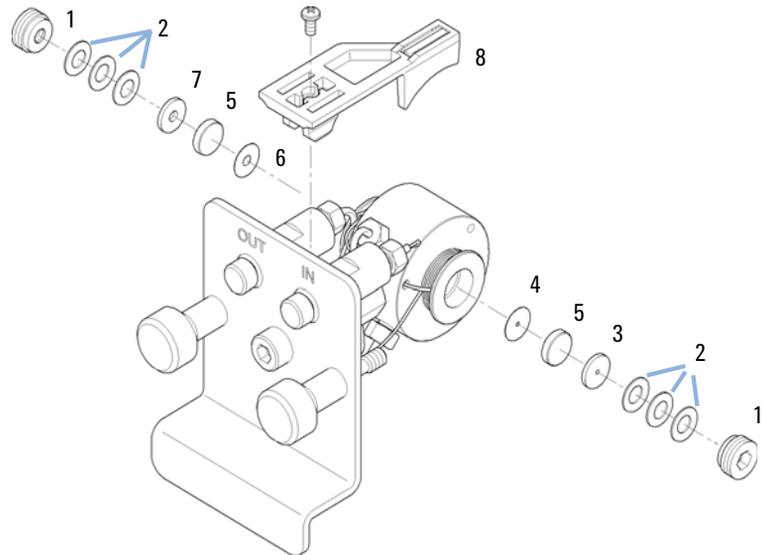

**Figure 46 Standard Flow Cell**

**1** Disassembling the Flow Cell.

- a** Unscrew the cell screw using a 4-mm hexagonal wrench.
- b** Remove the SST rings using a pair of tweezers.

**CAUTION**

Scatched window surfaces by tweezers

Window surfaces can easily be scratched by using tweezers for removing the windows.

→ Do not use tweezers to remove windows

- c** Use adhesive tape to remove the peek ring, the window and the gasket.
- d** Repeat step a through step c for the other window (keep the parts separate - otherwise they could be mixed!).

## 9 Maintenance and Repair

### Repairing the Flow Cells

#### 2 Cleaning the Flow Cell Parts

- a Pour isopropanol into the cell hole and wipe clean with a piece of lint-free cloth.
- b Clean the windows with ethanol or methanol. Dry it with a piece of lint-free cloth.

#### NOTE

Always use new gaskets.


#### 3 Reassembling the Flow Cell

- a Hold the flow cell cassette horizontally and place gasket in position. Ensure both cell holes can be seen through the holes of gasket.

#### NOTE

The semi-micro #1 and #2 gaskets (items 6 and 7, “Semi-micro Flow Cell 6 mm / 5  $\mu$ L” on page 156) look very similar. Do not mix them up.

- b Place the window on gasket.
- c Place the peek ring on the window.
- d Insert the conical springs. Make sure the conical springs point towards the window. Otherwise tightening the cell screw might break the window.

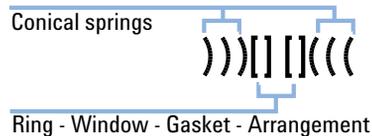


**Figure 47** Orientation of conical springs

- e Screw the cell screw into the flow cell and tighten the screw.
- 4 Repeat the procedure for the other cell side.
  - 5 Reconnect the capillaries.
  - 6 Perform a leak test. If OK, insert the flow cell.
  - 7 Perform “[Wavelength Verification-Calibration](#)” on page 120 to check the correct positioning of the flow cell.
  - 8 Replace the front cover.

## Using the Cuvette Holder

This cuvette holder can be placed instead of a flow cell in the variable wavelength detector. Standard cuvettes with standards in it, for example, National Institute of Standards & Technology (NIST) holmium oxide solution standard, can be fixed in it.

This can be used for wavelength verifications.

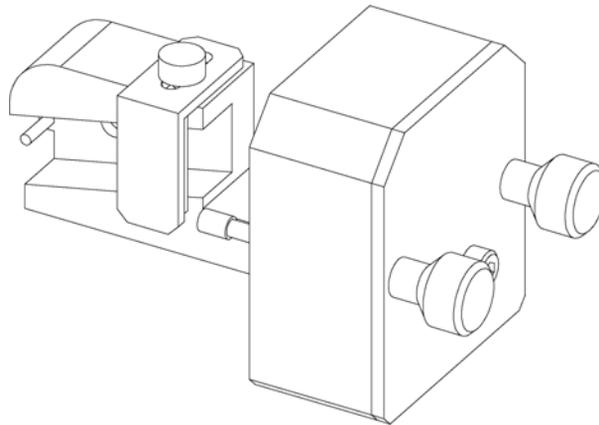

**When** If your own standard should be used to checkout the instrument.

Parts required	#	p/n	Description
	1	G1314-60200	Cuvette Holder
	1		Cuvette with the "standard", e.g. NIST certified holmium oxide sample


**Preparations**

- Remove the normal flow cell.
- Have cuvette with standard available.


## 9 Maintenance and Repair

### Using the Cuvette Holder


**1** Locate the cuvette holder on the desk.

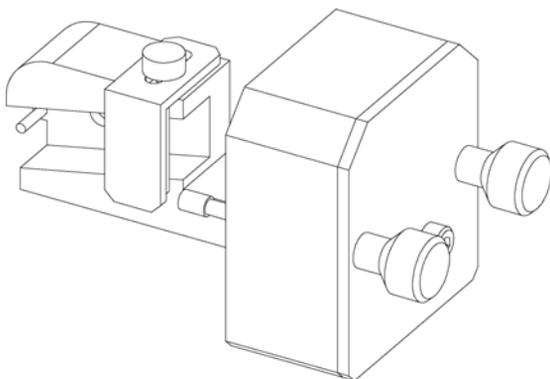


**2** Unscrew the bracket.

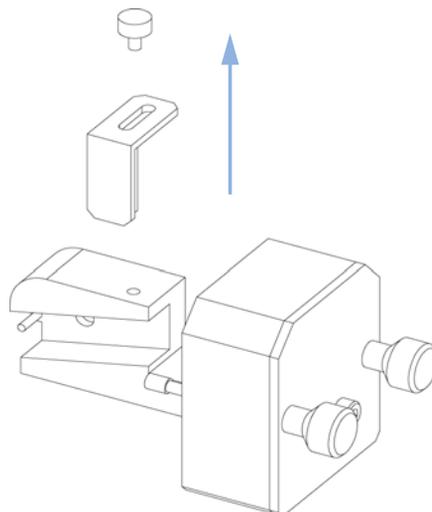

**3** Insert the cuvette with the sample into the holder. The clear side of the cuvette must be visible.

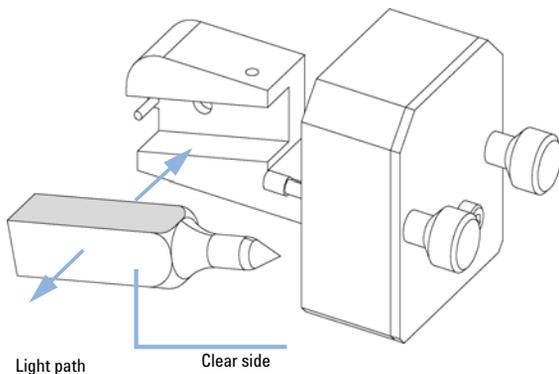

**4** Replace the bracket and fix the cuvette.

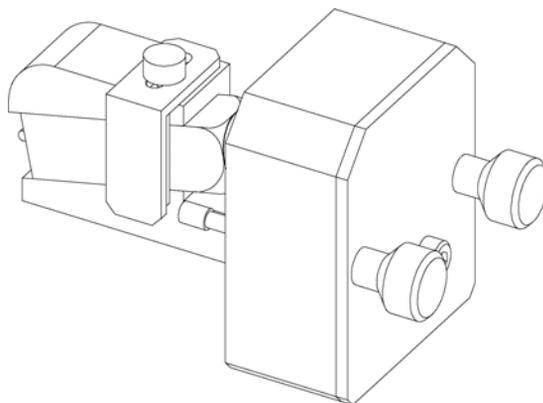

#### Next Steps:

- 5** Install the cuvette holder in the instrument.
- 6** Perform your Wavelength Verification/Calibration "[Wavelength Verification-Calibration](#)" on page 120 to check the correct position of the cuvette holder.

## Correcting Leaks

**When** If a leakage has occurred in the flow cell area or at the capillary connections.


**Tools required**

**Description**

Tissue

Wrench, 1/4 inch  
for capillary connections

- 1 Remove the front cover.
- 2 Use tissue to dry the leak sensor area.
- 3 Observe the capillary connections and the flow cell area for leaks and correct, if required.
- 4 Replace the front cover.

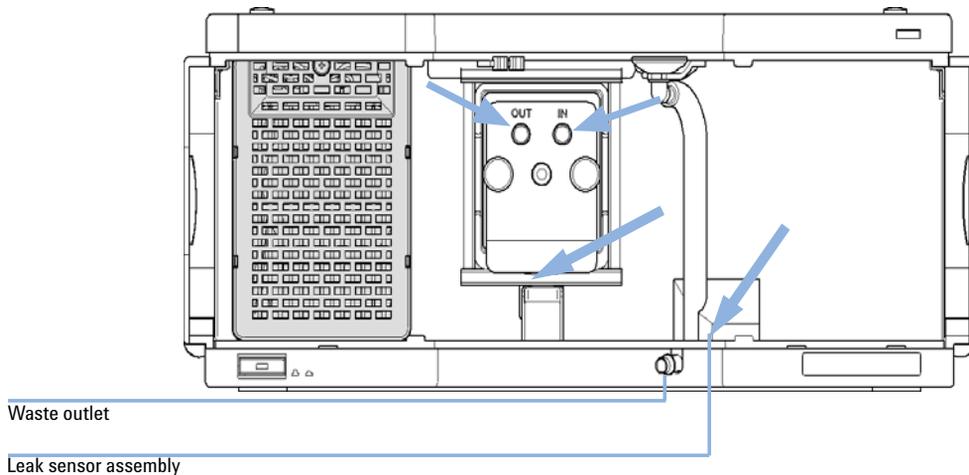

**Figure 48** Drying the Leak Sensor


## Replacing Leak Handling System Parts

**When** If the parts are corroded or broken.

**Tools required** None

<b>Parts required</b>	<b>#</b>	<b>p/n</b>	<b>Description</b>
	1	5041-8389	Leak funnel holder
	1	5041-8388	Leak funnel
	1	5062-2463	Corrugated tubing, PP, 6.5 mm id, 5 m

- 1 Remove the front cover to have access to the leak handling system.
- 2 Pull the leak funnel out of the leak funnel holder.
- 3 Pull the leak funnel with the tubing out of its location.
- 4 Replace the leak funnel and/or the tubing.
- 5 Insert the leak funnel with the tubing in its position.
- 6 Insert the leak funnel into the leak funnel holder.
- 7 Replace the front cover.

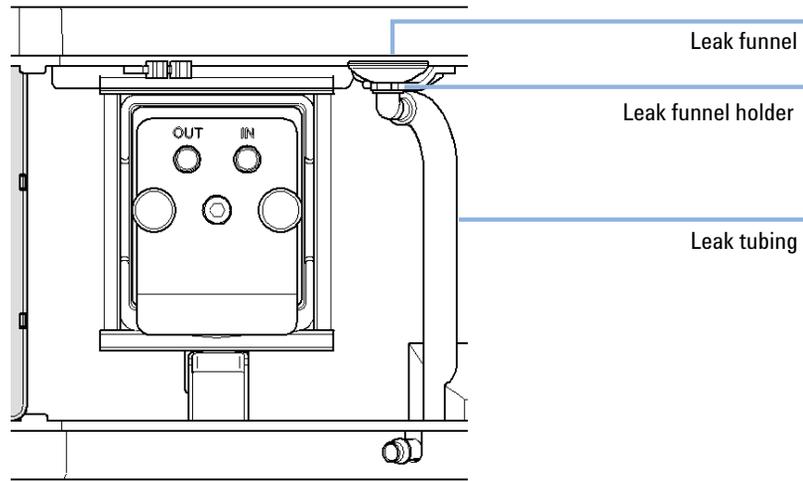

**Figure 49** Replacing Waste Handling System Parts

## Replacing the Module's Firmware

<b>When</b>	<p>The installation of newer firmware might be necessary</p> <ul style="list-style-type: none"> <li>• if a newer version solves problems of older versions or</li> <li>• to keep all systems on the same (validated) revision.</li> </ul> <p>The installation of older firmware might be necessary</p> <ul style="list-style-type: none"> <li>• to keep all systems on the same (validated) revision or</li> <li>• if a new module with newer firmware is added to a system or</li> <li>• if third party control software requires a special version.</li> </ul>
-------------	--

<b>Tools required</b>	<b>Description</b>
	LAN/RS-232 Firmware Update Tool
OR	Agilent Lab Advisor software
OR	Instant Pilot G4208A (only if supported by module)

<b>Parts required</b>	<b>#</b>	<b>Description</b>
	1	Firmware, tools and documentation from Agilent web site

**Preparations** Read update documentation provided with the Firmware Update Tool.

**To upgrade/downgrade the module's firmware carry out the following steps:**

- 1 Download the required module firmware, the latest LAN/RS-232 FW Update Tool and the documentation from the Agilent web.  
  - [http://www.chem.agilent.com/\\_layouts/agilent/downloadFirmware.aspx?whid=69761](http://www.chem.agilent.com/_layouts/agilent/downloadFirmware.aspx?whid=69761)
- 2 For loading the firmware into the module follow the instructions in the documentation.


## 9 Maintenance and Repair

### Replacing the Module's Firmware

#### *Module Specific Information*

**Table 17** Module Specific Information

	<b>G1314D</b>	<b>G1314E</b>	<b>G1314F</b>
Initial firmware	B.06.20	B.06.20	B.06.30
Compatibility with 1100 / 1200 series modules	When using the G1314D in a system, all other modules must have firmware revision A.06.10 or B.06.10 or above (main and resident). Otherwise the communication will not work.	When using the G1314E in a system, all other modules must have firmware revision A.06.10 or B.06.10 or above (main and resident). Otherwise the communication will not work.	When using the G1314F in a system, all other modules must have firmware revision A.06.30 or B.06.30 or above (main and resident). Otherwise the communication will not work.
Conversion to / emulation of G1314B or G1314C	Not possible due to different hardware and electronic platform		

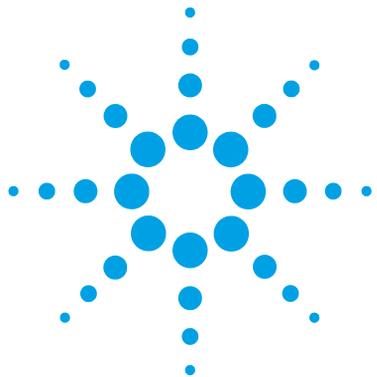

## 10 Parts and Materials for Maintenance

Overview of Maintenance Parts	150
Standard Flow Cell 10 mm / 14 $\mu$ L	152
Micro Flow Cell 3 mm / 2 $\mu$ L	154
Semi-micro Flow Cell 6 mm / 5 $\mu$ L	156
High Pressure Flow Cell 10 mm / 14 $\mu$ L	158
Cuvette Holder	160
Kits	151
Leak Parts	161

This chapter provides information on parts for maintenance.



## Overview of Maintenance Parts

<b>p/n</b>	<b>Description</b>
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m
G1314-60101	Deuterium lamp (with RFID tag)
G1314-60186	Standard flow cell 10 mm, 14 $\mu$ L, 40 bar (with RFID tag)
G1314-60187	Micro flow cell 3 mm, 2 $\mu$ L, 120 bar (with RFID tag)
G1314-60183	Semi-micro flow cell 6 mm, 5 $\mu$ L (with RFID tag)
G1314-60182	High pressure flow cell 10 mm, 14 $\mu$ L, 400 bar (with RFID tag)
G1314-60200	Cuvette Holder
5067-4691	Front Panel DAD/VWD/FLD (1260/1290)
5065-9982	Front Panel DAD/VWD/FLD (1200)

For details on flow cells, refer to

- [“Standard Flow Cell 10 mm / 14  \$\mu\$ L”](#) on page 152,
- [“Micro Flow Cell 3 mm / 2  \$\mu\$ L”](#) on page 154,
- [“Semi-micro Flow Cell 6 mm / 5  \$\mu\$ L”](#) on page 156 and
- [“High Pressure Flow Cell 10 mm / 14  \$\mu\$ L”](#) on page 158.

## Kits

### HPLC System Tool Kit

HPLC System Tool Kit (G4203-68708) contains some accessories and tools needed for installation and maintenance of the module.

### Accessory Kit

Accessory kit (G1314-68755) contains some accessories and tools needed for installation and repair of the module.

<b>p/n</b>	<b>Description</b>
0100-1516	Fitting male PEEK, 2/pk
5062-8535	Waste accessory kit, PEEK capillary 0.25 mm i.d., 1/16 o.d., 500 mm long plus 2 MT PTFE tubing i.d. 0.8 mm, 1/16 o.d.
5063-6527	Tubing assembly, i.d. 6 mm, o.d. 9 mm, 1.2 m (to waste)
5181-1516	CAN cable, Agilent module to module, 0.5 m

## 10 Parts and Materials for Maintenance

### Standard Flow Cell 10 mm / 14 $\mu$ L

## Standard Flow Cell 10 mm / 14 $\mu$ L

Item	p/n	Description
	G1314-60186	Standard flow cell 10 mm, 14 $\mu$ L, 40 bar (with RFID tag)
	5062-8522	Capillary column - detector PEEK 600 mm lg, 0.17 mm i.d., 1/16 inch o.d.
	G1314-65061	Cell Repair Kit, includes 2x Gasket #1, 2x Gasket #2, 2x Window Quartz
1	G1314-65062	Cell screw kit
2	79853-29100	Conical spring kit, 10/pk
3	G1314-65066	Ring #2 kit (IN small hole, i.d. 1 mm) PEEK, 2/pk
4	G1314-65064	Gaskets #2 IN (small hole i.d. 1 mm), KAPTON 10/pk
5	79853-68742	Window quartz kit, 2/pk
6	G1314-65063	Gasket #1 kit (OUT large hole, i.d. 2.4 mm) KAPTON, 2/pk
7	G1314-65065	Ring #1 kit (OUT large hole, i.d. 2.4 mm) PEEK, 2/pk
8	G1314-44010	Clip for RFI ID tag
9	0515-4780	Screw for Clip, M2.2, 4.5 mm long

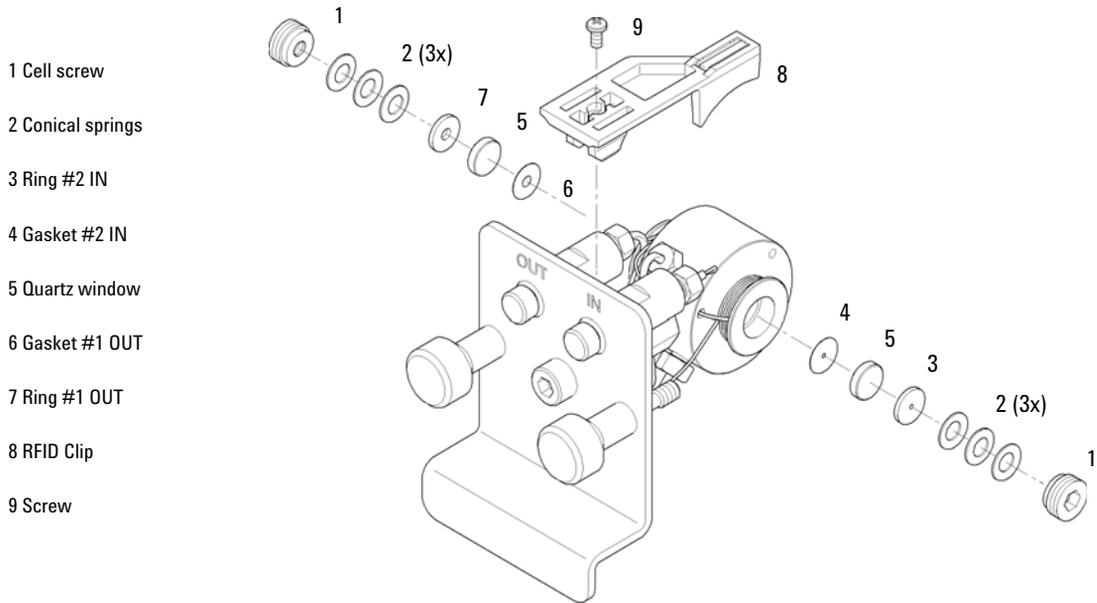


Figure 50 Standard Flow Cell

## 10 Parts and Materials for Maintenance

### Micro Flow Cell 3 mm / 2 $\mu$ L

## Micro Flow Cell 3 mm / 2 $\mu$ L

Item	p/n	Description
	G1314-60187	Micro flow cell 3 mm, 2 $\mu$ L, 120 bar (with RFID tag)
	5021-1823	Capillary column – detector SST 400 mm lg, 0.12 mm i.d.
1	79883-22402	Window screw
2	5062-8553	Washer kit (10/pk)
3	79883-28801	Compression washer
4	79883-22301	Window holder
5	1000-0488	Quartz window
6	G1315-68710	Gasket FRONT (PTFE), 1.3 mm hole, inlet side (12/pk)
7	79883-68702	Gasket BACK (PTFE), 1.8 mm hole, outlet side (12/pk)
8	G1314-44010	Clip for RFI ID tag
9	0515-4780	Screw for Clip, M2.2, 4.5 mm long
	G1314-87301	Capillary IN (0.12 mm, 310 mm lg)
	G1314-87302	Capillary OUT (0.17 mm, 120 mm lg)
	G1315-68713	Cell repair kit semi-micro, includes window screw kit, Gasket Kit BACK, Gasket Kit FRONT and 4 mm hexagonal wrench
	79883-68703	Window screw kit, includes 2 quartz windows, 2 compression washers, 2 window holders, 2 window screws and 10 washers

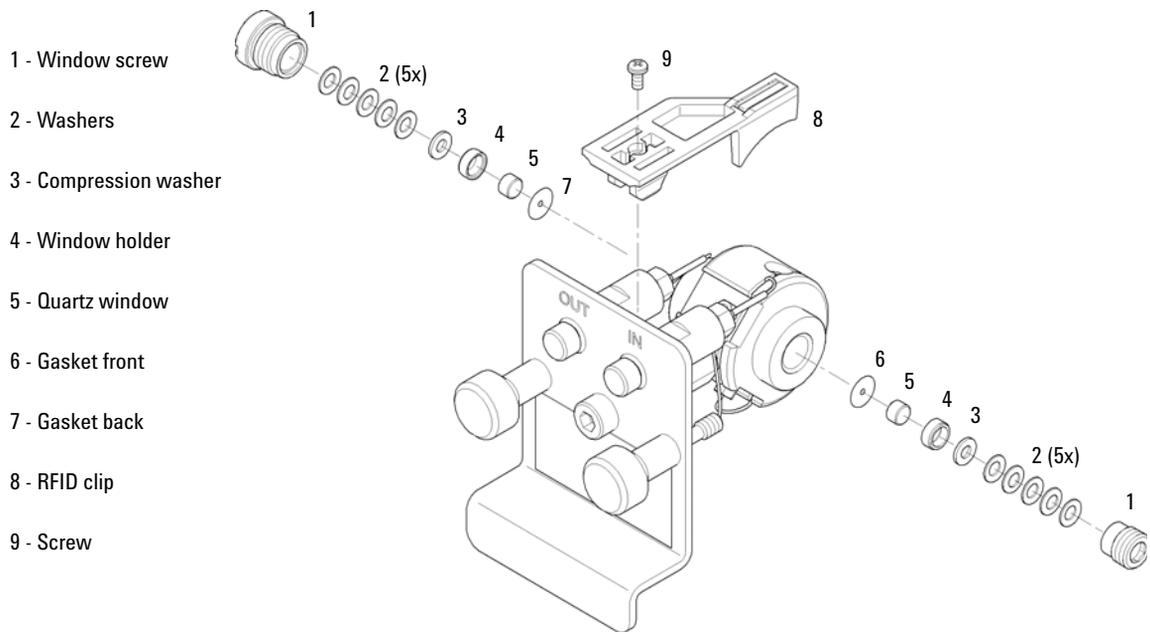


Figure 51 Micro Flow Cell

## 10 Parts and Materials for Maintenance


### Semi-micro Flow Cell 6 mm / 5 $\mu$ L

# Semi-micro Flow Cell 6 mm / 5 $\mu$ L

#### NOTE

The semi-micro #1 and #2 gaskets (items 6 and 7) look very similar. Do not mix them up.

Item	p/n	Description
	G1314-60183	Semi-micro flow cell 6 mm, 5 $\mu$ L (with RFID tag)
	5021-1823	Capillary column – detector SST 400 mm lg, 0.12 mm i.d.
1	G1314-20047	Cell screw
	G1314-65056	Semi-micro cell kit, includes two quartz windows, one gasket #1, one #2 and two PTFE gaskets.
2	79853-29100	Conical spring kit, 10/pk
3	79853-22500	Ring SST, 2/pk
4	79853-68743	PTFE gasket (round hole i.d. 2.5 mm, o.d. 8 mm), (10/pk)
5	79853-68742	Window quartz kit, 2/pk
6		Semi-micro #1 gasket (long hole 1.5 x 3.5 mm), PTFE
7		Semi-micro #2 gasket (long hole 2 x 4 mm), PTFE
8	G1314-44010	Clip for RFI ID tag
9	0515-4780	Screw for Clip, M2.2, 4.5 mm long

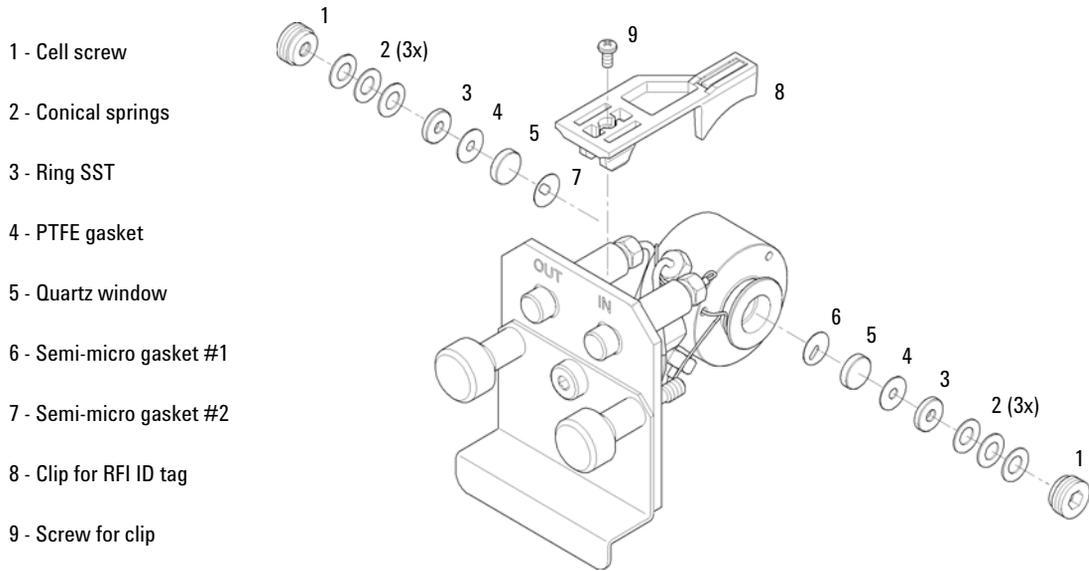


**Figure 52** Semi-micro Flow Cell

## 10 Parts and Materials for Maintenance

### High Pressure Flow Cell 10 mm / 14 $\mu$ L

## High Pressure Flow Cell 10 mm / 14 $\mu$ L

Item	p/n	Description
	G1314-60182	High pressure flow cell 10 mm, 14 $\mu$ L, 400 bar (with RFID tag)
	G1315-87311	Capillary ST 0.17 mm x 380 mm S/S
1	G1314-20047	Cell screw
	G1314-65054	Cell kit Agilent, comprises: two windows, two KAPTON gaskets and two PEEK rings
2		Ring PEEK kit
3		Window quartz kit
4		Gasket kit, KAPTON
5	G1314-44010	Clip for RFI ID tag
6	0515-4780	Screw for Clip, M2.2, 4.5 mm long

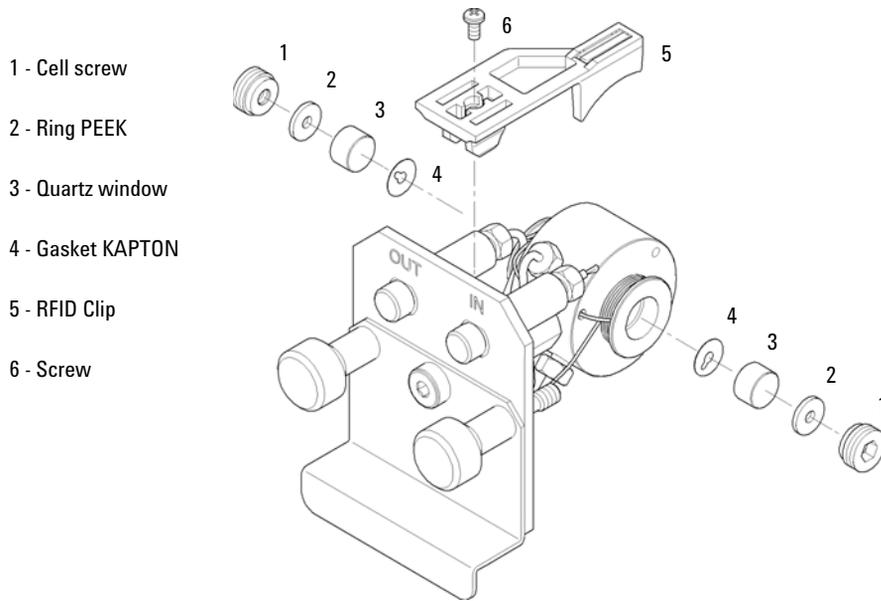


**Figure 53** High Pressure Flow Cell

## Cuvette Holder

For information the use of the cuvette holder, refer to “Using the Cuvette Holder” on page 143.


<b>p/n</b>	<b>Description</b>
G1314-60200	Cuvette Holder

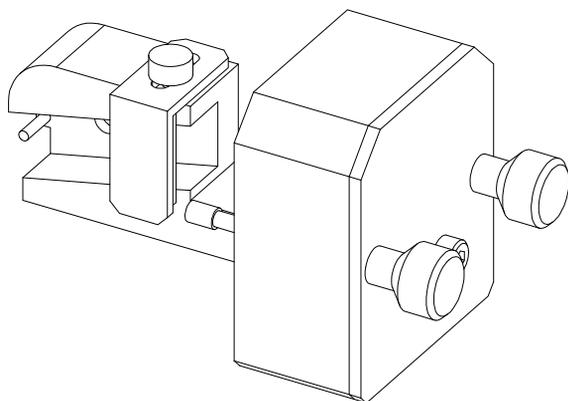

**Figure 54** Cuvette Holder

## Leak Parts


Item	p/n	Description
3	5041-8388	Leak funnel
4	5041-8389	Leak funnel holder
5	5041-8387	Tube clip
6	5062-2463	Corrugated tubing, PP, 6.5 mm id, 5 m
7	5062-2463	Corrugated tubing, PP, 6.5 mm id, 5 m

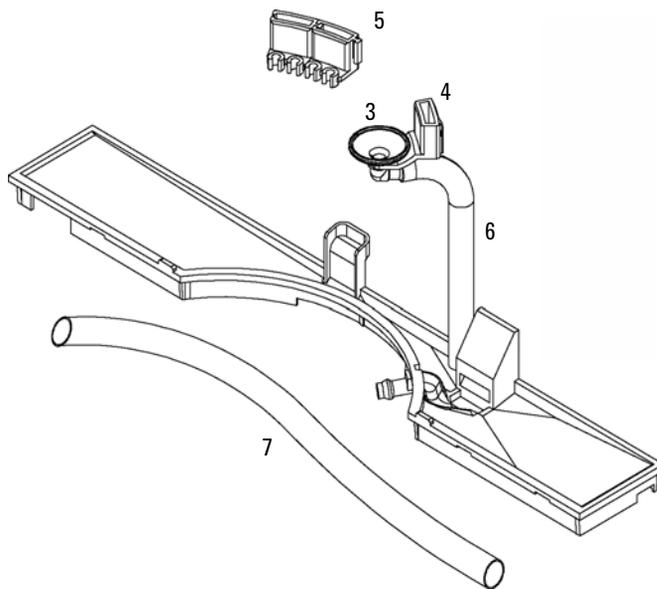

**Figure 55** Leak Parts

## **10 Parts and Materials for Maintenance**

### **Leak Parts**

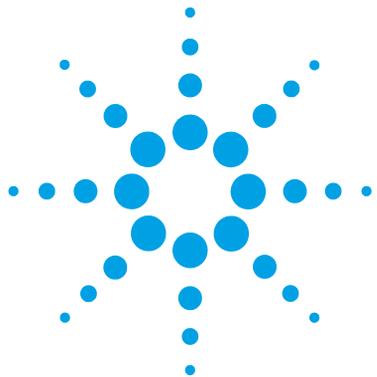

## 11 Identifying Cables

Cable Overview	164
Analog Cables	166
Remote Cables	168
BCD Cables	171
CAN/LAN Cables	173
RS-232 Cables	174

This chapter provides information on cables used with the Agilent 1200 Infinity Series modules.



# Cable Overview

**NOTE**

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

### Analog cables

p/n	Description
35900-60750	Agilent module to 3394/6 integrators
35900-60750	Agilent 35900A A/D converter
01046-60105	Analog cable (BNC to general purpose, spade lugs)

### Remote cables

p/n	Description
03394-60600	Agilent module to 3396A Series I integrators 3396 Series II / 3395A integrator, see details in section " <a href="#">Remote Cables</a> " on page 168
03396-61010	Agilent module to 3396 Series III / 3395B integrators
5061-3378	Remote Cable
01046-60201	Agilent module to general purpose

### BCD cables

p/n	Description
03396-60560	Agilent module to 3396 integrators
G1351-81600	Agilent module to general purpose

### CAN cables

p/n	Description
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m


### LAN cables

p/n	Description
5023-0203	Cross-over network cable, shielded, 3 m (for point to point connection)
5023-0202	Twisted pair network cable, shielded, 7 m (for point to point connection)

### RS-232 cables


p/n	Description
G1530-60600	RS-232 cable, 2 m
RS232-61601	RS-232 cable, 2.5 m Instrument to PC, 9-to-9 pin (female). This cable has special pin-out, and is not compatible with connecting printers and plotters. It's also called "Null Modem Cable" with full handshaking where the wiring is made between pins 1-1, 2-3, 3-2, 4-6, 5-5, 6-4, 7-8, 8-7, 9-9.
5181-1561	RS-232 cable, 8 m

## Analog Cables

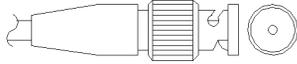


One end of these cables provides a BNC connector to be connected to Agilent modules. The other end depends on the instrument to which connection is being made.


### Agilent Module to 3394/6 Integrators

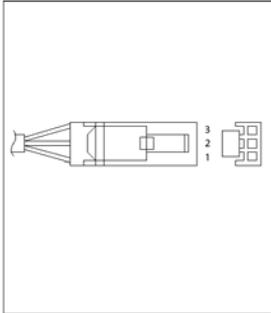p/n 35900-60750	Pin 3394/6	Pin Agilent module	Signal Name
	1		Not connected
	2	Shield	Analog -
	3	Center	Analog +

### Agilent Module to BNC Connector

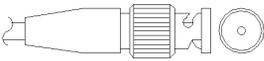p/n 8120-1840	Pin BNC	Pin Agilent module	Signal Name
	Shield	Shield	Analog -
	Center	Center	Analog +

### Agilent Module to General Purpose

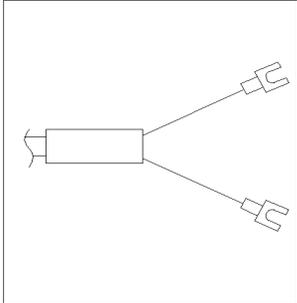p/n 01046-60105	Pin	Pin Agilent module	Signal Name
	1		Not connected
	2	Black	Analog -
	3	Red	Analog +

## Remote Cables



One end of these cables provides a Agilent Technologies APG (Analytical Products Group) remote connector to be connected to Agilent modules. The other end depends on the instrument to be connected to.


### Agilent Module to 3396A Integrators

p/n 03394-60600	Pin 3396A	Pin Agilent module	Signal Name	Active (TTL)
<p>A diagram of the Agilent module connector. It shows a vertical connector with pins labeled 1, 3, 5, 7, 9, 13, and 15. Pin 1 is at the bottom, and pin 15 is at the top. There are also pins 8 and 9 labeled on the right side of the connector.</p>	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	5,14	7 - Red	Ready	High
	1	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	


### Agilent Module to 3396 Series II / 3395A Integrators

Use the cable Agilent module to 3396A Series I integrators (03394-60600) and cut pin #5 on the integrator side. Otherwise the integrator prints START; not ready.

### Agilent Module to 3396 Series III / 3395B Integrators

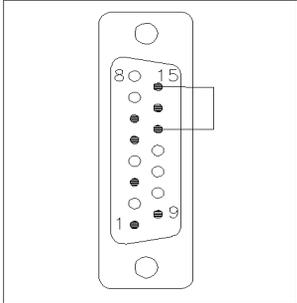p/n 03396-61010	Pin 33XX	Pin Agilent module	Signal Name	Active (TTL)
	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	14	7 - Red	Ready	High
	4	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	


### Agilent Module to Agilent 35900 A/D Converters

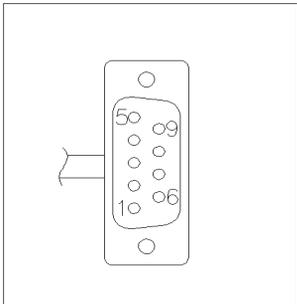p/n 5061-3378	Pin 35900 A/D	Pin Agilent module	Signal Name	Active (TTL)
	1 - White	1 - White	Digital ground	
	2 - Brown	2 - Brown	Prepare run	Low
	3 - Gray	3 - Gray	Start	Low
	4 - Blue	4 - Blue	Shut down	Low
	5 - Pink	5 - Pink	Not connected	
	6 - Yellow	6 - Yellow	Power on	High
	7 - Red	7 - Red	Ready	High
	8 - Green	8 - Green	Stop	Low
	9 - Black	9 - Black	Start request	Low

## 11 Identifying Cables

### Remote Cables

#### Agilent Module to General Purpose

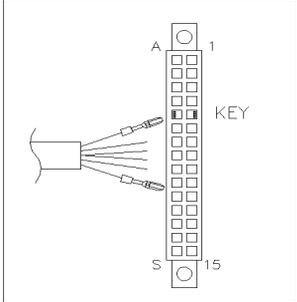
p/n 01046-60201	Wire Color	Pin Agilent module	Signal Name	Active (TTL)
	White	1	Digital ground	
	Brown	2	Prepare run	Low
	Gray	3	Start	Low
	Blue	4	Shut down	Low
	Pink	5	Not connected	
	Yellow	6	Power on	High
	Red	7	Ready	High
	Green	8	Stop	Low
	Black	9	Start request	Low

## BCD Cables



One end of these cables provides a 15-pin BCD connector to be connected to the Agilent modules. The other end depends on the instrument to be connected to


### Agilent Module to General Purpose

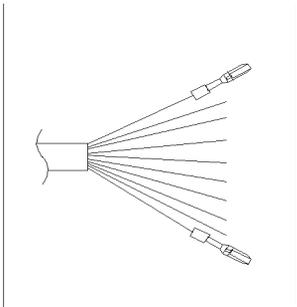p/n G1351-81600	Wire Color	Pin Agilent module	Signal Name	BCD Digit
	Green	1	BCD 5	20
	Violet	2	BCD 7	80
	Blue	3	BCD 6	40
	Yellow	4	BCD 4	10
	Black	5	BCD 0	1
	Orange	6	BCD 3	8
	Red	7	BCD 2	4
	Brown	8	BCD 1	2
	Gray	9	Digital ground	Gray
	Gray/pink	10	BCD 11	800
	Red/blue	11	BCD 10	400
	White/green	12	BCD 9	200
	Brown/green	13	BCD 8	100
	not connected	14		
	not connected	15	+ 5 V	Low


## 11 Identifying Cables

### BCD Cables

#### Agilent Module to 3396 Integrators

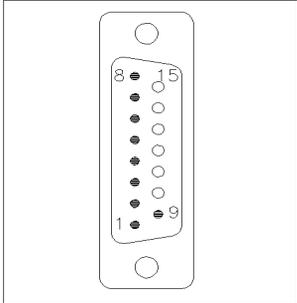p/n 03396-60560	Pin 3396	Pin Agilent module	Signal Name	BCD Digit
	1	1	BCD 5	20
	2	2	BCD 7	80
	3	3	BCD 6	40
	4	4	BCD 4	10
	5	5	BCD0	1
	6	6	BCD 3	8
	7	7	BCD 2	4
	8	8	BCD 1	2
	9	9	Digital ground	
	NC	15	+ 5 V	Low

## CAN/LAN Cables

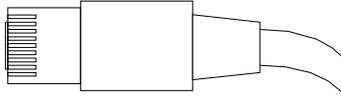

Both ends of this cable provide a modular plug to be connected to Agilent modules CAN or LAN connectors.

### CAN Cables


p/n	Description
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m

### LAN Cables

p/n	Description
5023-0203	Cross-over network cable, shielded, 3 m (for point to point connection)
5023-0202	Twisted pair network cable, shielded, 7 m (for point to point connection)

## RS-232 Cables

<b>p/n</b>	<b>Description</b>
G1530-60600	RS-232 cable, 2 m
RS232-61601	RS-232 cable, 2.5 m Instrument to PC, 9-to-9 pin (female). This cable has special pin-out, and is not compatible with connecting printers and plotters. It's also called "Null Modem Cable" with full handshaking where the wiring is made between pins 1-1, 2-3, 3-2, 4-6, 5-5, 6-4, 7-8, 8-7, 9-9.
5181-1561	RS-232 cable, 8 m

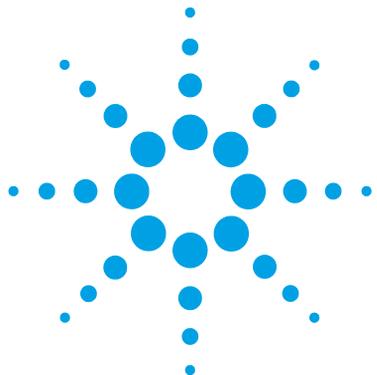

## 12 Hardware Information

Firmware Description	176
Electrical Connections	179
Rear View of the Module	180
Information on Instrument Serial Number	181
Interfaces	182
Overview Interfaces	185
Setting the 8-bit Configuration Switch	189
Special Settings	191
Instrument Layout	193
Early Maintenance Feedback (EMF)	194
EMF Counter	194
Using the EMF Counters	195

This chapter describes the detector in more detail on hardware and electronics.



## Firmware Description

The firmware of the instrument consists of two independent sections:

- a non-instrument specific section, called *resident system*
- an instrument specific section, called *main system*

### Resident System

This resident section of the firmware is identical for all Agilent 1100/1200/1220/1260/1290 series modules. Its properties are:

- the complete communication capabilities (CAN, LAN and RS-232C)
- memory management
- ability to update the firmware of the 'main system'

### Main System

Its properties are:

- the complete communication capabilities (CAN, LAN and RS-232C)
- memory management
- ability to update the firmware of the 'resident system'

In addition the main system comprises the instrument functions that are divided into common functions like

- run synchronization through APG remote,
- error handling,
- diagnostic functions,
- or module specific functions like
  - internal events such as lamp control, filter movements,
  - raw data collection and conversion to absorbance.

## Firmware Updates

Firmware updates can be done using your user interface:

- PC and Firmware Update Tool with local files on the hard disk
- Instant Pilot (G4208A) with files from a USB Flash Disk
- Agilent Lab Advisor software B.01.03 and above

The file naming conventions are:

PPPP\_RVVV\_XXX.dlb, where

PPPP is the product number, for example, 1315AB for the G1315A/B DAD,

R the firmware revision, for example, A for G1315B or B for the G1315C DAD,

VVV is the revision number, for example 102 is revision 1.02,


XXX is the build number of the firmware.

For instructions on firmware updates refer to section *Replacing Firmware* in chapter "Maintenance" or use the documentation provided with the *Firmware Update Tools*.

### NOTE

Update of main system can be done in the resident system only. Update of the resident system can be done in the main system only.

Main and resident firmware must be from the same set.

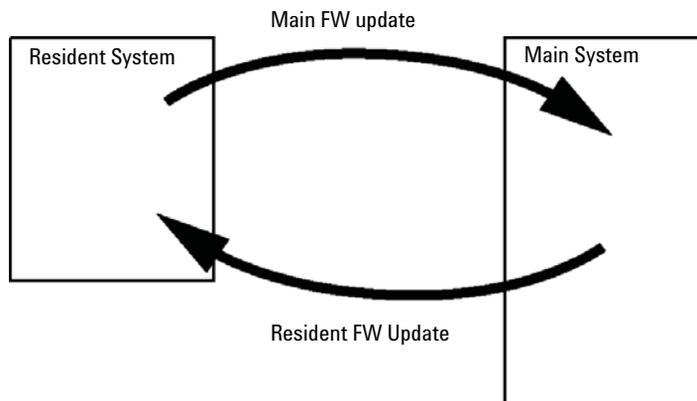

**Figure 56** Firmware Update Mechanism

## 12 Hardware Information

### Firmware Description

#### NOTE

Some modules are limited in downgrading due to their main board version or their initial firmware revision. For example, a G1315C DAD SL cannot be downgraded below firmware revision B.01.02 or to a A.xx.xx.

Some modules can be re-branded (e.g. G1314C to G1314B) to allow operation in specific control software environments. In this case the feature set of the target type are use and the feature set of the original are lost. After re-branding (e.g. from G1314B to G1314C), the original feature set is available again.

All these specific informations are described in the documentation provided with the firmware update tools.

---

The firmware update tools, firmware and documentation are available from the Agilent web.

- [http://www.chem.agilent.com/\\_layouts/agilent/downloadFirmware.aspx?whid=69761](http://www.chem.agilent.com/_layouts/agilent/downloadFirmware.aspx?whid=69761)

## Electrical Connections


- The CAN bus is a serial bus with high speed data transfer. The two connectors for the CAN bus are used for internal module data transfer and synchronization.
- One analog output provides signals for integrators or data handling systems.
- The REMOTE connector may be used in combination with other analytical instruments from Agilent Technologies if you want to use features such as start, stop, common shut down, prepare, and so on.
- With the appropriate software, the RS-232C connector may be used to control the module from a computer through a RS-232C connection. This connector is activated and can be configured with the configuration switch.
- The power input socket accepts a line voltage of 100 – 240 VAC  $\pm$  10 % with a line frequency of 50 or 60 Hz. Maximum power consumption varies by module. There is no voltage selector on your module because the power supply has wide-ranging capability. There are no externally accessible fuses, because automatic electronic fuses are implemented in the power supply.

### NOTE

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

---

## Rear View of the Module

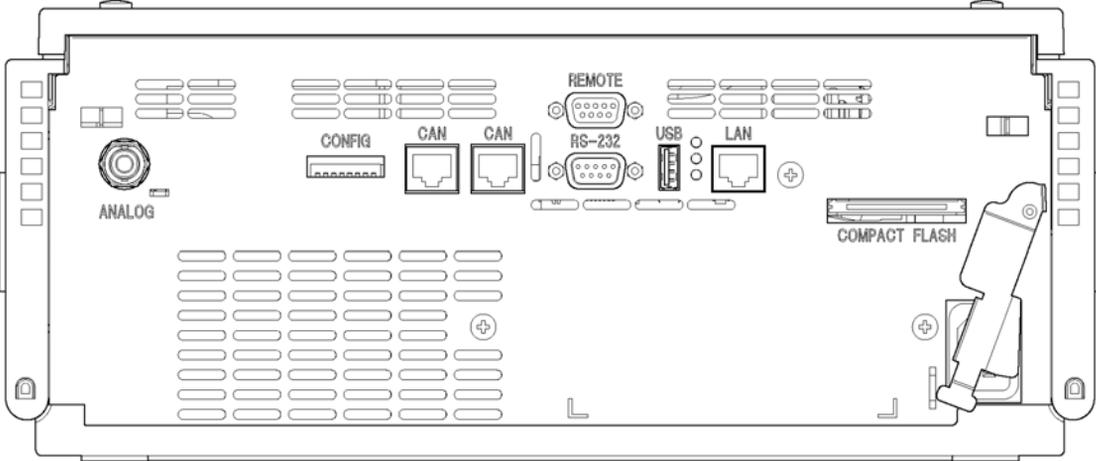

**Figure 57** Rear View of Detector

**NOTE**

The Compact Flash card slot is used for the G1314E VWD only.

## Information on Instrument Serial Number

### Serial Number Information 1200 Series and 1290 Infinity

The serial number information on the instrument labels provide the following information:

CCYWWSSSSS	Format
CC	country of manufacturing <ul style="list-style-type: none"> <li>• DE = Germany</li> <li>• JP = Japan</li> <li>• CN = China</li> </ul>
YWW	year and week of last major manufacturing change, e.g. 820 could be week 20 of 1998 or 2008
SSSSS	real serial number

### Serial Number Information 1260 Infinity

The serial number information on the instrument labels provide the following information:

CCXZZ00000	Format
CC	Country of manufacturing <ul style="list-style-type: none"> <li>• DE = Germany</li> <li>• JP = Japan</li> <li>• CN = China</li> </ul>
X	Alphabetic character A-Z (used by manufacturing)
ZZ	Alpha-numeric code 0-9, A-Z, where each combination unambiguously denotes a module (there can be more than one code for the same module)
00000	Serial number

## Interfaces

The Agilent 1200 Infinity Series modules provide the following interfaces:

**Table 18** Agilent 1200 Infinity Series Interfaces

Module	CAN	LAN/BCD (optional)	LAN (on-board)	RS-232	Analog	APG Remote	Special
<b>Pumps</b>							
G1310B Iso Pump G1311B Quat Pump G1311C Quat Pump VL G1312B Bin Pump K1312B Bin Pump Clinical Ed. G1312C Bin Pump VL 1376A Cap Pump G2226A Nano Pump G5611A Bio-inert Quat Pump	2	Yes	No	Yes	1	Yes	
G4220A/B Bin Pump G4204A Quat Pump	2	No	Yes	Yes	No	Yes	CAN-DC- OUT for CAN slaves
G1361A Prep Pump	2	Yes	No	Yes	No	Yes	CAN-DC- OUT for CAN slaves
<b>Samplers</b>							
G1329B ALS G2260A Prep ALS	2	Yes	No	Yes	No	Yes	THERMOSTAT for G1330B/K1330B
G1364B FC-PS G1364C FC-AS G1364D FC- $\mu$ S G1367E HiP ALS K1367E HiP ALS Clinical Ed. G1377A HiP micro ALS G2258A DL ALS G5664A Bio-inert FC-AS G5667A Bio-inert Autosampler	2	Yes	No	Yes	No	Yes	THERMOSTAT for G1330B/K1330B CAN-DC- OUT for CAN slaves
G4226A ALS	2	Yes	No	Yes	No	Yes	

**Table 18** Agilent 1200 Infinity Series Interfaces

Module	CAN	LAN/BCD (optional)	LAN (on-board)	RS-232	Analog	APG Remote	Special
<b>Detectors</b>							
G1314B VWD VL G1314C VWD VL+	2	Yes	No	Yes	1	Yes	
G1314E/F VWD K1314F Clinical Ed.	2	No	Yes	Yes	1	Yes	
G4212A/B DAD K4212B DAD Clinical Ed.	2	No	Yes	Yes	1	Yes	
G1315C DAD VL+ G1365C MWD G1315D DAD VL G1365D MWD VL	2	No	Yes	Yes	2	Yes	
G1321B FLD K1321B FLD Clinical Ed. G1321C FLD	2	Yes	No	Yes	2	Yes	
G1362A RID	2	Yes	No	Yes	1	Yes	
G4280A ELSD	No	No	No	Yes	Yes	Yes	EXT Contact AUTOZERO
<b>Others</b>							
G1170A Valve Drive	2	No	No	No	No	No	1
G1316A/C TCC K1316C TCC Clinical Ed.	2	No	No	Yes	No	Yes	
G1322A DEG K1322A DEG Clinical Ed.	No	No	No	No	No	Yes	AUX
G1379B DEG	No	No	No	Yes	No	Yes	
G4225A DEG K4225A DEG Clinical Ed.	No	No	No	Yes	No	Yes	

## 12 Hardware Information

### Interfaces

**Table 18** Agilent 1200 Infinity Series Interfaces

Module	CAN	LAN/BCD (optional)	LAN (on-board)	RS-232	Analog	APG Remote	Special
G4227A Flex Cube	2	No	No	No	No	No	CAN-DC- OUT for CAN slaves 1
G4240A CHIP CUBE	2	Yes	No	Yes	No	Yes	CAN-DC- OUT for CAN slaves THERMOSTAT for G1330A/B (NOT USED), K1330B

<sup>1</sup> Requires a HOST module with on-board LAN (e.g. G4212A or G4220A with minimum firmware B.06.40 or C.06.40) or with additional G1369C LAN Card

#### NOTE

The detector (DAD/MWD/FLD/VWD/RID) is the preferred access point for control via LAN. The inter-module communication is done via CAN.

- CAN connectors as interface to other modules
- LAN connector as interface to the control software
- RS-232C as interface to a computer
- REMOTE connector as interface to other Agilent products
- Analog output connector(s) for signal output

## Overview Interfaces

### CAN

The CAN is inter-module communication interface. It is a 2-wire serial bus system supporting high speed data communication and real-time requirement.

### LAN

The modules have either an interface slot for an LAN card (e.g. Agilent G1369B/C LAN Interface) or they have an on-board LAN interface (e.g. detectors G1315C/D DAD and G1365C/D MWD). This interface allows the control of the module/system via a PC with the appropriate control software. Some modules have neither on-board LAN nor an interface slot for a LAN card (e.g. G1170A Valve Drive or G4227A Flex Cube). These are hosted modules and require a Host module with firmware B.06.40 or later or with additional G1369C LAN Card.

#### NOTE

If an Agilent detector (DAD/MWD/FLD/VWD/RID) is in the system, the LAN should be connected to the DAD/MWD/FLD/VWD/RID (due to higher data load). If no Agilent detector is part of the system, the LAN interface should be installed in the pump or autosampler.

### RS-232C (Serial)

The RS-232C connector is used to control the module from a computer through RS-232C connection, using the appropriate software. This connector can be configured with the configuration switch module at the rear of the module. Refer to *Communication Settings for RS-232C*.

#### NOTE

There is no configuration possible on main boards with on-board LAN. These are pre-configured for

- 19200 baud,
- 8 data bit with no parity and
- one start bit and one stop bit are always used (not selectable).


The RS-232C is designed as DCE (data communication equipment) with a 9-pin male SUB-D type connector. The pins are defined as:

## 12 Hardware Information

### Interfaces

**Table 19** RS-232C Connection Table

Pin	Direction	Function
1	In	DCD
2	In	RxD
3	Out	TxD
4	Out	DTR
5		Ground
6	In	DSR
7	Out	RTS
8	In	CTS
9	In	RI

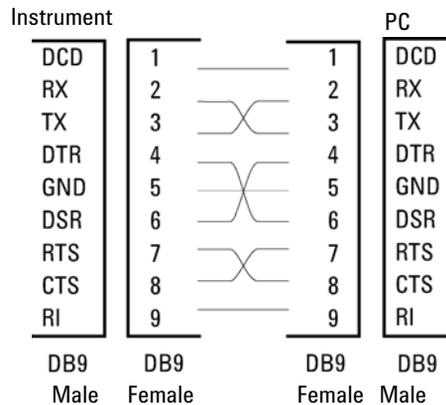

**Figure 58** RS-232 Cable

### Analog Signal Output

The analog signal output can be distributed to a recording device. For details refer to the description of the module's main board.

## APG Remote

The APG Remote connector may be used in combination with other analytical instruments from Agilent Technologies if you want to use features as common shut down, prepare, and so on.

Remote control allows easy connection between single instruments or systems to ensure coordinated analysis with simple coupling requirements.

The subminiature D connector is used. The module provides one remote connector which is inputs/outputs (wired- or technique).

To provide maximum safety within a distributed analysis system, one line is dedicated to **SHUT DOWN** the system's critical parts in case any module detects a serious problem. To detect whether all participating modules are switched on or properly powered, one line is defined to summarize the **POWER ON** state of all connected modules. Control of analysis is maintained by signal readiness **READY** for next analysis, followed by **START** of run and optional **STOP** of run triggered on the respective lines. In addition **PREPARE** and **START REQUEST** may be issued. The signal levels are defined as:

- standard TTL levels (0 V is logic true, + 5.0 V is false),
- fan-out is 10 ,
- input load is 2.2 kOhm against + 5.0 V, and
- output are open collector type, inputs/outputs (wired- or technique).

### NOTE

All common TTL circuits operate with a 5 V power supply. A TTL signal is defined as "low" or L when between 0 V and 0.8 V and "high" or H when between 2.0 V and 5.0 V (with respect to the ground terminal).

**Table 20** Remote Signal Distribution

Pin	Signal	Description
1	DGND	Digital ground
2	PREPARE	(L) Request to prepare for analysis (for example, calibration, detector lamp on). Receiver is any module performing pre-analysis activities.
3	START	(L) Request to start run / timetable. Receiver is any module performing run-time controlled activities.
4	SHUT DOWN	(L) System has serious problem (for example, leak: stops pump). Receiver is any module capable to reduce safety risk.
5		Not used
6	POWER ON	(H) All modules connected to system are switched on. Receiver is any module relying on operation of others.
7	READY	(H) System is ready for next analysis. Receiver is any sequence controller.
8	STOP	(L) Request to reach system ready state as soon as possible (for example, stop run, abort or finish and stop injection). Receiver is any module performing run-time controlled activities.
9	START REQUEST	(L) Request to start injection cycle (for example, by start key on any module). Receiver is the autosampler.

### Special Interfaces

There is no special interface for this module.

## Setting the 8-bit Configuration Switch


The 8-bit configuration switch is located at the rear of the module. Switch settings provide configuration parameters for LAN, serial communication protocol and instrument specific initialization procedures.

All modules with on-board LAN:

- Default is ALL switches DOWN (best settings).
  - Bootp mode for LAN and
  - 19200 baud, 8 data bit / 1 stop bit with no parity for RS-232
- For specific LAN modes switches 3-8 must be set as required.
- For boot/test modes switches 1+2 must be UP plus required mode.

### NOTE

For normal operation use the default (best) settings.

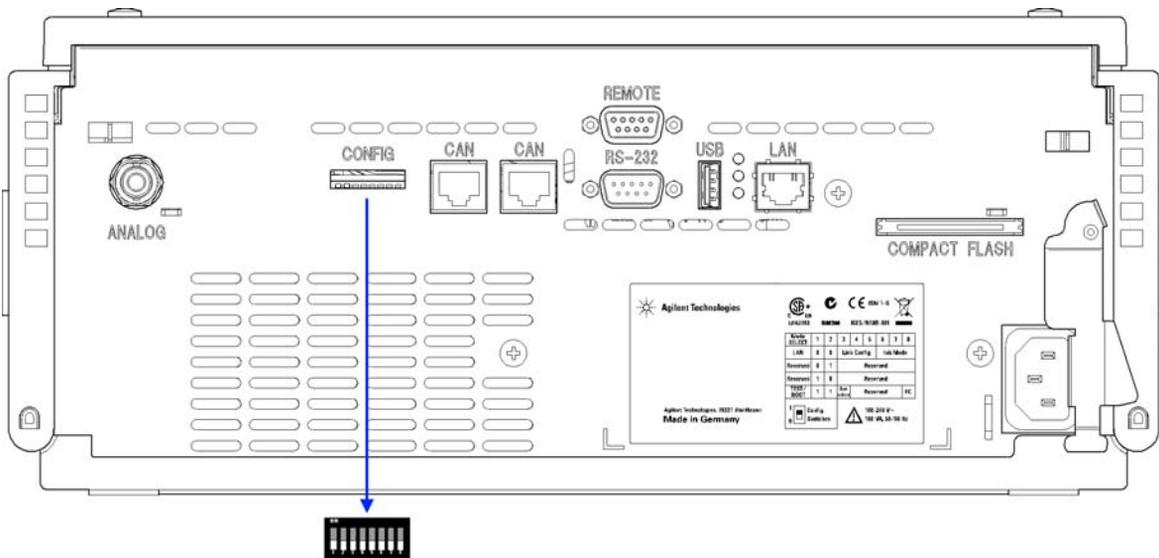

**Figure 59** Location of Configuration Switch (example shows a G4212A DAD)

## 12 Hardware Information

### Setting the 8-bit Configuration Switch

#### NOTE

To perform any LAN configuration, SW1 and SW2 must be set to OFF. For details on the LAN settings/configuration refer to chapter LAN Configuration.

**Table 21** 8-bit Configuration Switch (with on-board LAN)

	Mode		Function					
	SW 1	SW 2	SW 3	SW 4	SW 5	SW 6	SW 7	SW 8
LAN	0	0	Link Configuration			Init Mode Selection		
Auto-negotiation			0	x	x	x	x	x
10 MBit, half-duplex			1	0	0	x	x	x
10 MBit, full-duplex			1	0	1	x	x	x
100 MBit, half-duplex			1	1	0	x	x	x
100 MBit, full-duplex			1	1	1	x	x	x
Bootp			x	x	x	0	0	0
Bootp & Store			x	x	x	0	0	1
Using Stored			x	x	x	0	1	0
DHCP			x	x	x	1	0	0
Using Default			x	x	x	0	1	1
TEST	1	1	System					NVRAM
Boot Resident System			1					x
Revert to Default Data (Coldstart)			x	x	x			1

#### Legend:

0 (switch down), 1 (switch up), x (any position)

#### NOTE

When selecting the mode TEST, the LAN settings are: Auto-Negotiation & Using Stored.

#### NOTE

For explanation of "Boot Resident System" and "Revert to Default Data (Coldstart)" refer to "Special Settings" on page 191.

## Special Settings

The special settings are required for specific actions (normally in a service case).

**NOTE**

The tables include both settings for modules – with on-board LAN and without on-board LAN. They are identified as LAN and no LAN.

### Boot-Resident

Firmware update procedures may require this mode in case of firmware loading errors (main firmware part).

If you use the following switch settings and power the instrument up again, the instrument firmware stays in the resident mode. It is not operable as a module. It only uses basic functions of the operating system for example, for communication. In this mode the main firmware can be loaded (using update utilities).

**Table 22** Boot Resident Settings (On-board LAN)

Mode Select	SW1	SW2	SW3	SW4	SW5	SW6	SW7	SW8
TEST/BOOT	1	1	1	0	0	0	0	0

## 12 Hardware Information

### Setting the 8-bit Configuration Switch

#### Forced Cold Start

A forced cold start can be used to bring the module into a defined mode with default parameter settings.

#### CAUTION

Loss of data

Forced cold start erases all methods and data stored in the non-volatile memory. Exceptions are calibration settings, diagnosis and repair log books which will not be erased.

→ Save your methods and data before executing a forced cold start.

If you use the following switch settings and power the instrument up again, a forced cold start has been completed.

**Table 23** Forced Cold Start Settings (On-board LAN)

Mode Select	SW1	SW2	SW3	SW4	SW5	SW6	SW7	SW8
TEST/BOOT	1	1	0	0	0	0	0	1

## Instrument Layout

The industrial design of the module incorporates several innovative features. It uses Agilent's E-PAC concept for the packaging of electronics and mechanical assemblies. This concept is based upon the use of expanded polypropylene (EPP) layers of foam plastic spacers in which the mechanical and electronic boards components of the module are placed. This pack is then housed in a metal inner cabinet which is enclosed by a plastic external cabinet. The advantages of this packaging technology are:

- virtual elimination of fixing screws, bolts or ties, reducing the number of components and increasing the speed of assembly/disassembly,
- the plastic layers have air channels molded into them so that cooling air can be guided exactly to the required locations,
- the plastic layers help cushion the electronic and mechanical parts from physical shock, and
- the metal inner cabinet shields the internal electronics from electromagnetic interference and also helps to reduce or eliminate radio frequency emissions from the instrument itself.

## Early Maintenance Feedback (EMF)

Maintenance requires the exchange of components which are subject to wear or stress. Ideally, the frequency at which components are exchanged should be based on the intensity of usage of the instrument and the analytical conditions, and not on a predefined time interval. The early maintenance feedback (EMF) feature monitors the usage of specific components in the instrument, and provides feedback when the user-selectable limits have been exceeded. The visual feedback in the user interface provides an indication that maintenance procedures should be scheduled.

### EMF Counter

The detector module provides an EMF counter for the lamp. The counter increments with lamp use, and can be assigned a maximum limit which provides visual feedback in the user interface when the limit is exceeded. Dependig on the lamp type the counter can be reset to zero after the lamp is exchanged.

Lamp Type	Counter Reset	Comment
lamp with RFID tag	NO	
lamp without RFID tag	YES	via Lab Advisor or Instant Pilot

The detector provides the following EMF counters:

- Deuterium Lamp On-Time
- Number of UV lamp ignitions

## Using the EMF Counters

The user-settable EMF limits for the EMF counters enable the early maintenance feedback to be adapted to specific user requirements. The useful lamp burn time is dependent on the requirements for the analysis (high or low sensitivity analysis, wavelength etc.), therefore, the definition of the maximum limits need to be determined based on the specific operating conditions of the instrument.

### Setting the EMF Limits

The setting of the EMF limits must be optimized over one or two maintenance cycles. Initially, no EMF limit should be set. When instrument performance indicates maintenance is necessary, take note of the values displayed by lamp counters. Enter these values (or values slightly less than the displayed values) as EMF limits, and then reset the EMF counters to zero. The next time the EMF counters exceed the new EMF limits, the EMF flag will be displayed, providing a reminder that maintenance needs to be scheduled.


**NOTE**

This function is only available via Agilent Lab Advisor or Instant Pilot.

---

## **12 Hardware Information**

### **Early Maintenance Feedback (EMF)**

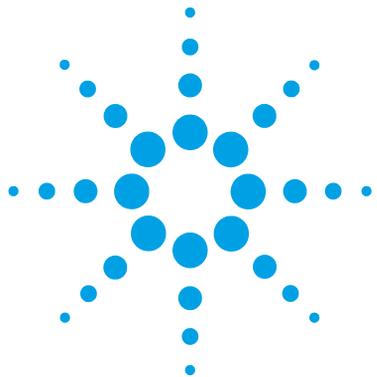

## 13 LAN Configuration

What you have to do first	198
TCP-IP parameter configuration	199
Configuration Switches	200
Initialization mode selection	201
Dynamic Host Configuration Protocol (DHCP)	205
General Information (DHCP)	205
Setup (DHCP)	206
Link configuration selection	208
Automatic configuration with Bootp	209
About Agilent BootP Service	209
How BootP Service Works	210
Situation: Cannot Establish LAN Communication	210
Installation of BootP Service	211
Two Methods to Determine the MAC Address	213
Assigning IP Addresses Using the Agilent BootP Service	214
Changing the IP Address of an Instrument Using the Agilent BootP Service	217
Manual Configuration	219
With Telnet	220
With the Instant Pilot (G4208A)	224
PC and User Interface Software Setup	225
PC Setup for Local Configuration	225
User Interface Software Setup	226

This chapter provides information on connecting the detector to the Agilent ChemStation PC.




## 13 LAN Configuration

What you have to do first

### What you have to do first

The module has an on-board LAN communication interface.

- 1 Note the MAC (Media Access Control) address for further reference. The MAC or hardware address of the LAN interfaces is a world wide unique identifier. No other network device will have the same hardware address. The MAC address can be found on a label at the rear of the module underneath the configuration switch.

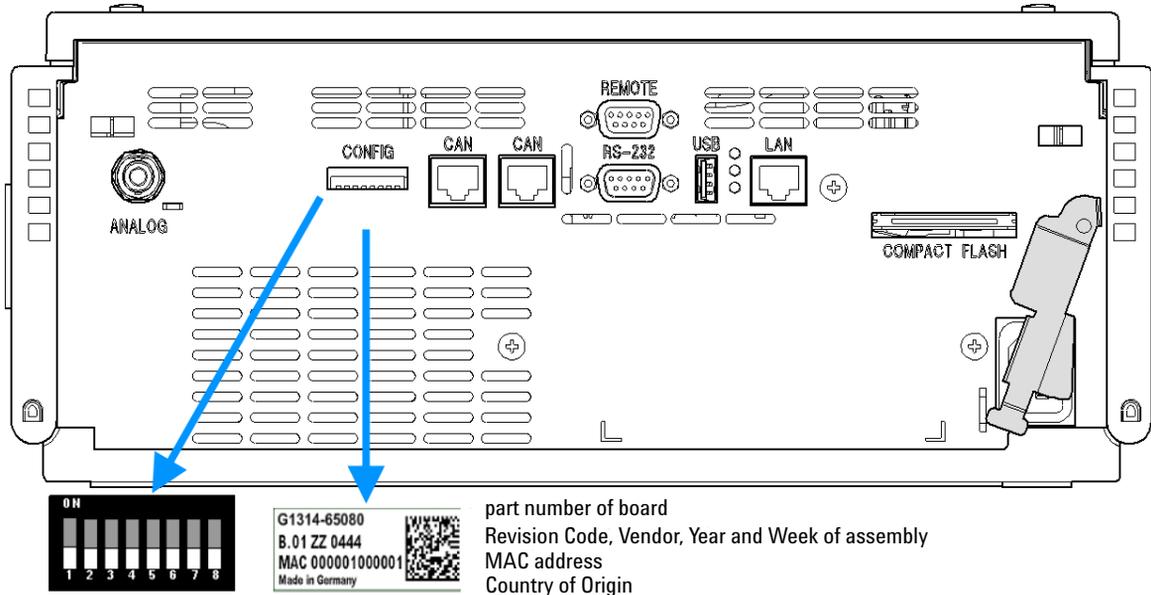

**Figure 60** Location of Configuration Switch and MAC Label

- 2 Connect the instrument's LAN interface to
  - the PC network card using a crossover network cable (point-to-point) or
  - a hub or switch using a standard LAN cable.

## TCP-IP parameter configuration

To operate properly in a network environment, the LAN interface must be configured with valid TCP/IP network parameters. These parameters are:

- IP address
- Subnet Mask
- Default Gateway


The TCP/IP parameters can be configured by the following methods:

- by automatically requesting the parameters from a network-based BOOTP Server (using the so-called Bootstrap Protocol)
- by automatically requesting the parameters from a network-based DHCP Server (using the so-called Dynamic Host Configuration Protocol). This mode requires a LAN-onboard Module or a G1369C LAN Interface card, see “[Setup \(DHCP\)](#)” on page 206
- by manually setting the parameters using Telnet
- by manually setting the parameters using the Instant Pilot (G4208A)

The LAN interface differentiates between several initialization modes. The initialization mode (short form ‘init mode’) defines how to determine the active TCP/IP parameters after power-on. The parameters may be derived from a Bootp cycle, non-volatile memory or initialized with known default values. The initialization mode is selected by the configuration switch, see [Table 25](#) on page 201.

## Configuration Switches

The configuration switch can be accessed at the rear of the module, see figure below.

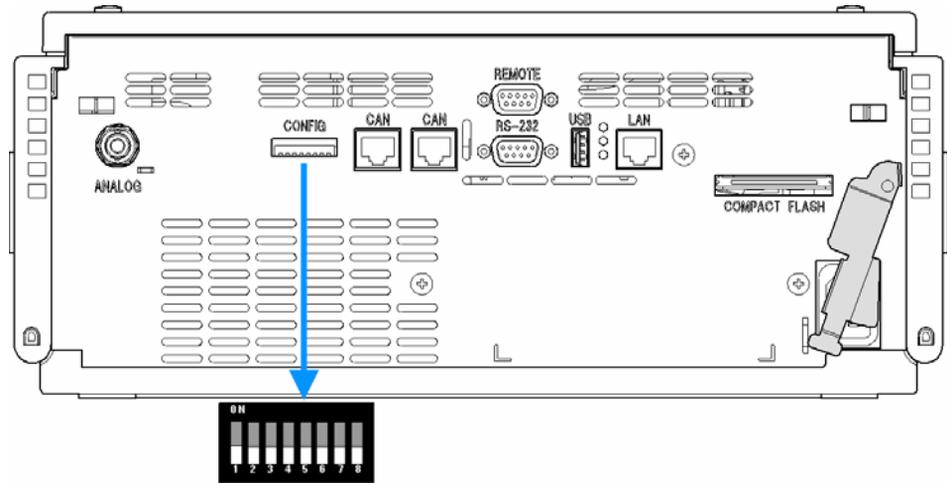

**Figure 61** Location of Configuration Switch

The module is shipped with all switches set to OFF, as shown above.

**NOTE**

To perform any LAN configuration, SW1 and SW2 must be set to OFF.


**Table 24** Factory Default Settings

Initialization ('Init') Mode	Bootp, all switches down. For details see <a href="#">Figure 62</a> on page 201
Link Configuration	speed and duplex mode determined by auto-negotiation, for details see " <a href="#">Link configuration selection</a> " on page 208

## Initialization mode selection

The following initialization (init) modes are selectable:


**Table 25** Initialization Mode Switches

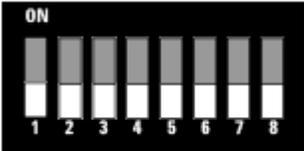	SW 6	SW 7	SW 8	Init Mode
	OFF	OFF	OFF	Bootp
	OFF	OFF	ON	Bootp & Store
	OFF	ON	OFF	Using Stored
	OFF	ON	ON	Using Default
	ON	OFF	OFF	DHCP <sup>1</sup>

<sup>1</sup> Requires firmware B.06.40 or above. Modules without LAN on board, see G1369C LAN Interface Card

### Bootp

When the initialization mode **Bootp** is selected, the module tries to download the parameters from a **Bootp** Server. The parameters obtained become the active parameters immediately. They are not stored to the non-volatile memory of the module. Therefore, the parameters are lost with the next power cycle of the module.

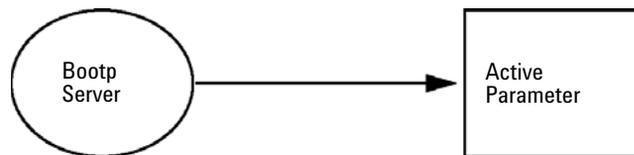


**Figure 62** Bootp (Principle)

## Bootp & Store

When **Bootp & Store** is selected, the parameters obtained from a **Bootp** Server become the active parameters immediately. In addition, they are stored to the non-volatile memory of the module. Thus, after a power cycle they are still available. This enables a kind of bootp once configuration of the module.

*Example:* The user may not want to have a **Bootp** Server be active in his network all the time. But on the other side, he may not have any other configuration method than **Bootp**. In this case he starts the **Bootp** Server temporarily, powers on the module using the initialization mode **Bootp & Store**, waits for the **Bootp** cycle to be completed, closes the **Bootp** Server and powers off the module. Then he selects the initialization mode Using Stored and powers on the module again. From now on, he is able to establish the TCP/IP connection to the module with the parameters obtained in that single **Bootp** cycle.

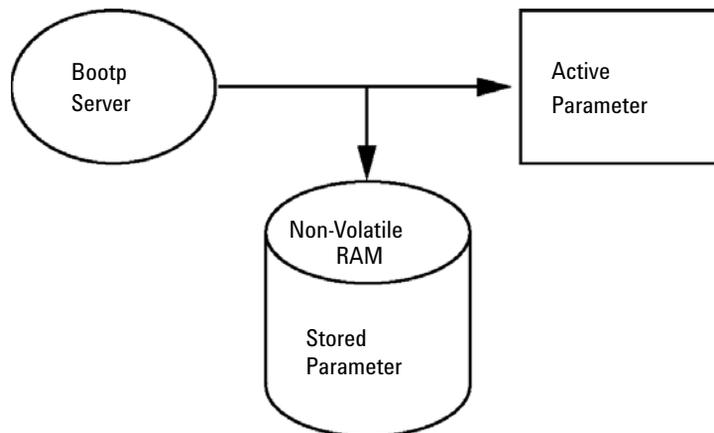


**Figure 63** Bootp & Store (Principle)

### NOTE

Use the initialization mode **Bootp & Store** carefully, because writing to the non-volatile memory takes time. Therefore, when the module shall obtain its parameters from a **Bootp** Server every time it is powered on, the recommended initialization mode is **Bootp**!

### Using Stored


When initialization mode **Using Stored** is selected, the parameters are taken from the non-volatile memory of the module. The TCP/IP connection will be established using these parameters. The parameters were configured previously by one of the described methods.

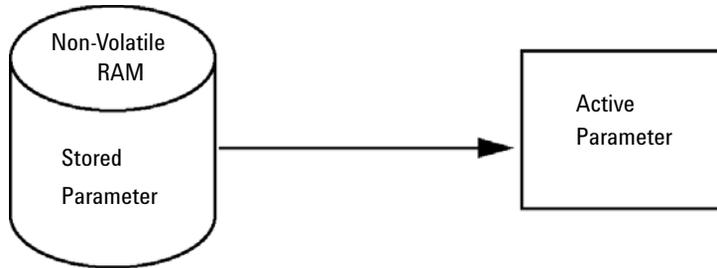

**Figure 64** Using Stored (Principle)

### Using Default

When **Using Default** is selected, the factory default parameters are taken instead. These parameters enable a TCP/IP connection to the LAN interface without further configuration, see [Table 26](#) on page 204.

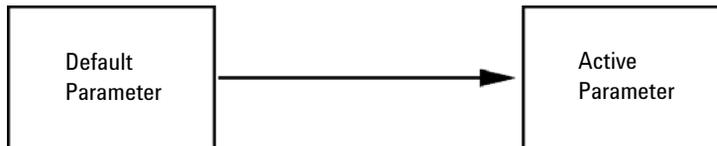

**Figure 65** Using Default (Principle)

**NOTE**

Using the default address in your local area network may result in network problems. Take care and change it to a valid address immediately.

## 13 LAN Configuration

### Initialization mode selection

**Table 26** Using Default Parameters

IP address:	192.168.254.11
Subnet Mask:	255.255.255.0
Default Gateway	not specified

Since the default IP address is a so-called local address, it will not be routed by any network device. Thus, the PC and the module must reside in the same subnet.

The user may open a Telnet session using the default IP address and change the parameters stored in the non-volatile memory of the module. He may then close the session, select the initialization mode Using Stored, power-on again and establish the TCP/IP connection using the new parameters.

When the module is wired to the PC directly (e.g. using a cross-over cable or a local hub), separated from the local area network, the user may simply keep the default parameters to establish the TCP/IP connection.

#### NOTE

In the **Using Default** mode, the parameters stored in the memory of the module are not cleared automatically. If not changed by the user, they are still available, when switching back to the mode Using Stored.


# Dynamic Host Configuration Protocol (DHCP)

## General Information (DHCP)

The Dynamic Host Configuration Protocol (DHCP) is an auto configuration protocol used on IP networks. The DHCP functionality is available on all Agilent HPLC modules with on-board LAN Interface or LAN Interface Card, and “B”-firmware (B.06.40 or above).

When the initialization mode “DHCP” is selected, the card tries to download the parameters from a DHCP Server. The parameters obtained become the active parameters immediately. They are not stored to the non-volatile memory of the card.

Besides requesting the network parameters, the card also submits its hostname to the DHCP Server. The hostname equals the MAC address of the card, e.g. *0030d3177321*. It is the DHCP server's responsibility to forward the hostname/address information to the Domain Name Server. The card does not offer any services for hostname resolution (e.g. NetBIOS).

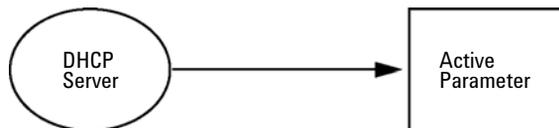

**Figure 66** DHCP (Principle)

### NOTE

- 1 It may take some time until the DHCP server has updated the DNS server with the hostname information.
- 2 It may be necessary to fully qualify the hostname with the DNS suffix, e.g. *0030d3177321.country.company.com*.
- 3 The DHCP server may reject the hostname proposed by the card and assign a name following local naming conventions.

## 13 LAN Configuration


### Dynamic Host Configuration Protocol (DHCP)

## Setup (DHCP)

**Software required** The modules in the stack must have at least firmware from set A.06.34 and the above mentioned modules B.06.40 or above (must from the same firmware set).

- 1 Note the MAC address of the LAN interface (provided with G1369C LAN Interface Card or Main Board). This MAC address is on a label on the card or at the rear of the main board, e.g. *0030d3177321*.

On the Instant Pilot the MAC address can be found under **Details** in the LAN section.

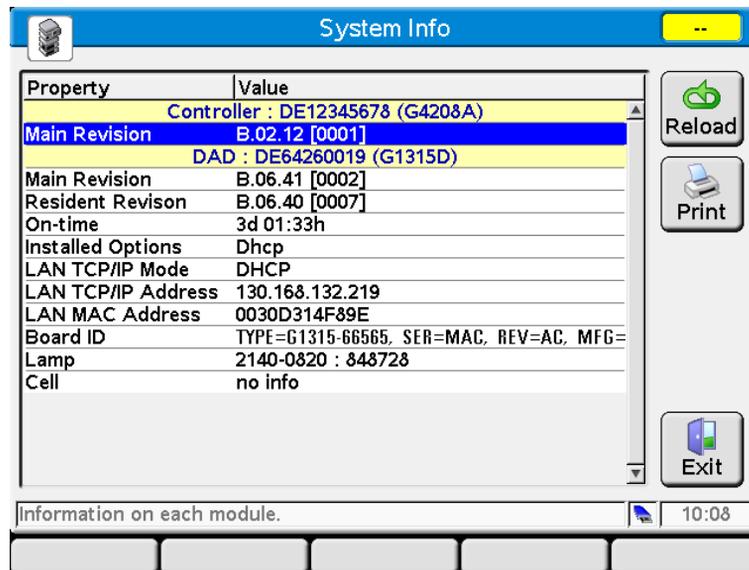

**Figure 67** LAN Setting on Instant Pilot

- 2 Set the Configuration Switch to DHCP either on the G1369C LAN Interface Card or the main board of above mentioned modules.

**Table 27** G1369C LAN Interface Card (configuration switch on the card)

SW 4	SW 5	SW 6	SW 7	SW 8	Initialization Mode
ON	OFF	OFF	OFF	OFF	DHCP

**Table 28** LC Modules inclusive 1120/1220 (configuration switch at rear of the instrument)

SW 6	SW 7	SW 8	Initialization Mode
ON	OFF	OFF	DHCP

- 3 Turn on the module that hosts the LAN interface.
- 4 Configure your Control Software (e.g. Agilent ChemStation, Lab Advisor, Firmware Update Tool) and use MAC address as host name, e.g. *0030d3177321*.

The LC system should become visible in the control software (see Note in section “[General Information \(DHCP\)](#)” on page 205).

## 13 LAN Configuration

### Link configuration selection

## Link configuration selection

The LAN interface supports 10 or 100 Mbps operation in full- or half-duplex modes. In most cases, full-duplex is supported when the connecting network device - such as a network switch or hub - supports IEEE 802.3u auto-negotiation specifications.

When connecting to network devices that do not support auto-negotiation, the LAN interface will configure itself for 10- or 100-Mbps half-duplex operation.

For example, when connected to a non-negotiating 10-Mbps hub, the LAN interface will be automatically set to operate at 10-Mbps half-duplex.

If the module is not able to connect to the network through auto-negotiation, you can manually set the link operating mode using link configuration switches on the module.

**Table 29** Link Configuration Switches

	SW 3	SW 4	SW 5	Link Configuration
	OFF	-	-	speed and duplex mode determined by auto-negotiation
	ON	OFF	OFF	manually set to 10 Mbps, half-duplex
	ON	OFF	ON	manually set to 10 Mbps, full-duplex
	ON	ON	OFF	manually set to 100 Mbps, half-duplex
	ON	ON	ON	manually set to 100 Mbps, full-duplex

## Automatic configuration with Bootp

### NOTE

All examples shown in this chapter will not work in your environment. You need your own IP-, Subnet-Mask- and Gateway addresses.

---

### NOTE

Assure that the detector configuration switch is set properly. The setting should be either **BootP** or **BootP & Store**, see [Table 25](#) on page 201.

---

### NOTE

Assure that the detector connected to the network is powered off.

---

### NOTE

If the Agilent BootP Service program is not already installed on your PC, then install it from your Agilent ChemStation DVD, located in folder **BootP**.

---

## About Agilent BootP Service

The Agilent BootP Service is used to assign the LAN Interface with an IP address.

The Agilent BootP Service is provided on the ChemStation DVD. The Agilent BootP Service is installed on a server or PC on the LAN to provide central administration of IP addresses for Agilent instruments on a LAN. The BootP service must be running TCP/IP network protocol and cannot run a DHCP server.

## How BootP Service Works

When an instrument is powered on, an LAN Interface in the instrument broadcasts a request for an IP address or host name and provides its hardware MAC address as an identifier. The Agilent BootP Service answers this request and passes a previously defined IP address and host name associated with the hardware MAC address to the requesting instrument.

The instrument receives its IP address and host name and maintains the IP address as long as it is powered on. Powering down the instrument causes it to lose its IP address, so the Agilent BootP Service must be running every time the instrument powers up. If the Agilent BootP Service runs in the background, the instrument will receive its IP address on power-up.

The Agilent LAN Interface can be set to store the IP address and will not lose the IP address if power cycled.

## Situation: Cannot Establish LAN Communication

If a LAN communication with BootP service cannot be established, check the following on the PC:

- Is the BootP service started? During installation of BootP, the service is not started automatically.
- Does the Firewall block the BootP service? Add the BootP service as an exception.
- Is the LAN Interface using the BootP-mode instead of "Using Stored" or "Using Default" modes?

## Installation of BootP Service


Before installing and configuring the Agilent BootP Service, be sure to have the IP addresses of the computer and instruments on hand.

- 1 Log on as Administrator or other user with Administrator privileges.
- 2 Close all Windows programs.
- 3 Insert the Agilent ChemStation software DVD into the drive. If the setup program starts automatically, click **Cancel** to stop it.
- 4 Open Windows Explorer.
- 5 Go to the BootP directory on the Agilent ChemStation DVD and double-click **BootPPackage.msi**.
- 6 If necessary, click the **Agilent BootP Service...** icon in the task bar.
- 7 The **Welcome** screen of the **Agilent BootP Service Setup Wizard** appears. Click **Next**.
- 8 The **End-User License Agreement** screen appears. Read the terms, indicate acceptance, then click **Next**.
- 9 The **Destination Folder** selection screen appears. Install BootP to the default folder or click **Browse** to choose another location. Click **Next**.  
The default location for installation is:  
C:\Program Files\Agilent\BootPService\  
10 Click **Install** to begin installation.

## 13 LAN Configuration

### Automatic configuration with Bootp

11 Files load; when finished, the **BootP Settings** screen appears.

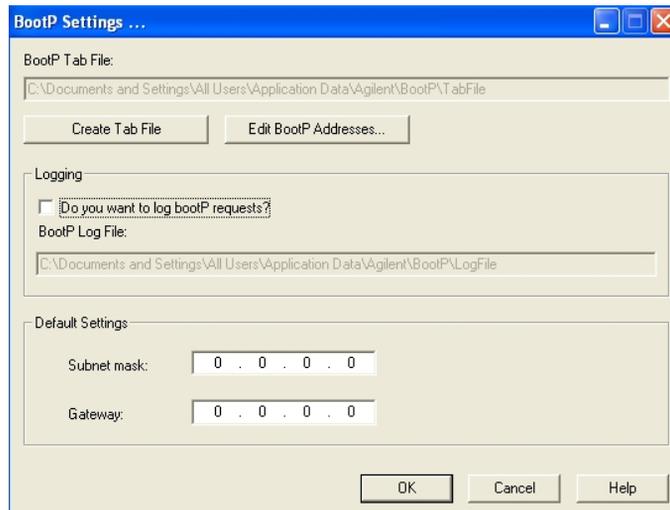

**Figure 68** BootP Settings screen

12 In the **Default Settings** part of the screen, if known, you can enter the subnet mask and gateway.

Defaults can be used:

- The default subnet mask is 255.255.255.0
- The default gateway is 192.168.254.11

13 On the **BootP Settings** screen, click **OK**. The **Agilent BootP Service Setup** screen indicates completion.

14 Click **Finish** to exit the **Agilent BootP Service Setup** screen.

15 Remove the DVD from the drive.

This completes installation.

16 Start BootP Service in the Windows® services: On the Windows® desktop click right on **Computer** icon, select **Manage > Services and Applications > Services**. Select the **Agilent BootP Service** and click **Start**.

## Two Methods to Determine the MAC Address

### Enabling logging to discover the MAC address using BootP

If you want to see the MAC address, select the **Do you want to log BootP requests?** check box.

- 1 Open BootP Settings from **Start > All Programs > Agilent BootP Service > EditBootPSettings**.
- 2 In **BootP Settings...** check **Do you want to log BootP requests?** to enable logging.



**Figure 69** Enable BootP logging

The log file is located in

C:\Documents and Settings\All Users\Application Data\Agilent\BootP\LogFile

It contains a MAC address entry for each device that requests configuration information from BootP.

- 3 Click **OK** to save the values or **Cancel** to discard them. The editing ends.
- 4 After each modification of the BootP settings (i.e. **EditBootPSettings**) a stop or start of the BootP service is required for the BootP service to accept changes. See “[Stopping the Agilent BootP Service](#)” on page 217 or “[Restarting the Agilent BootP Service](#)” on page 218.
- 5 Uncheck the **Do you want to log BootP requests?** box after configuring instruments; otherwise, the log file will quickly fill up disk space.

### Determining the MAC address directly from the LAN Interface card label

- 1 Turn off the instrument.
- 2 Read the MAC address from the label and record it.  
The MAC address is printed on a label on the rear of the module. It is the number below the barcode and after the colon (:) and usually begins with the letters AD, see [Figure 60](#) on page 198 .
- 3 Turn on the instrument.

## Assigning IP Addresses Using the Agilent BootP Service

The Agilent BootP Service assigns the Hardware MAC address of the instrument to an IP address.

### Determining the MAC address of the instrument using BootP Service

- 1 Power cycle the Instrument.
- 2 After the instrument completes self-test, open the log file of the BootP Service using Notepad.
  - The default location for the logfile is C:\Documents and Settings\All Users\Application Data\Agilent\BootP\LogFile.
  - The logfile will not be updated if it is open.

The contents will be similar to the following:

02/25/10 15:30:49 PM

Status: BootP Request received at outermost layer

Status: BootP Request received from hardware address: 0010835675AC

Error: Hardware address not found in BootPTAB: 0010835675AC

Status: BootP Request finished processing at outermost layer


- 3 Record the hardware (MAC) address (for example, 0010835675AC).
- 4 The Error means the MAC address has not been assigned an IP address and the Tab File does not have this entry. The MAC address is saved to the Tab File when an IP address is assigned.
- 5 Close the log file before turning on another instrument.
- 6 Uncheck the **Do you want to log BootP requests?** box after configuring instruments to avoid having the logfile use up excessive disk space.

## Adding each instrument to the network using BootP

- 1 Follow **Start > All Programs > Agilent BootP Service** and select **Edit BootP Settings**. The BootP Settings screen appears.
- 2 Uncheck the **Do you want to log BootP requests?** once all instruments have been added.

The **Do you want to log BootP requests?** box must be unchecked when you have finished configuring instruments; otherwise, the log file will quickly fill up disk space.

- 3 Click **Edit BootP Addresses...** The **Edit BootP Addresses** screen appears.
- 4 Click **Add...** The **Add BootP Entry** screen appears.

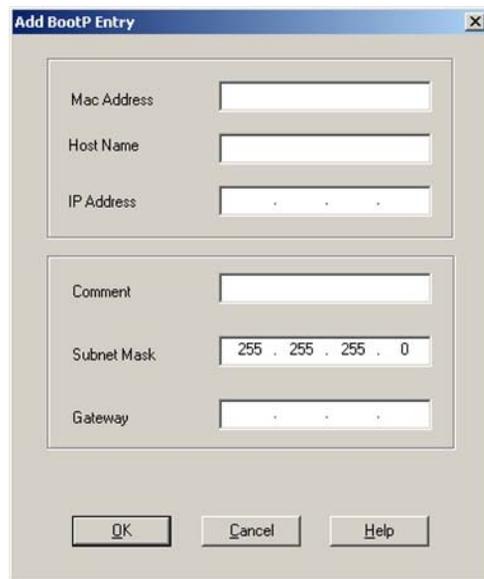

**Figure 70** Enable BootP logging

- 5 Make these entries for the instrument:
  - MAC address
  - Host name, Enter a Hostname of your choice.  
The Host Name must begin with "alpha" characters (i.e. LC1260)
  - IP address
  - Comment (optional)

## 13 LAN Configuration

### Automatic configuration with Bootp

- Subnet mask
- Gateway address (optional)

The configuration information entered is saved in the Tab File.

**6** Click **OK**.

**7** Leave **Edit BootP Addresses** by pressing **Close**.

**8** Exit **BootP Settings** by pressing **OK**.

**9** After each modification of the BootP settings (i.e. EditBootPSettings) a stop or start of the BootP service is required for the BootP service to accept changes. See “[Stopping the Agilent BootP Service](#)” on page 217 or “[Restarting the Agilent BootP Service](#)” on page 218.

**10** Power cycle the Instrument.

OR

If you changed the IP address, power cycle the instrument for the changes to take effect.

**11** Use the PING utility to verify connectivity by opening a command window and typing:

Ping 192.168.254.11 for example.

The Tab File is located at


C:\Documents and Settings\All Users\Application Data\Agilent\BootP\TabFile

## Changing the IP Address of an Instrument Using the Agilent BootP Service

Agilent BootP Service starts automatically when your PC reboots. To change Agilent BootP Service settings, you must stop the service, make the changes, and then restart the service.

### Stopping the Agilent BootP Service

- 1 From the Windows control panel, select **Administrative Tools > Services**. The **Services** screen appears.

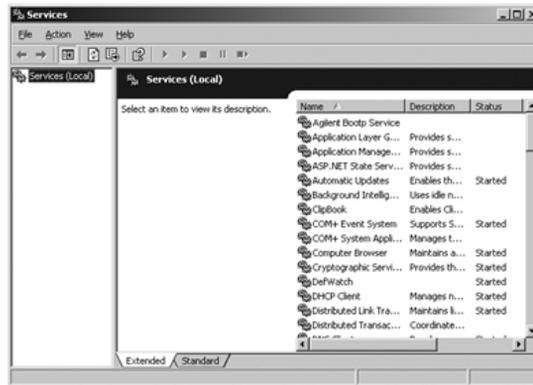

**Figure 71** Windows Services screen


- 2 Right-click **Agilent BootP Service**.
- 3 Select **Stop**.
- 4 Close the **Services and Administrative Tools** screen.

## 13 LAN Configuration

### Automatic configuration with Bootp

#### Editing the IP address and other parameters in EditBootPSettings

- 1 Select **Start > All Programs > Agilent BootP Service** and select **Edit BootP Settings**. The **BootP Settings** screen appears.
- 2 When the **BootP Settings** screen is first opened, it shows the default settings from installation.
- 3 Press **Edit BootP Addresses...** to edit the Tab File.

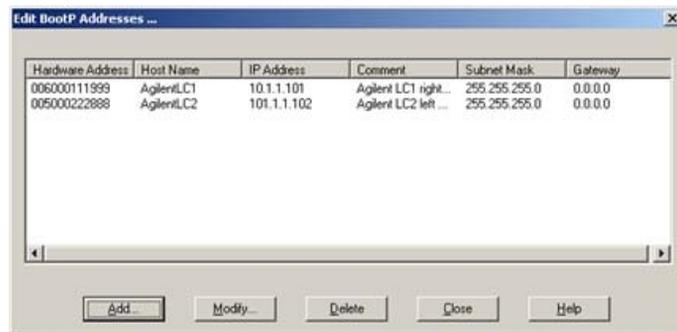

**Figure 72** Edit BootP Adresses screen

- 4 In the **Edit BootP Addresses...** screen press **Add...** to create a new entry or select an existing line from the table and press **Modify...** or **Delete** to change the IP address, comment, subnet mask, for example, in the Tab File.

If you change the IP address, it will be necessary to power cycle the instrument for the changes to take effect.


- 5 Leave **Edit BootP Addresses...** by pressing **Close**.
- 6 Exit **BootP Settings** by pressing **OK**.

#### Restarting the Agilent BootP Service

- 1 In the Windows control panel, select **Administrative Tools > Services**. The **Services** screen appears, see [Figure 71](#) on page 217.
- 2 Right-click **Agilent BootP Service** and select **Start**.
- 3 Close the **Services and Administrative Tools** screens.

## Manual Configuration

Manual configuration only alters the set of parameters stored in the non-volatile memory of the module. It never affects the currently active parameters. Therefore, manual configuration can be done at any time. A power cycle is mandatory to make the stored parameters become the active parameters, given that the initialization mode selection switches are allowing it.

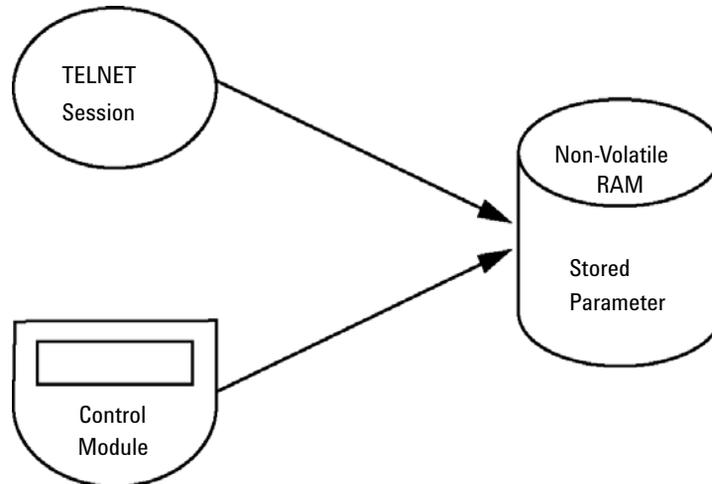


**Figure 73** Manual Configuration (Principle)

## With Telnet

Whenever a TCP/IP connection to the module is possible (TCP/IP parameters set by any method), the parameters may be altered by opening a Telnet session.

- 1 Open the system (DOS) prompt window by clicking on Windows **START** button and select “**Run...**”. Type “cmd” and press OK.
- 2 Type the following at the system (DOS) prompt:
  - `c:\>telnet <IP address>` or
  - `c:\>telnet <host name>`

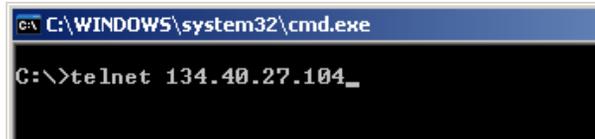


```
C:\WINDOWS\system32\cmd.exe
C:\>telnet 134.40.27.104_
```

Figure 74 Telnet - Starting a session

where <IP address> may be the assigned address from a Bootp cycle, a configuration session with the Local Controller, or the default IP address (see “[Configuration Switches](#)” on page 200).


When the connection was established successfully, the module responds with the following:

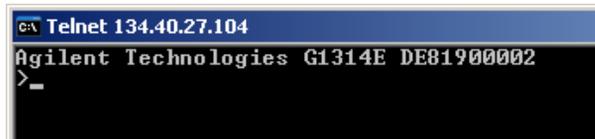

```
C:\ Telnet 134.40.27.104
Agilent Technologies G1314E DE81900002
>_
```

Figure 75 A connection to the module is made

- 3 Type ? and press enter to see the available commands.

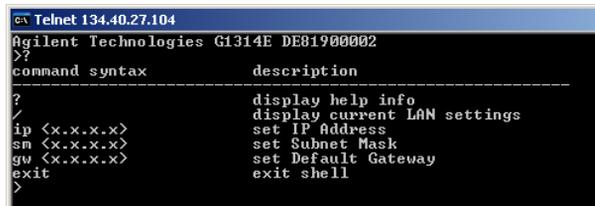

```
C:\ Telnet 134.40.27.104
Agilent Technologies G1314E DE81900002
>?
command syntax      description
-----
?                    display help info
/                    display current LAN settings
ip <x.x.x.x>          set IP Address
sm <x.x.x.x>          set Subnet Mask
gw <x.x.x.x>          set Default Gateway
exit                 exit shell
>
```

Figure 76 Telnet Commands

**Table 30** Telnet Commands

Value	Description
?	displays syntax and descriptions of commands
/	displays current LAN settings
ip <x.x.x.x>	sets new ip address
sm <x.x.x.x>	sets new subnet mask
gw <x.x.x.x>	sets new default gateway
exit	exits shell and saves all changes

**4** To change a parameter follows the style:

- parameter value, for example:  
**ip 134.40.27.230**

Then press [Enter], where parameter refers to the configuration parameter you are defining, and value refers to the definitions you are assigning to that parameter. Each parameter entry is followed by a carriage return.

## 13 LAN Configuration

### Manual Configuration

- 5 Use the “/” and press Enter to list the current settings.

```
c:\ Telnet 134.40.27.104
Agilent Technologies G1314E DE81900002
>/
LAN Status Page
-----
MAC Address      : 0030D30611BD
Init Mode        : Using Stored
-----
TCP/IP Properties
- active -
IP Address       : 134.40.27.104
Subnet Mask      : 255.255.248.0
Def. Gateway     : 134.40.24.1
-----
TCP/IP Status   : Ready
-----
Controllers     : no connections
>_
```

information about the LAN interface

MAC address, initialization mode

Initialization mode is Using Stored

active TCP/IP settings

TCP/IP status - here ready

connected to PC with controller software

here not connected

**Figure 77** Telnet - Current settings in “Using Stored” mode

- 6 Change the IP address (in this example 134.40.27.99) and type “/” to list current settings.

```
c:\ Telnet 134.40.27.104
>ip 134.40.27.99
>/
LAN Status Page
-----
MAC Address      : 0030D30611BD
Init Mode        : Using Stored
-----
TCP/IP Properties
- active -
IP Address       : 134.40.27.104
Subnet Mask      : 255.255.248.0
Def. Gateway     : 134.40.24.1
- stored -
IP Address       : 134.40.27.99
Subnet Mask      : 255.255.248.0
Def. Gateway     : 134.40.24.1
-----
TCP/IP Status   : Ready
-----
Controllers     : no connections
>_
```

change of IP setting to

Initialization mode is Using Stored

active TCP/IP settings

stored TCP/IP settings in non-volatile memory


TCP/IP status - here ready

connected to PC with controller software

here not connected

**Figure 78** Telnet - Change IP settings

- 7 When you have finished typing the configuration parameters, type **exit** and press [Enter] to exit with storing parameters.

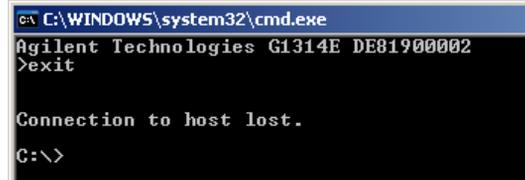

```
C:\WINDOWS\system32\cmd.exe
Agilent Technologies G1314E DE81900002
>exit

Connection to host lost.
C:\>
```

**Figure 79** Closing the Telnet Session


**NOTE**

If the Initialization Mode Switch is changed now to “Using Stored” mode, the instrument will take the stored settings when the module is re-booted.

## With the Instant Pilot (G4208A)

To configure the TCP/IP parameters before connecting the detector to the network, the Instant Pilot (G4208A) can be used.

- 1 From the Welcome screen press the **More** button.
- 2 Select **Configure**.
- 3 Press the **VWD** button.
- 4 Scroll down to the LAN settings.

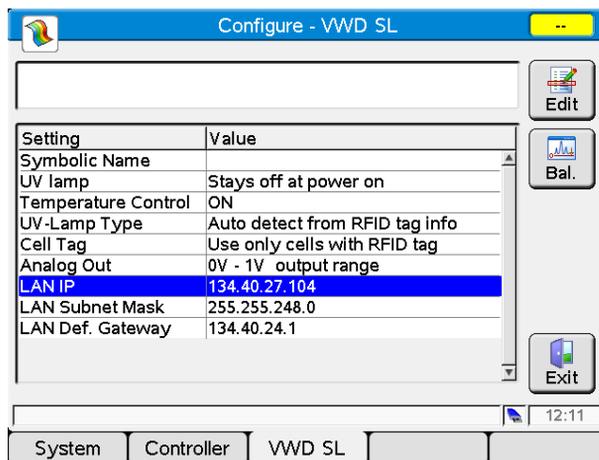

**Figure 80** Instant Pilot - LAN Configuration


- 5 Press the **Edit** button (only visible if not in Edit mode), perform the required changes and press the **Done** button.
- 6 Leave the screen by clicking **Exit**.

# PC and User Interface Software Setup

## PC Setup for Local Configuration

This procedure describes the change of the TCP/IP settings on your PC to match the module’s default parameters in a local configuration (see also “Initialization mode selection” on page 201).

FIXED IP address

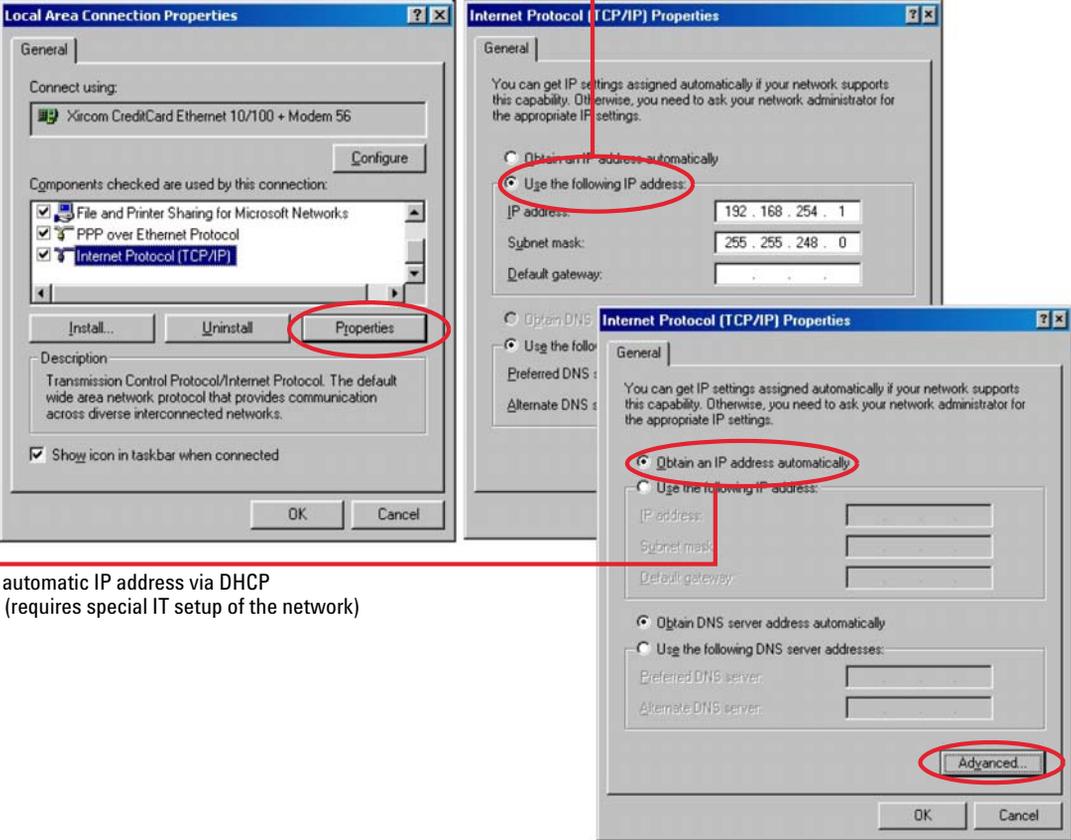

automatic IP address via DHCP  
(requires special IT setup of the network)


Figure 81 Changing the TCP/IP settings of the PC

## 13 LAN Configuration

### PC and User Interface Software Setup

#### User Interface Software Setup

Install you user interface software according the provided *User Interface Software Setup Guide*.

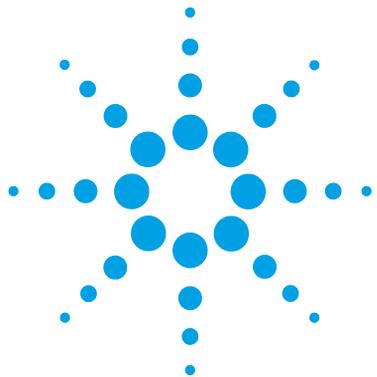

## 14 Appendix

General Safety Information	228
Safety Symbols	228
General Safety Information	229
Safety Standards	229
Operation	229
Radio Interference	231
Sound Emission	232
UV Radiation	233
Solvent Information	234
Declaration of Conformity for HOX2 Filter	236
Agilent Technologies on Internet	237

This chapter provides addition information on safety, legal and web.



## General Safety Information

### Safety Symbols

Table 31 Safety Symbols

Symbol	Description
	The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.
	Indicates dangerous voltages.
	Indicates a protected ground terminal.
	Indicates eye damage may result from directly viewing the light produced by the deuterium lamp used in this product.
	The apparatus is marked with this symbol when hot surfaces are available and the user should not touch it when heated up.

#### WARNING

##### A WARNING

alerts you to situations that could cause physical injury or death.

- Do not proceed beyond a warning until you have fully understood and met the indicated conditions.

#### CAUTION

##### A CAUTION

alerts you to situations that could cause loss of data, or damage of equipment.

- Do not proceed beyond a caution until you have fully understood and met the indicated conditions.

## General Safety Information

The following general safety precautions must be observed during all phases of operation, service, and repair of this instrument. Failure to comply with these precautions or with specific warnings elsewhere in this manual violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements.

### WARNING

**Ensure the proper usage of the equipment.**

**The protection provided by the equipment may be impaired.**

→ The operator of this instrument is advised to use the equipment in a manner as specified in this manual.

---

## Safety Standards

This is a Safety Class I instrument (provided with terminal for protective earthing) and has been manufactured and tested according to international safety standards.

## Operation

Before applying power, comply with the installation section. Additionally the following must be observed.

Do not remove instrument covers when operating. Before the instrument is switched on, all protective earth terminals, extension cords, auto-transformers, and devices connected to it must be connected to a protective earth via a ground socket. Any interruption of the protective earth grounding will cause a potential shock hazard that could result in serious personal injury. Whenever it is likely that the protection has been impaired, the instrument must be made inoperative and be secured against any intended operation.

Make sure that only fuses with the required rated current and of the specified type (normal blow, time delay, and so on) are used for

replacement. The use of repaired fuses and the short-circuiting of fuse holders must be avoided.

Some adjustments described in the manual, are made with power supplied to the instrument, and protective covers removed. Energy available at many points may, if contacted, result in personal injury.

Any adjustment, maintenance, and repair of the opened instrument under voltage should be avoided whenever possible. When inevitable, this has to be carried out by a skilled person who is aware of the hazard involved. Do not attempt internal service or adjustment unless another person, capable of rendering first aid and resuscitation, is present. Do not replace components with power cable connected.

Do not operate the instrument in the presence of flammable gases or fumes. Operation of any electrical instrument in such an environment constitutes a definite safety hazard.

Do not install substitute parts or make any unauthorized modification to the instrument.

Capacitors inside the instrument may still be charged, even though the instrument has been disconnected from its source of supply. Dangerous voltages, capable of causing serious personal injury, are present in this instrument. Use extreme caution when handling, testing and adjusting.

When working with solvents, observe appropriate safety procedures (for example, goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet by the solvent vendor, especially when toxic or hazardous solvents are used.

## Radio Interference

Cables supplied by Agilent Technologies are screened to provide optimized protection against radio interference. All cables are in compliance with safety or EMC regulations.

### Test and Measurement

If test and measurement equipment is operated with unscreened cables, or used for measurements on open set-ups, the user has to assure that under operating conditions the radio interference limits are still met within the premises.

## Sound Emission

### Manufacturer's Declaration

This statement is provided to comply with the requirements of the German Sound Emission Directive of 18 January 1991.

This product has a sound pressure emission (at the operator position) < 70 dB.

- Sound Pressure  $L_p < 70$  dB (A)
- At Operator Position
- Normal Operation
- According to ISO 7779:1988/EN 27779/1991 (Type Test)

## UV Radiation

Emissions of ultraviolet radiation (200 – 315 nm) from this product is limited such that radiant exposure incident upon the unprotected skin or eye of operator or service personnel is limited to the following TLVs (Threshold Limit Values) according to the American Conference of Governmental Industrial Hygienists:

**Table 32** UV radiation limits

Exposure/day	Effective irradiance
8 h	0.1 $\mu\text{W}/\text{cm}^2$
10 min	5.0 $\mu\text{W}/\text{cm}^2$

Typically the radiation values are much smaller than these limits:

**Table 33** UV radiation typical values

Position	Effective irradiance
Lamp installed, 50 cm distance	average 0.016 $\mu\text{W}/\text{cm}^2$
Lamp installed, 50 cm distance	maximum 0.14 $\mu\text{W}/\text{cm}^2$

## Solvent Information

Observe the following recommendations on the use of solvents.

### Flow Cell

Avoid the use of alkaline solutions (pH > 9.5 ) which can attack quartz and thus impair the optical properties of the flow cell.

Prevent any crystallization of buffer solutions. This will lead into a blockage/damage of the flow cell.

If the flow cell is transported while temperatures are below 5 °C, it must be assured that the cell is filled with alcohol.


Aqueous solvents in the flow cell can built up algae. Therefore do not leave aqueous solvents sitting in the flow cell. Add small % of organic solvents (for example, acetonitrile or methanol ~5 %).

### Solvents

Brown glass ware can avoid growth of algae.

Always filter solvents, small particles can permanently block the capillaries. Avoid the use of the following steel-corrosive solvents:

- Solutions of alkali halides and their respective acids (for example, lithium iodide, potassium chloride, and so on).
- High concentrations of inorganic acids like nitric acid, sulfuric acid especially at higher temperatures (replace, if your chromatography method allows, by phosphoric acid or phosphate buffer which are less corrosive against stainless steel).
- Halogenated solvents or mixtures which form radicals and/or acids, for example:

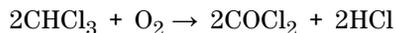

This reaction, in which stainless steel probably acts as a catalyst, occurs quickly with dried chloroform if the drying process removes the stabilizing alcohol.

- Chromatographic grade ethers, which can contain peroxides (for example, THF, dioxane, di-isopropylether) such ethers should be filtered through dry aluminium oxide which adsorbs the peroxides.
- Solutions of organic acids (acetic acid, formic acid, and so on) in organic solvents. For example, a 1 % solution of acetic acid in methanol will attack steel.
- Solutions containing strong complexing agents (for example, EDTA, ethylene diamine tetra-acetic acid).
- Mixtures of carbon tetrachloride with 2-propanol or THF.

# Declaration of Conformity for HOX2 Filter

Declaration of Conformity				
We herewith inform you that the				
<b>Holmium Oxide Glass Filter</b>				
used in Agilent's absorbance detectors listed in the table below meets the requirements of National Institute of Standards and Technology (NIST) to be applied as certified wavelength standard.				
According to the publication of NIST in J. Res. Natl. Inst. Stand. Technol. 112, 303-306 (2007) the holmium oxide glass filters are inherently stable with respect to the wavelength scale and need no recertification. The expanded uncertainty of the certified wavelength values is 0.2 nm.				
Agilent Technologies guarantees, as required by NIST, that the material of the filters is holmium oxide glass representing the inherently existent holmium oxide absorption bands.				
Test wavelengths:				
Product Number	Series	Measured Wavelength *	Wavelength Accuracy	Optical Bandwidth
79883A	1090	361.0 nm	+/- 1 nm	2 nm
79854A	1050	418.9 nm		
G1306A	1050	453.7 nm		
G1315A, G1365A	1100	536.7 nm		
G1315B/C, G1365B/C	1100 / 1200 / 1260			
G1600A, G7100A	CE			
79853C	1050	360.8nm 418.5nm 536.4nm	+/- 2 nm	6 nm
G1314A/B/C	1100 / 1200 / 1260	360.8nm 418.5nm	+/- 1 nm	6 nm
G1314D/E/F		418.5nm		
G4286..... 90A/B/C	1120 / 1220	536.4nm		
*) The variation in Measured Wavelength depends on the different Optical Bandwidth.				
May 19, 2010				
----- (Date)				
 ----- (R&D Manager)		 ----- (Quality Manager)		
P/N 89550-90501 		Revision: H Effective by: May 19, 2010		

## Agilent Technologies on Internet

For the latest information on products and services visit our worldwide web site on the Internet at:

<http://www.agilent.com>

# Index

## 8

8-bit configuration switch  
on-board LAN 189

## A

absorbance  
Lambert-Beer 86  
Agilent Lab Advisor software 94  
Agilent Lab Advisor 94  
Agilent  
on internet 237  
User Interface Software Setup 225  
alga information 234  
ambient non-operating temperature 23  
ambient operating temperature 23  
analog out  
put 31, 28, 25  
analog signal 186  
analog  
cable 166  
output range 75  
output settings 75  
apg remote 187  
ASTM Drift 122  
automatic configuration with Bootp 209

## B

band width 6.5 nm 24, 27, 30  
BCD  
cable 171  
beam splitter 15  
Beer-Lambert  
absorbance 86

bench space 22  
boards  
photodiode boards (ADC) 15  
BootP service  
installation 211  
restart 218  
settings 218  
stop 217  
Bootp  
& Store 202  
automatic configuration 209  
initialization modes 201  
using default 203  
using stored 203

## C

cable  
analog 166  
BCD 171  
CAN 173  
connecting APG remote 42  
connecting CAN 42  
connecting LAN 42  
connecting the ChemStation 42  
connecting the power 42  
LAN 173  
remote 168  
RS-232 174  
cables  
analog 164  
BCD 164  
CAN 165  
LAN 165  
overview 164  
remote 164

RS-232 165  
CAN  
cable 173  
cautions and warnings 131  
cell test 118  
chromatogram 60  
cleaning 134  
communications 25, 28, 31  
compensation sensor open 102  
compensation sensor short 102  
condensation 22  
configuration switch  
location 200  
configuration  
one stack 37, 37, 39  
RFID tag (lamp and cell) 72  
stack 36  
temperature control 72  
two stack front 43  
two stack rear 44  
two stack 41, 43  
control and data evaluation 25, 28, 31  
correction factors for flow cells 87  
cutoff filter 14  
cuvette holder 143  
parts 160  
**D**  
data rate  
sampling rate 30, 27, 24  
data recovery  
DRC 80  
declaration of conformity 236  
defect on arrival 34

- delivery checklist 35
  - detection type 24, 27, 30
  - detector error messages 104
  - DHCP
    - general information 205
    - setup 206
  - diagnostic
    - signals 90
    - test functions 115
  - dimensions 23
  - DRC
    - run recovery 80
  - drift 24, 27, 30, 63
    - initial 68
- ## E
- electrical connections
    - descriptions of 179
  - EMF
    - early maintenance feedback 194
    - setting limits 195
    - using counters 195
  - entrance slit assembly 13
  - environment 22
  - error message
    - ADC hardware error 110
  - error messages
    - calibration failed 106
    - compensation sensor open 102
    - compensation sensor short 102
    - detector 104
    - fan failed 103
    - filter check failed 109
    - grating/filter motor defective 108
    - heater current missing 106
    - heater failed 111
    - heater power at limit 112
    - holmium oxide test failed 107
    - ignition without cover 103, 103
- ## F
- fan failed 103
  - features
    - GLP 31, 28, 25
    - safety and maintenance 31, 28, 25
  - firmware
    - description 176
    - main system 176
    - resident system 176
    - update tool 177
    - updates 177, 147, 147
    - upgrade/downgrade 147
    - upgrade/downgrade 147
  - flow cell
    - configuration 72
    - correction factors 87
    - high pressure (parts) 158
    - micro (parts) 154
    - semimicro (parts) 156
    - standard (parts) 152
    - types and data 30, 27, 24
    - with RFID tag 12
  - flow cells
    - matching for application 85
  - illegal temperature value from sensor
    - at air inlet 111
  - illegal value from sensor on main board 110
  - lamp current missing 104
  - lamp ignition failed 105
  - lamp voltage missing 105
  - leak sensor open 101
  - leak sensor short 101
  - leak 100
  - lost CAN partner 100
  - no run data available in device 112
  - remote timeout 99
  - shutdown 98
  - timeout 98
  - wavelength check failed 109
- ## G
- general error messages 98
  - grating
    - assembly 15
- ## H
- holmium oxide
    - declaration of conformity 236
    - filter 14
    - test 126
  - humidity 23
- ## I
- information
    - on cuvette holder 143
    - on solvents 234
    - on sound emission 232
    - on UV radiation 233
  - initialization mode selection 201
  - installation
    - bench space 22
    - of flow connections 52
    - of the detector 49
  - Instant Pilot G4208A 10
  - instrument layout 193
  - interfaces 182
  - internet 237
  - introduction 10
    - to optical system 11
- ## L
- lamp
    - configuration 72
    - initial drift 68
    - intensity test 116
    - type 30, 27, 24
    - with RFID tag 13

## Index

### LAN

- automatic configuration with
  - Bootp 209
  - Bootp & Store 202
  - Bootp 201
- cable 173
- configuration switches 200
- configuration 197
- initialization mode selection 201
- link configuration selection 208
- manual configuration with telnet 220
- manual configuration 219
- PC and User Interface Software Setup 225
- TCP/IP parameter configuration 199
  - using default 203
  - using stored 203
- what you have to do first 198
- leak sensor open 101
- leak sensor short 101
- leak
  - correcting 145
  - parts 161
- line frequency 23
- line voltage 23
- linearity 24, 27, 30
- link configuration selection 208
- lost CAN partner 100

### M

- MAC address
  - determine 214
- maintenance
  - correcting leaks 145
  - definition of 130
  - exchanging flow cell 138
  - exchanging lamps 135
  - exchanging leak handling system 146

- for parts see 'parts for maintenance' 149
- introduction 129
- overview 133
- replacing firmware 147, 147
- standard flow cell 140
- using the cuvette holder 143
- manual configuration
  - of LAN 219
- message
  - ADC hardware error 110
  - calibration failed 106
  - calibration lost 109
  - filter check failed 109
  - grating/filter motor defective 108
  - heater current missing 106
  - heater failed 111
  - heater power at limit 112
  - holmium oxide test failed 107
  - ignition without cover 103, 103
  - illegal temperature value from sensor at air inlet 111
  - illegal value from sensor on main board 110
  - lamp current missing 104
  - lamp ignition failed 105
  - lamp voltage missing 105
  - no run data available in device 112
  - remote timeout 99
  - wavelength check failed 109

### method

- load 64

### mirror

- assemblies 15

### N

- Noise and Linearity
  - specifications 32, 29, 26
- Noise Test 122
- noise, short term 24, 27, 30

- non-operating altitude 23
- non-operating temperature 23

### O

- online
  - plot 68
  - spectra 73
- operating Altitude 23
- operating temperature 23
- optical unit
  - beam splitter assembly 15
  - entrance slit assembly 13
  - filter assembly 14
  - filter 14
  - flow cell 12
  - grating assembly 15
  - lamp 13
  - mirrors 15
  - photodiode assemblies 15
  - photodiode boards 15
  - source lens assembly 13
  - stepper motor 15
- optimization
  - detector performance 84
  - of the system 61
- overview
  - optical path 11
  - optical system 11
  - system overview 11

### P

- packaging
  - damaged 34
- parameters
  - detector 65
- parts for maintenance
  - cuvette holder 160
  - high pressure flow cell 158
  - leak parts 161
  - micro flow cell 154

- semimicro flow cell 156
  - standard flow cell 152
  - parts
    - and materials for maintenance 149
  - PC and User Interface Software Setup 225
  - peakwidth
    - settings 76
  - performance
    - optimization 84
    - specifications 24
  - photodiode
    - assemblies 15
    - boards 15
  - photometric accuracy 87
  - physical specifications 23
  - physical
    - specifications 23
  - power consumption 23
  - power cords 21
  - power
    - considerations 20
  - preparing the HPLC system 61
- Q**
- quick noise test 123
- R**
- Radio Frequency Identification
    - flow cell and lamp 10
  - radio interference 231
  - recalibration of wavelength 90
  - remote
    - cable 168
  - repairs
    - cautions and warnings 131
    - introduction 129
    - overview of simple repairs 133
    - replacing firmware 147, 147
  - RFID
    - Radio Frequency Identification 10
  - RS-232C
    - cable 174
  - run recovery
    - automatic 81
    - manual 82
    - no run data available in 112
  - running the sample 70
- S**
- safety class I 229
  - safety
    - general information 229
    - standards 23
    - symbols 228
  - sample info 69
  - sampling rate
    - data rate 30, 27, 24
  - scanning 74
  - serial number
    - information 181, 181
  - setting up an analysis 57
  - settings
    - analog output settings 75
    - peakwidth 76
  - setup of detector 65
  - shutdown 98
  - signal
    - diagnostic 90
    - plot 67
  - site requirements
    - environment 22
    - power considerations 20
    - power cords 21
  - solvent information 234
  - sound emission 232
  - source lens assembly 13
  - special interfaces 188
  - special setpoints 76
  - special settings
    - boot-resident 191
    - forced cold start 192
  - specification
    - physical 23
  - specifications 26, 29, 32
    - noise and linearity 32, 29, 26
    - performance 24
    - physical 23
  - spectra
    - online 73
    - tools 31, 28, 25
  - stable baseline 63
  - stack configuration
    - front view 41
    - rear view 42
  - status indicators 90, 91
  - status indicator 92
  - stepper motor 15
  - stop-flow condition 73
- T**
- TCP/IP parameter configuration 199
  - telnet
    - configuration 220
  - temperature control
    - configuration 72
  - temperature sensor 100
  - test functions 90, 115
  - tests
    - available tests vs interface 93
    - holmium oxide 126
    - intensity of deuterium lamp 116
    - wavelength calibration 120
  - timeout 98
  - tool kit
    - hplc system 151
  - troubleshooting

## Index

- available tests vs interface 93
- diagnostic signals 90
- error messages 90, 97
- overview 90
- status indicators 90, 91
- test functions 115, 90

## U

- unpacking 34
- using
  - analog output settings 75
  - control settings 71
  - detector parameters 65
  - drift 63
  - EMF 194
  - load method 64
  - online plot 68
  - online spectra 73
  - peakwidth settings 76
  - preparing the HPLC system 61
  - priming and purging the system 57
  - requirements and conditions 59
  - running the sample 70
  - sample info 69
  - scanning 74
  - setting up an analysis 57
  - setup of detector 65
  - signal plot 67
  - special setpoints 76
  - special settings 71
  - stable baseline 63
  - stop-flow condition 73
  - the cuvette holder 143
  - the detector 55
  - turn on 62
  - typical chromatogram 60
  - warm up 63
- UV radiation 233

## V

- voltage range 23

## W

- warm-up
  - of detector 63
- warnings and cautions 131
- wavelength
  - accuracy 30, 27, 24
  - calibration 120
  - range 190-600 nm 30, 27, 24
  - recalibration 90
- weight 23



## In This Book

This manual contains technical reference information about the Agilent 1290 Infinity Variable Wavelength Detector (G1314E), the Agilent 1260 Infinity Variable Wavelength Detector (G1314F) and the Agilent 1200 series Variable Wavelength Detector (G1314D) (obsolete).

The manual describes the following:

- introduction and specifications,
- installation,
- using and optimizing,
- troubleshooting and diagnose,
- maintenance and repair,
- parts identification,
- hardware information,
- safety and related information.

© Agilent Technologies 2008, 2010-2012, 2013

Printed in Germany  
11/2013



G1314-90034 Rev. B