



Air

Fast separation of oxygen and nitrogen

Application Note

Energy & Fuels

Authors

Agilent Technologies, Inc.

Introduction

Gas chromatography with an Agilent CP-Molsieve 5Å column and Agilent 490 Micro GC separates oxygen and nitrogen in 40 seconds.



Agilent Technologies


Conditions

Technique : Micro-GC
GC-Channel : Agilent CP-Molsieve 5Å
Temperature : 100 °C
Carrier Gas : Helium, 26 psig
Injector : 100 ms, 110 °C
BackFlush : after 7 s
Detector : chip TCD

Courtesy : Jim Luong, Dow Chemical Canada

Peak identification

1. oxygen 21%
2. nitrogen 79%

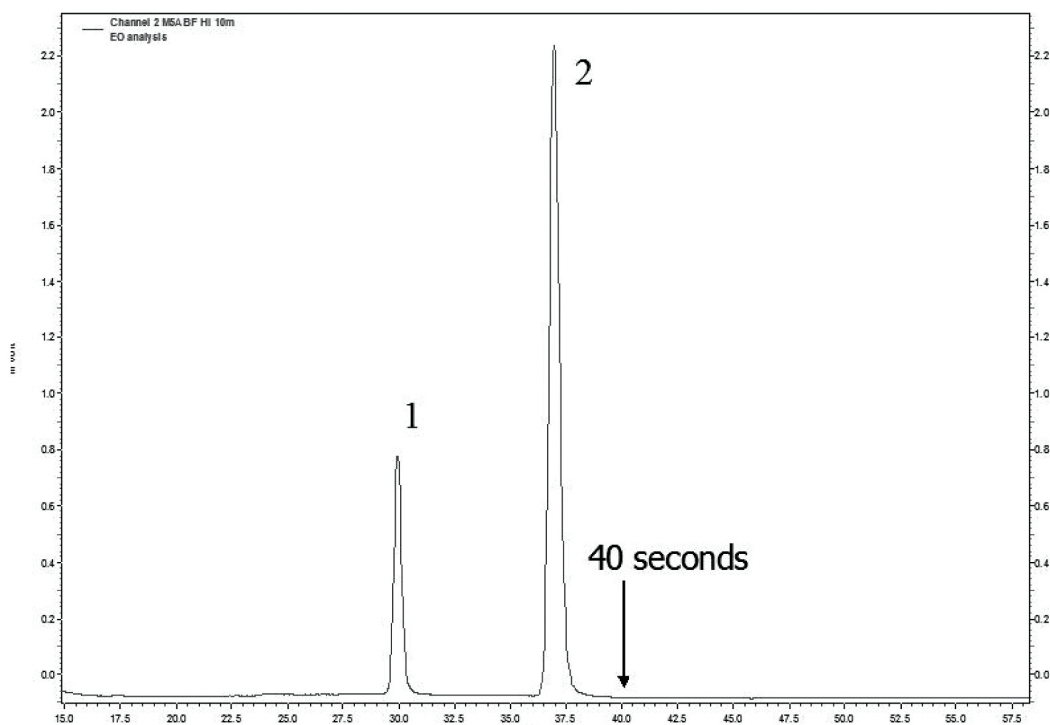

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A02061



Agilent Technologies